

SIERRA ANALYTICS, INC.

Version 3.0.0

PolymerixTM
Software
User Manual

VERSION 3.0.0 – MARCH 2013

Polymerix Software User Manual



© Copyright 2010 to 2013
Sierra Analytics, Inc.
5815 Stoddard Road, Suite 601
Modesto, CA 95356 USA
+1-209-545-8506

Polymerix@MassSpec.com
<http://www.MassSpec.com>

Table of Contents

Introduction	1
How this manual is organized	1
Some conventions used in this manual.....	1
Sierra's Web site	2
Getting Started	3
Goals for this chapter	3
Installing Polymerix	3
Registration.....	9
Installing the license key	16
What was covered in this chapter	19
In the next chapter	19
Polymerix Workflow Model and User Interface Overview	20
Goals for this chapter.....	20
The workflow model	20
Introducing the Polymerix user interface	22
What was covered in this chapter	27
In the next chapter	27
Data Import, Spectrum Selection, and Processing	28
Goals for this chapter.....	28
Importing data.....	29
Restricting the processing mass range	33
Detecting peaks	35
Mass adjustment.....	38
Mass mode	40
Deisotoping	41
Saving and loading parameter template files	46
What was covered in this chapter	47
In the next chapter	48
Find Homopolymer Series.....	49
Goals for this chapter	49
Overview	49

Simplest use case example	50
Understanding reference peaks	51
How Polymerix handles reference peak selection.....	52
Automatic reference peak selection	52
Find Homopolymer Series page.....	53
Selecting the repeat unit	54
Selecting end groups and adducts	55
Setting constraints on chemical composition	56
Troubleshooting Tips.....	59
What was covered in this chapter	59
In the next chapter	59
Homopolymer Analysis.....	60
Goals for this chapter	60
Molecular weight distribution statistics	60
The Homopolymer Analysis page user interface.....	62
Setting homopolymer processing parameters.....	63
Defining homopolymer series.....	65
Homopolymer molecular weight statistics	69
Visualizing homopolymer series assignments.....	70
Assignments	70
Filtered Assignments	70
Filtered Residuals	71
Details	71
Mass Ranges	71
Distribution Plots.....	72
Assignment and matching details	73
The simple matching algorithm.....	73
Cluster matching.....	73
Centroid cluster matching	74
Centroid cluster matching algorithm	74
What was covered in this chapter	74
In the next chapter	75
Copolymer Analysis	76
Goals for this chapter	76
Copolymer definition	76
The Copolymer Analysis page user interface.....	77
Setting copolymer processing parameters	78
Defining copolymer series.....	79
Copolymer molecular weight statistics	80

Visualizing copolymer series assignments	80
Assignments	80
Filtered Assignments	81
Filtered Residuals	81
Details	82
Mass Ranges	82
AB Distribution	83
Repeat A Statistics.....	84
Repeat B Statistics.....	85
Repeat A Plot.....	85
Repeat B Plot.....	86
What was covered in this chapter	86
In the next chapter	87
The Relative Area Tool.....	88
Goals for this chapter	88
The Relative Area page	88
Specifying ranges	89
Relative Area results	89
In the next chapter	90
Autocorrelation	91
Goals for this chapter	91
Accessing the Autocorrelation page.....	91
The autocorrelation spectrum	91
The end group spectrum	93
User interface and parameters.....	94
What was covered in this chapter	97
In the next chapter	97
Creating and Customizing Output	98
Goals for this chapter	98
Polymerix options.....	98
Copying to the clipboard	100
Saving and exporting results.....	101
Printing.....	103
What was covered in this chapter	106
In the next chapter	106
Release Notes.....	107
Version 3.0.....	107
Major new features	107
Updates, bug fixes, and new features	107

Version 2.0	108
Updates, bug fixes, and new features	108
Example Report	110

Introduction

Some tips on using this manual effectively

Welcome to Polymerix. We're looking forward to helping you get the most out of your data with a minimum of headache. But before getting started, we would like to explain some of the conventions we use in this manual which we hope will make it easier for you to understand how to use the manual and software to maximum effectiveness.

How this manual is organized

Our goal in writing this manual is to guide you from initial installation and registration to expert use of Polymerix by presenting the material in a way that models your workflow. Instead of a programmer's point of view ("Here's the **File** menu. It has five submenus..."), we try to take a task-oriented approach, in which each major task is broken into a series of smaller subtasks. We use a minimum of explanation and a maximum of illustration to help make the tasks as clear and easy to follow as possible.

So, the first part of this manual is a description of the tasks to get you started in using Polymerix. In many cases, the steps will be sufficient to solve most problems. However, Polymerix is complex software, with many parameters and options to handle a wide range of sample types and analytical approaches. The second part of the manual goes into much greater detail to describe this complexity, explaining each setting and its effect on the analysis. Finally, because no single software product can accomplish all things for all people, the last part of the manual details the export format used to store all of Polymerix's parameters and results and gives some examples of how to extract information for use in other software.

Some conventions used in this manual

We have standardized on certain typographical conventions to help distinguish normal text from labels, prompts, or other text you see onscreen, to highlight important items, and to point out areas where you must pay attention to avoid mistakes that could affect the quality or accuracy of your results.

INTRODUCTION

- Text that appears in **boldface** normal font is used to display menu items, labels, or other prompts that appear onscreen.
- Text that appears in **boldface Arial font** is used to emphasize some point in this manual.
- An icon is displayed to indicate an important point or warning:
 - ⇒ An arrow sets off an item of importance.
 - ! An exclamation point indicates a warning.
 - ⌚ An hourglass indicates a step that might require lengthy processing.
- In the later parts of the manual, text in `computer font` is used to present items from Polymerix's parameter and results formats, or to indicate text you would write in scripting files used to extract and process these formats.

ICON KEY

⇒ Valuable information

! Warning

⌚ Lengthy process

Sierra's Web site

Visit our Web site at <http://www.MassSpec.com> for more Polymerix information, application notes, publications, and other useful information. We plan to update the site frequently with new content.

Getting Started

How to install Polymerix, submit registration details, and install a license key

Polymerix is a registered and licensed product. Depending on the details of your purchase, you may have a limited or a perpetual term license. In either case, we will issue a license key valid for the term of your license which must be installed on your PC before the software will function. At the end of a limited license term, the software will cease to run until you renew your license and obtain a new key.

Goals for this chapter

In this chapter, we will explain how to:

- Install Polymerix from the distribution media or through a Web download
- Register the software and obtain a license key
- Install the license key and begin using the software

Installing Polymerix

We deliver Polymerix software in one of two ways: physically, on a CD-ROM, or electronically via an Internet download from Sierra's Web site.


Starting the installation program

If you received Polymerix on a CD-ROM, insert the disk into your CD-ROM drive. The installation program should start automatically. If it does not start, browse to the CD-ROM folder and double-click the **Polymerix_3.0_FullSetup** EXE file to start the installation program.

If you received an e-mail with a Web link, double-click the link in the e-mail (or copy the link to your Web browser if your e-mail program does not permit opening links), and follow the instructions on the Web page to download an appropriate setup program, which differs depending upon the data importers which are provided. Save


INSTALLATION, LICENSING, AND REGISTRATION

the downloaded file to your disk, unzip the distribution, and then double-click the **Polymerix_3.0_FullSetup.exe** file (or the **Polymerix_3.0_*Setup.exe** file, where * represents a vendor name) to start the installation program.

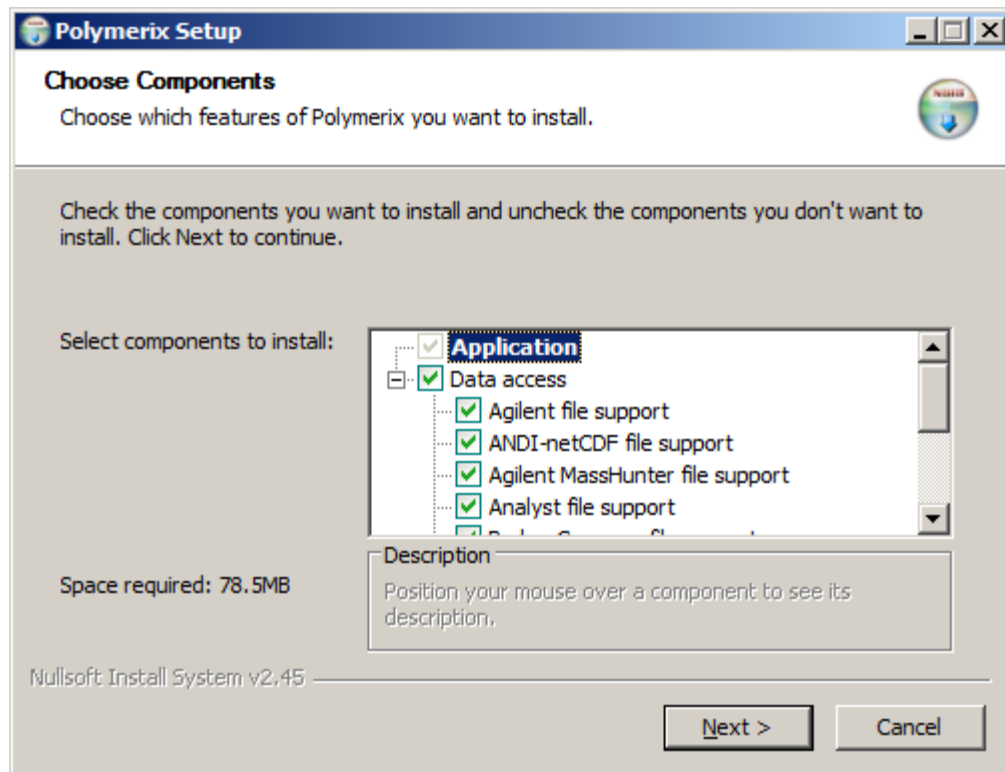


The installation program is a dialog-based “wizard” that guides you through the setup steps

1. The first page of the setup wizard is shown above. The item labeled **Application** is already selected. In order to be able to read raw data directly from mass spectrometry vendor format files, you must also select one or more of the formats listed under the **Data access** section. In this example, the full set of supported formats is available and has been selected. Depending on the details of your purchase, you may have a different set of data access selections. If your copy of Polymerix was purchased through an MS instrument vendor, only that vendor’s formats will be available.

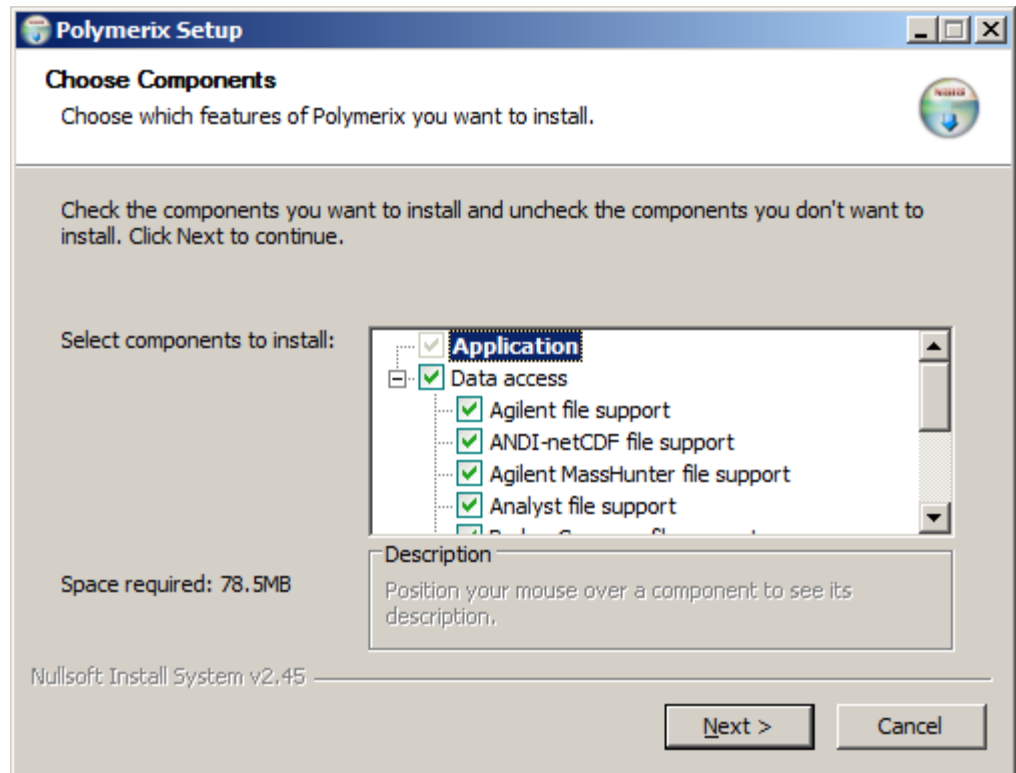


2. Select the desired data access option(s), then click the **Next >** button. At any time, you may click the **Cancel** button to stop the installation and remove any installed files from your computer.


You **must** select at least one of the data access items if you wish to read raw data from vendor format files.

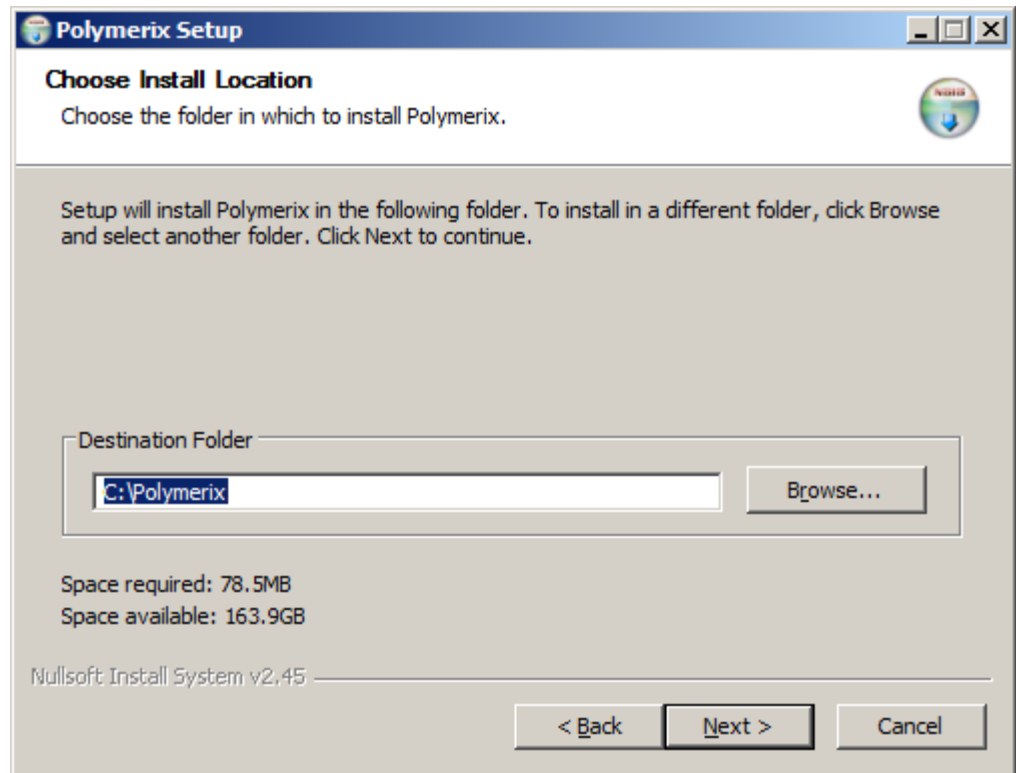


3. Next, choose the installation location. By default, Polymerix will be installed in the folder **C:\Polymerix**. If this folder does not exist, it will be created. If you wish to install Polymerix in a different folder, click the **Browse...** button to choose an alternative. Click **Next >** to continue.

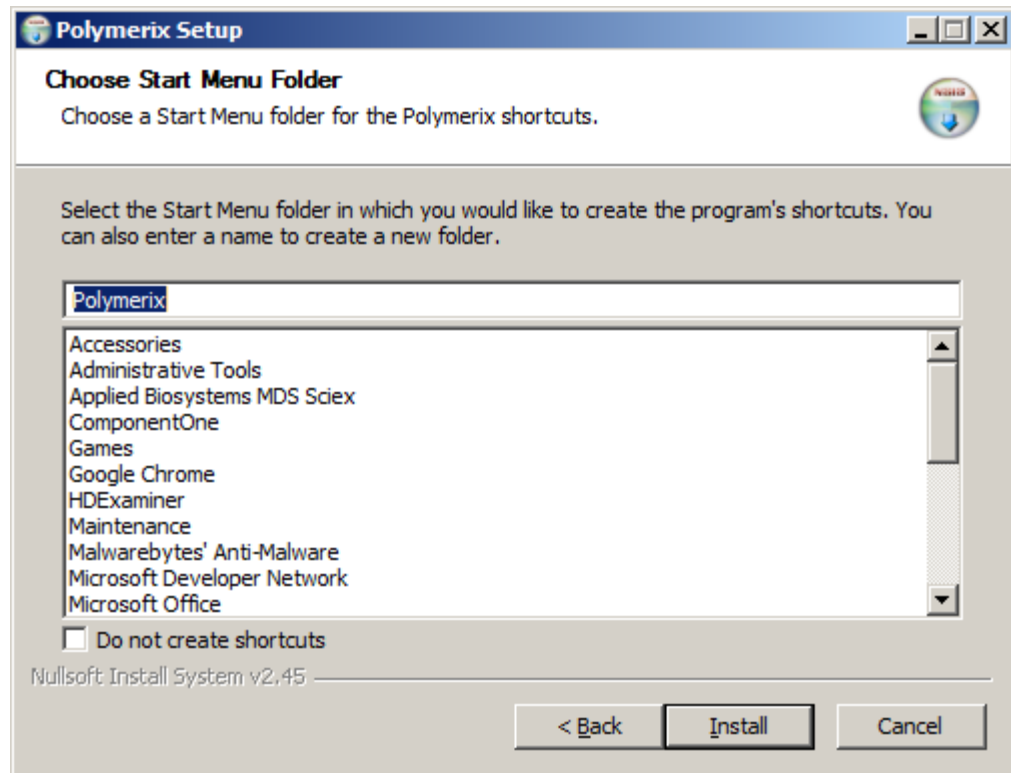

4. Choose the name for the Start Menu folder. If you do not wish to have an entry made on the Start Menu and a Polymerix icon placed on your desktop, select the **Do not create shortcuts** option. Click **Install** to begin copying files to the installation location.

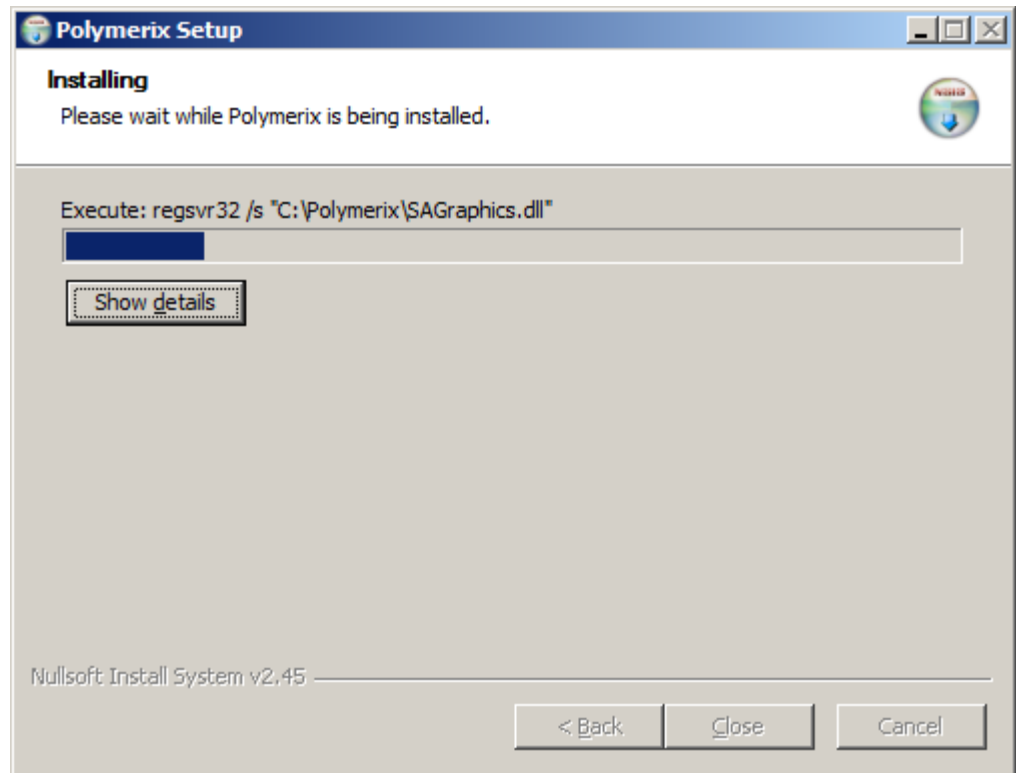

5. As the installation program extracts and copies files to the destination folder, the progress is displayed. To see details of this process, click the **Show details** button.

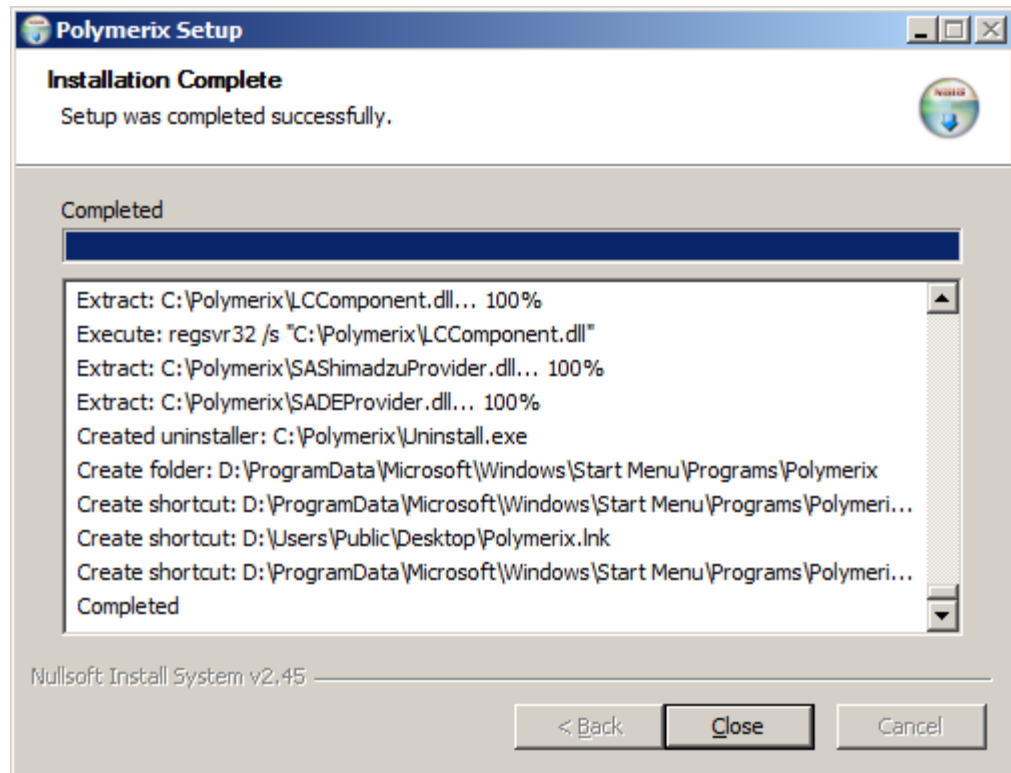

- Once installation has completed, click the **Close** button to end the setup program. This screenshot shows the installation wizard after the **Show details** button was clicked. The details shown for your installation might be different.

Registration

After installation, the next step is to register the software and submit a request for a license key. Registration is required. The software will not function until registration has been completed and a valid license key has been entered.

Two options for licensing

You may license your installation of Polymerix using either a fixed license or a floating license.

- A fixed license locks your copy of Polymerix to a specific PC. The registration procedure reads information from the PC hardware creates a license key request. Using this request and the term of your license, we create a license key that permits Polymerix to run only on the chosen PC.
- A floating license allows installation and registration of Polymerix on up to 4 PCs, using a combination of each PC's hardware information and a USB memory stick. Once licensed, the software may be run on any of the PCs, but only one at a time. The USB memory stick **must** be inserted at all times while

the software is in use. If the USB device is removed, the software will cease working until the device is reinserted.



You must choose a license type when you first register Polymerix. Once you have submitted a registration request and have installed a license key, we may not permit you to change the license type until the license is renewed at the end of the license term. We may issue a new license key under special circumstances, such as a lost USB memory stick or an upgrade to a new PC, but we may require that you first un-register the software and send us proof via e-mail.

A summary of the registration process

These are the steps involved in registration:

1. Run the **Registration Wizard** and enter information used to identify the type of license, USB device if applicable, and your contact details.
2. E-mail the registration information to us. The Wizard will automatically create the e-mail for you.
3. Wait for our e-mail in reply. It will contain your license key.


Using the Registration Wizard to request a license key

When you start Polymerix for the first time, it detects that the software is not registered and invokes the Registration Wizard. The wizard is a dialog-based procedure that guides you stepwise through the registration process.




1. Double-click the Polymerix icon on your desktop or select **Programs → Polymerix → Polymerix** from the **Start** menu to start the Polymerix application. The Polymerix “splash screen” will appear for a few seconds, and then the first page of the Registration Wizard will be shown.


INSTALLATION, LICENSING, AND REGISTRATION

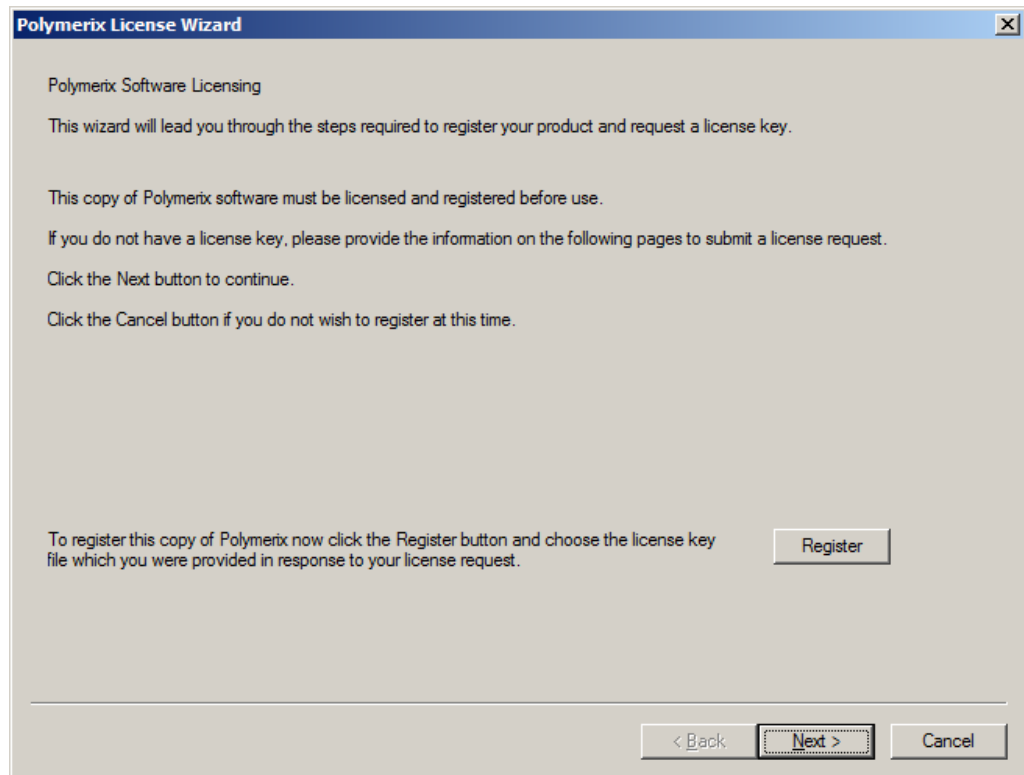

2. Since you do not yet have a license key, click the **Next >** button to continue.

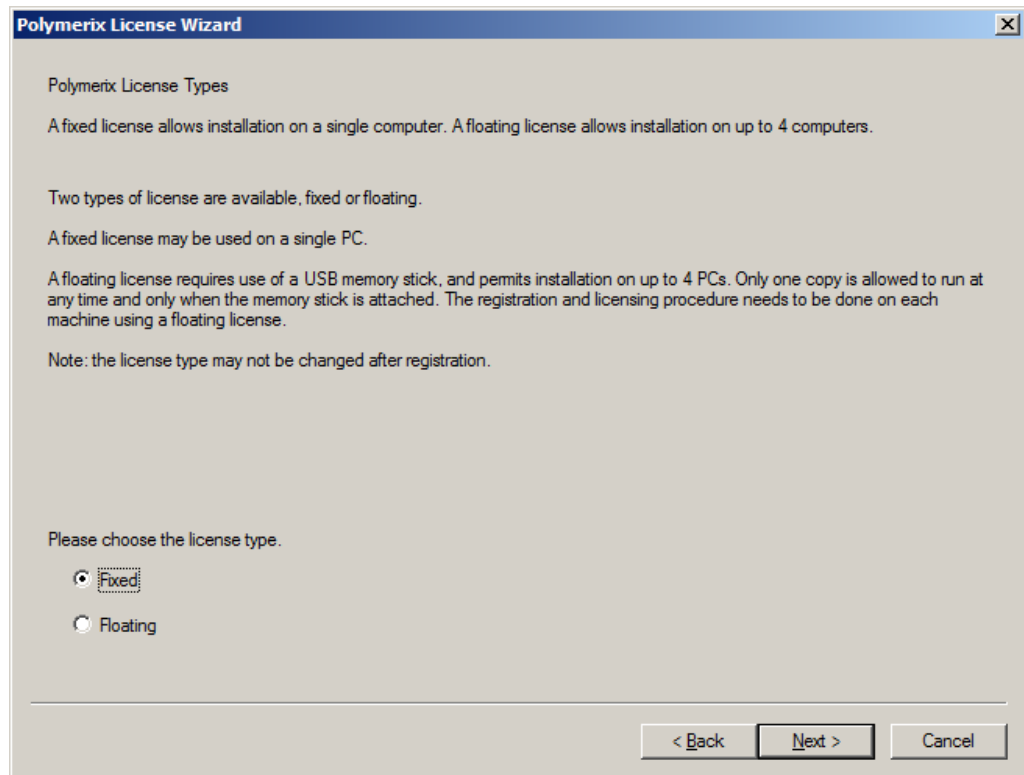

3. On this second page you choose the type of license, fixed or floating. If you wish to lock the software to a single PC, choose the **Fixed** option, otherwise choose the **Floating** option. A floating license permits installation on up to 4 PCs, but requires use of a USB device that you can transfer from PC to PC when you wish to run Polymerix. The USB device must be installed in the PC during the entire duration of your Polymerix session. Usually, a USB “memory stick” is the most convenient way to do this. Any kind of USB device may be used, as long as it has a unique name and serial number that can be detected by the software. Click the **Next >** button to continue.

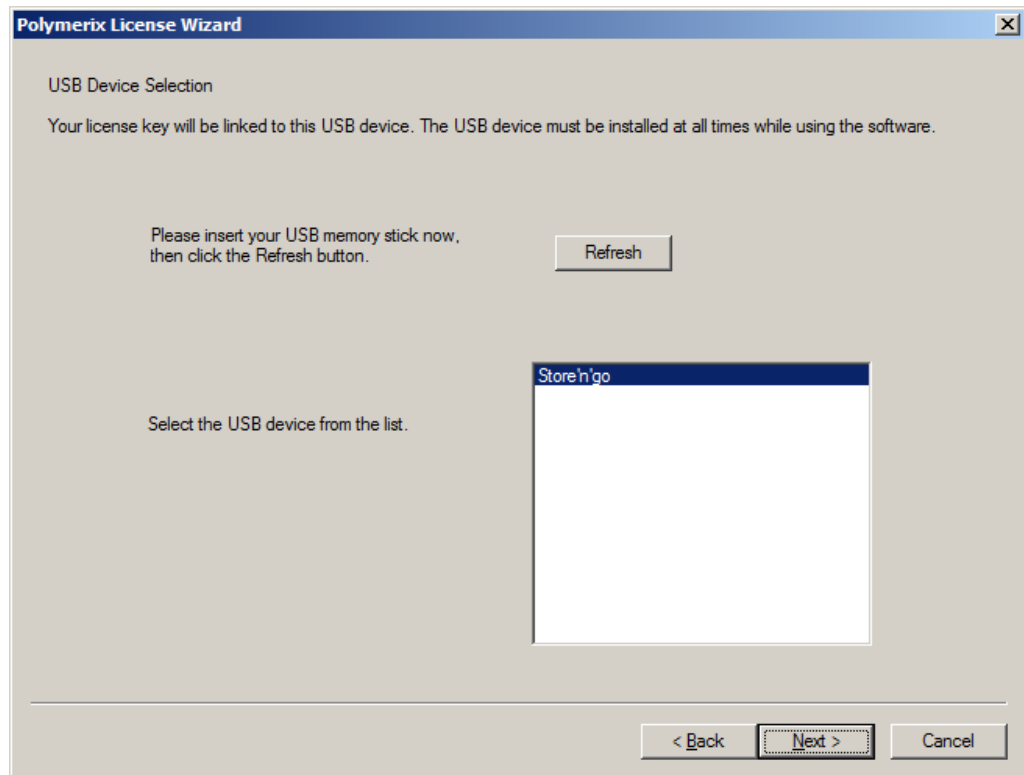

4. If you chose the floating license option, on this page you will select the USB device used for the license key. If the USB device is not inserted, install it now, and then click the **Refresh** button until it appears in the list of available devices. **Note: If you are registering the software on more than one PC, you must use the same USB device on each one.** We will not issue license keys based on different USB devices if you have only purchased a single license to the software. Select the USB device from the list and click the **Next >** button to continue.

INSTALLATION, LICENSING, AND REGISTRATION

Polymerix License Wizard

License Registration Information

This information is required to create a license key for your product.

Please provide the information below. Items labeled with * are required.

Name *	<input type="text" value="Ignatius Jaques Reilly"/>		
E-mail *	<input type="text" value="Ignatius@MassSpec.com"/>	Telephone *	<input type="text" value="504-555-5555"/>
Institution *	<input type="text" value="Sierra Analytics"/>		
Street	<input type="text"/>		
	<input type="text"/>		
City	<input type="text"/>	State / Region	<input type="text"/>
Postal Code	<input type="text"/>	Country	<input type="text"/>

This information will be used only to register the software and to enable us to provide software support. Will will not make this information available to third parties, and will not use it to send spam email.

< Back Next > Cancel

5. Please enter your contact details on this page. We would appreciate complete details, including mailing address. We collect this information so that we can provide personalized support and updates for your software. We do not share this information with anyone. Click the **Next >** button to continue.

INSTALLATION, LICENSING, AND REGISTRATION

The screenshot shows a window titled "Polymerix License Wizard" with a close button in the top right corner. The main content area is titled "License Request Submission" and contains the following text:

License Request Submission

License request submission consists of saving registration information to a license request file

Please check this information carefully. Use the Back button to make any changes.


The information to be submitted is displayed in a text box:

```
License type : fixed
User name : Ignatius Jaques Reilly
Institution : Sierra Analytics
Address 1 :
Address 2 :
City :
State / Region :
Country :
Postal code :
E-mail : Ignatius@MassSpec.com
Telephone : 504-555-5555
```

Below the text box, it says: "To create the license request file click the Save button."

At the bottom of the dialog, there are three buttons: "< Back", "Next >", and "Cancel". The "Next >" button is highlighted with a dashed border.

6. This final page shows the information that will be submitted as your license request. Please review it carefully. If you need to make changes, click the < **Back** button to go to the appropriate page. If everything is correct, click the **Save...** button to create the registration request file.

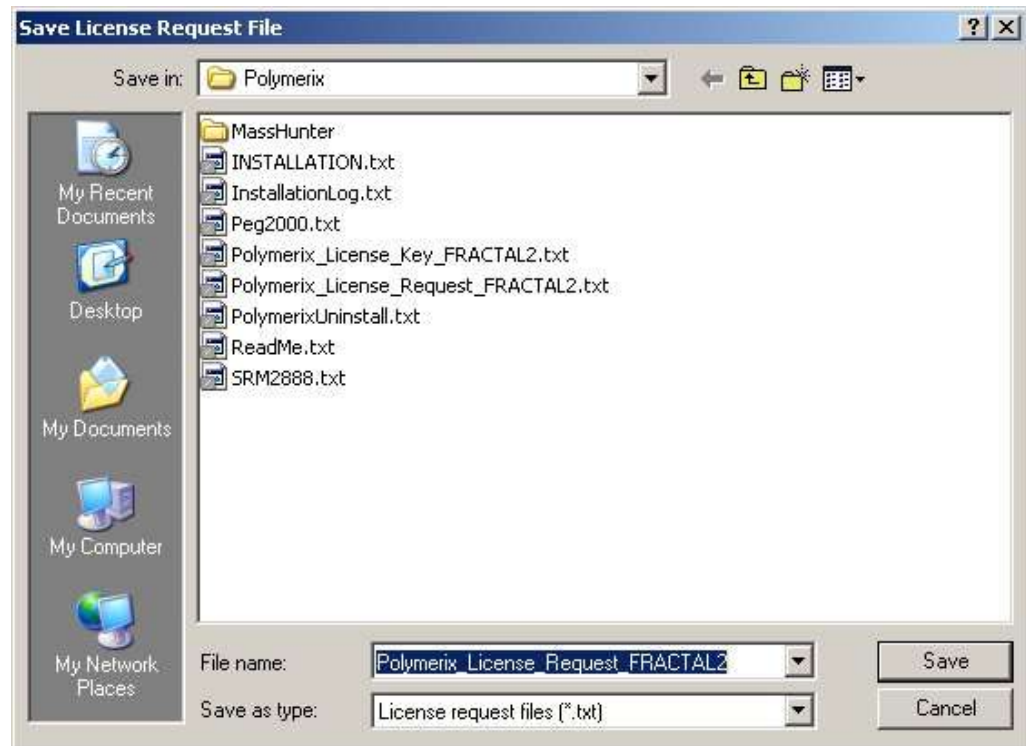

7. Once the license request file is saved, continue with the Polymerix redeem download instructions. They will allow you to retrieve a license key file, which you will use in the next section.

Installing the license key

If registration was successful, we will send you a license key by reply to your e-mail. You will use the Registration Wizard again, this time to copy the license key into the software and license the copy.

A summary of the license installation process

These are the steps involved in installing your license key:


1. Open the saved file with the license key.
2. Run the **Registration Wizard** and load the license key file into the wizard.
3. Start using Polymerix!

Using the Registration Wizard to install a license key


Until you have installed a valid license key, each time you start the Polymerix software it will run the Registration Wizard.



1. Double-click the Polymerix icon on your desktop or select **Programs** → **Polymerix** → **Polymerix** from the **Start** menu to start the Polymerix application. The Polymerix “splash screen” will appear for a few seconds, and then the first page of the Registration Wizard will be shown.

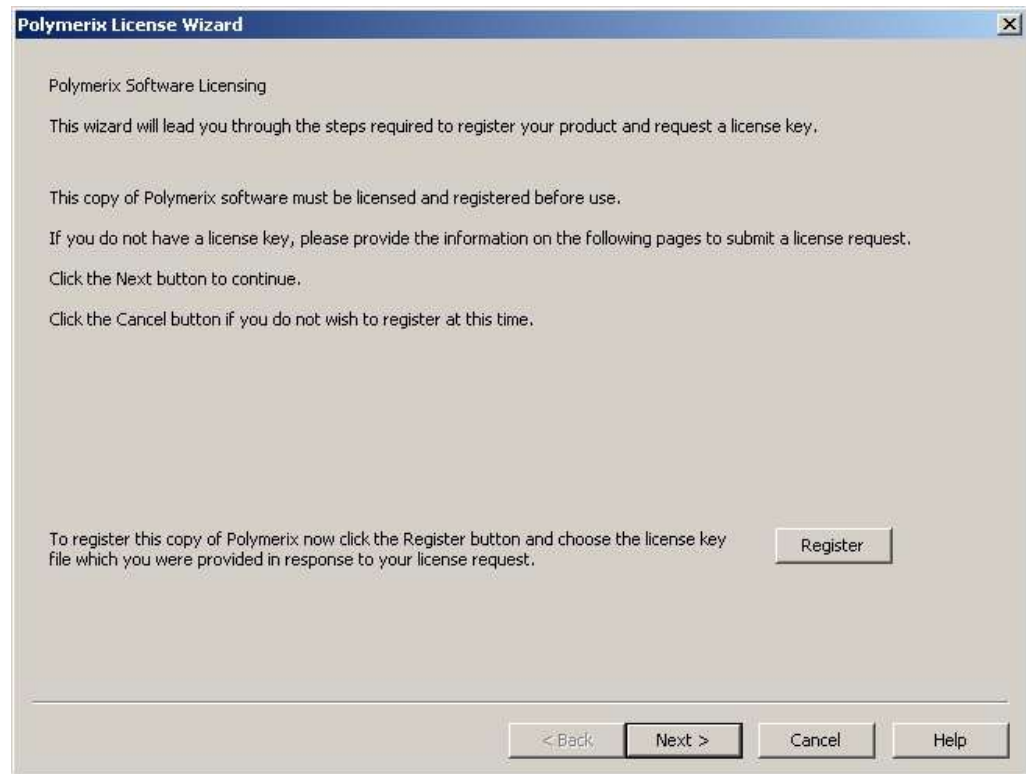

2. On the first page of the wizard, click the **Register** button. This will open a file selection dialog. Choose the registration key file you saved from online registration process..

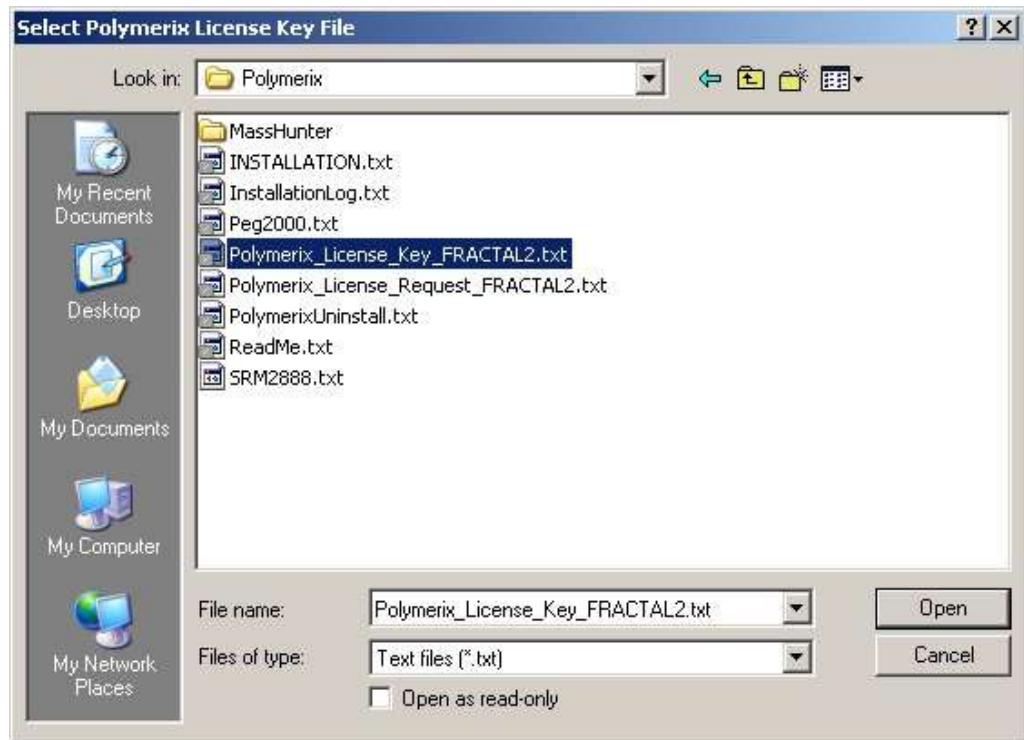


3. Select the license key file and click the **Open** button. The Registration Wizard will read the file contents and verify them.



4. If everything is OK, the “Thank you” dialog will be displayed. Click the **OK** button to close this dialog, then the **Finish** button on the Registration Wizard.
5. The main Polymerix page will then be displayed. If the registration failed for any reason, a message will be displayed with that information, and the registration wizard will remain open. If you used a USB memory device for the registration request, **be sure it is inserted** and recognized by your PC. If everything seems OK, but the licensing still fails, please contact us at **Polymerix@MassSpec.com** and we will try to work through the problem with you.
6. Please note that registration keys are specific to the PC that was used to generate the license request. You cannot use the same file to license Polymerix on a different PC. If you are using a floating registration, **you**

must create a registration request and receive a license key file for each PC, using the same USB device each time.

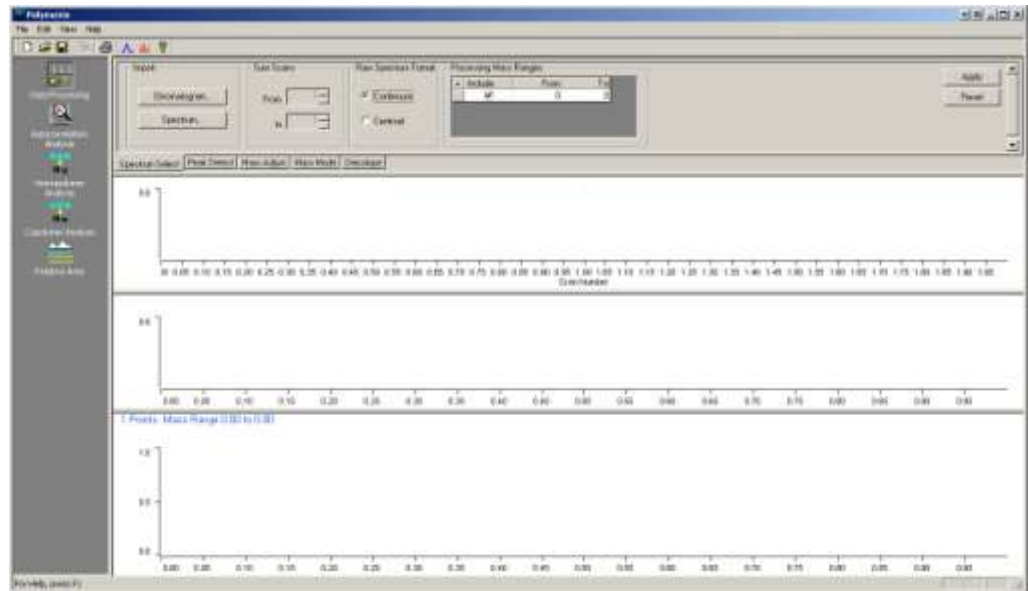

7. Time to get to some real work!

What was covered in this chapter

In this chapter, we described how to:

- Install the Polymerix software from the distribution.
- Create a registration request and submit it by e-mail to obtain a license key.
- Install the license key in the software

In the next chapter

We will describe the workflow model used in the Polymerix software, and show how that workflow is implemented in the main parts of the Polymerix user interface.

Polymerix Workflow Model and User Interface Overview

How Polymerix models the data reduction, analysis, and reporting workflow and the principal elements of the user interface

Polymerix's workflow model is designed to simulate the same processes and steps you would use if you were to perform this data analysis and reduction manually. The user interface implements this workflow model in an intuitive and graphical fashion.

Goals for this chapter

In this chapter we will cover the following material:

- We will describe the model used in Polymerix for the data reduction, analysis, and reporting workflow.
- We will present the major parts of the Polymerix user interface and how they fit the workflow model.

The workflow model

A typical polymeric or related sample consists of one or more homologous series, each consisting of one (for a *homopolymer*) or two (for a *copolymer*) repeating units (the *monomers*), with *alpha* and *omega* (or *total*) *end groups*. When analyzed by mass spectrometry, the resulting spectrum may likewise show spectral features for each of these series and possibly more, dependent on the sampling and ionization method and resolving power of the instrument. The goals of polymer analysis are to assign a unique chemical composition and abundance to as many of the detected features as possible, and to compute the molecular weight distribution statistics for each series.

Goals of polymer analysis

Steps in the workflow

The workflow for this analytical process can be broken down into a series of dependent steps:

1. Data import and reduction of the profile (continuum) mode spectrum to an m/z and abundance peak list, followed by optional mass adjustment of the peak list to assign highly accurate m/z values and/or deisotoping to merge higher isotopomers into the monoisotopic peak.
2. Creation of one or more homologous series definitions to permit assignment of the sets of peaks in the spectrum.
3. Visualization of sample composition and series assignments through the use of graphs and tables.
4. Creation of a report of the results of the analysis.

Each of these steps depends on successful completion of the previous step. Failure to accurately peak detect the profile spectrum can lead to missed peaks or false additional peaks, which will result in the failure of subsequent steps to produce correct results. Likewise, bad mass assignment will result in inaccurate correction of peak m/z values, which in turn leads to incorrect or missed series assignments, and so forth.

**Mapping workflow
to the user
interface**

Mapping this workflow onto an intuitive user interface is a major goal of the software. The user interface must permit specification of the relevant parameters for each processing stage, and then allow browsing the processing results to validate their correctness. Corresponding to the workflow described above, the user interface is organized into pages each of which corresponds to a step in the workflow:


1. A data reduction and processing page, which displays the total ion chromatogram (TIC) if present, the input profile spectrum, and the fully processed spectrum that will be used for series assignment.
2. A page with graphical tools for helping to determine from spectral data the chemical compositions of the units which construct a homopolymer. The resultant chemical compositions can be transferred to the homopolymer analysis page.
3. A homopolymer analysis page used to view the quantitative results of performing homopolymer series assignments. Homopolymer series definitions can be specified here.
4. A copolymer analysis page, used to setup copolymer series definitions and to view the assignments and decomposition into spectral subsets by repeat unit count.
5. A relative area page, where relative areas of multimodal distributions can be selected and measured.
6. An optional autocorrelation page, which can assist in determining repeat units and total end groups for an unknown polymer. The functionality of

this page has largely been superseded by the Find Homopolymer Series page in item two above.

In support of these pages are many dialogs and other user interface items for data import and export, setting parameters, displaying tables and graphics, and switching between the various pages and data views.


Introducing the Polymerix user interface

The general layout of the Polymerix user interface is similar to that used in many popular software applications. There are menu and tool bars, a set of main workspace pages, and a control panel for rapid switching between these pages. These are indicated in the screenshot below.

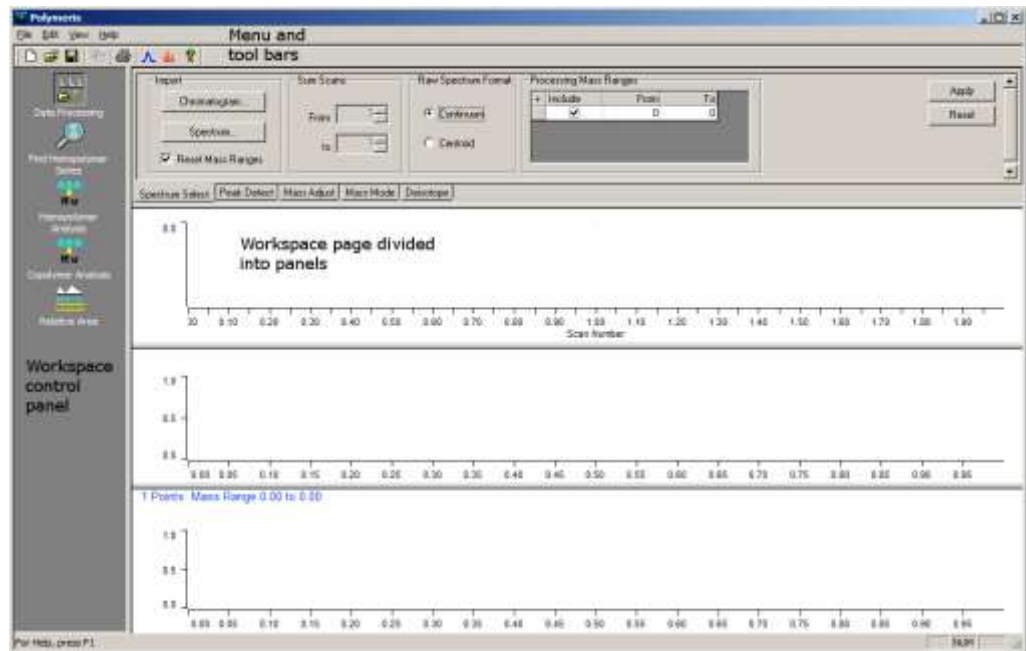

Processing Page

The buttons in the workspace control panel are used to switch between the top-level pages on the right. In almost all cases, one of the first steps in your workflow will be to import raw data, perform peak detection and other processing, and review the results to ensure that the processing is correct. The Processing page, shown below, is used for that purpose.

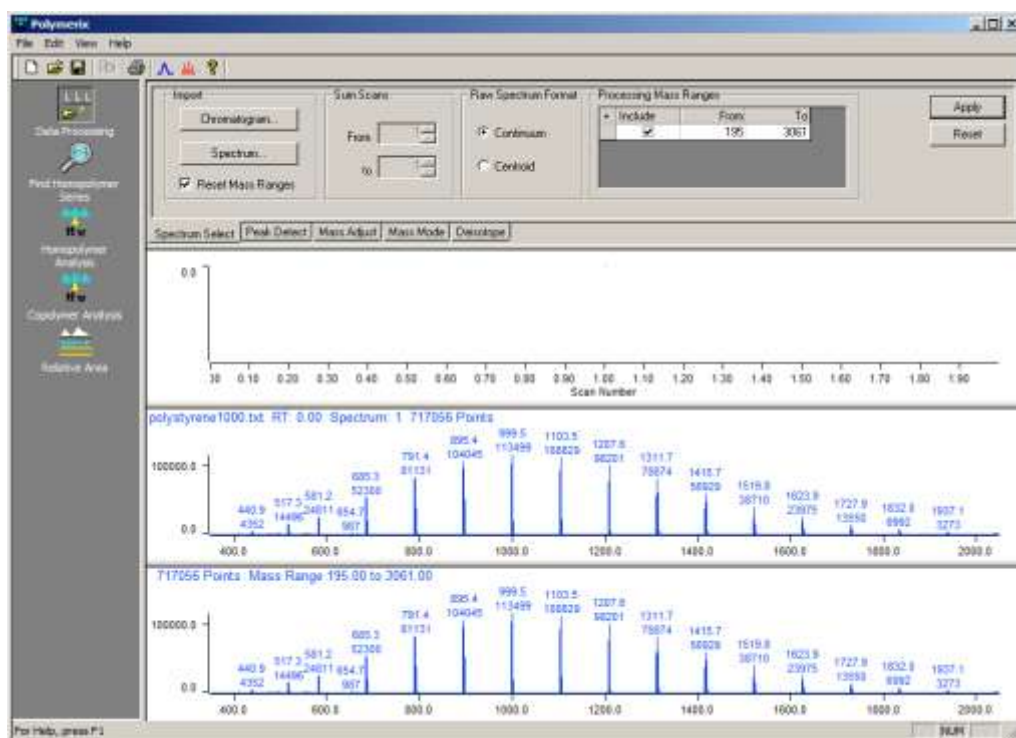


The topmost pane uses tabs to choose between different data import and processing parameter specification editors. In the second pane is the TIC (or equivalent for multi-scan infusions or other sampling). The third pane shows the raw data as imported, prior to any processing. In the bottom pane shows the final processed spectrum after all processing parameters have been applied. The m/z axes of the spectrum plots are linked, and zooming or panning in one by changes the other to the same axis range.

Autocorrelation Page

The **Autocorrelation Analysis** button is not displayed by default, because **Find Homopolymer Series** (see below) supersedes the functionality of autocorrelation analysis in most cases. To enable and show the button, select **View Autocorrelation** from the Menu bar and then exit and restart Polymerix..

Clicking the **Autocorrelation Analysis** button on the control panel changes the right side to the **Autocorrelation Page**. For an unknown polymer, autocorrelation analysis may provide insights into the nature of the repeat units and total end group. This page provides a way to review the autocorrelation and end group spectra and manipulate their settings.

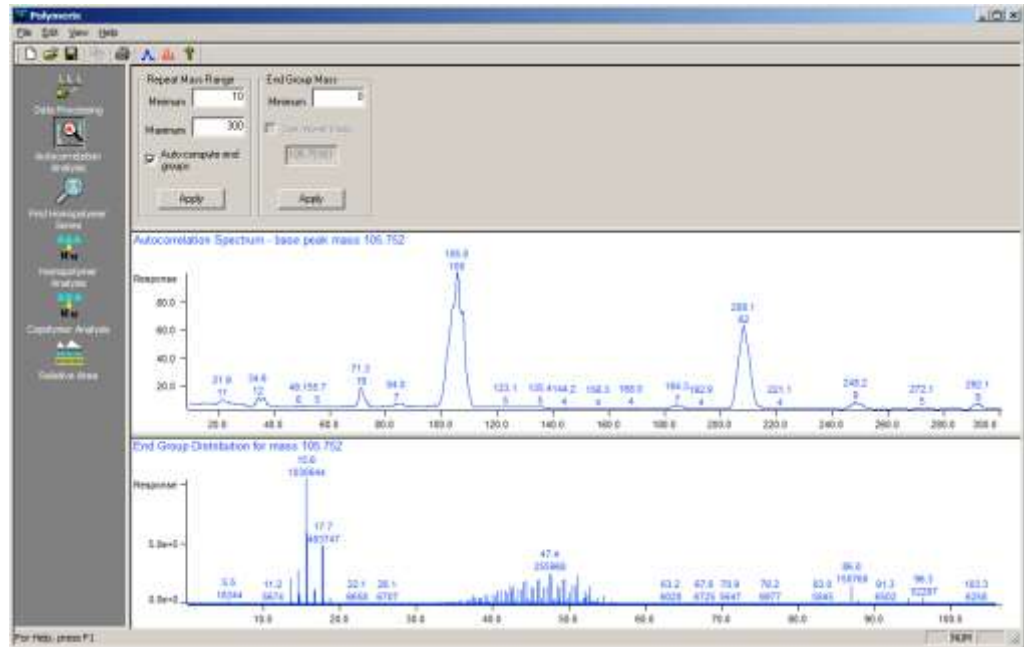


In the topmost pane the autocorrelation and end group parameters may be adjusted. In the middle pane the autocorrelation spectrum is displayed. In the bottommost pane the end group spectrum is shown. Each spectrum is manipulated independently.

Find Homopolymer Series Page

Clicking the **Find Homopolymer Series** button on the control panel changes the right side to the **Find Homopolymer Series** Page. This page assists the user in determining the chemical formulas of a centroided spectrum's homopolymer repeat units, end groups, and adducts. Once determined the chemical formulas can be transferred to the **Homopolymer Analysis** Page for quantitative analysis.


POLYMERIX WORKFLOW AND USER INTERFACE

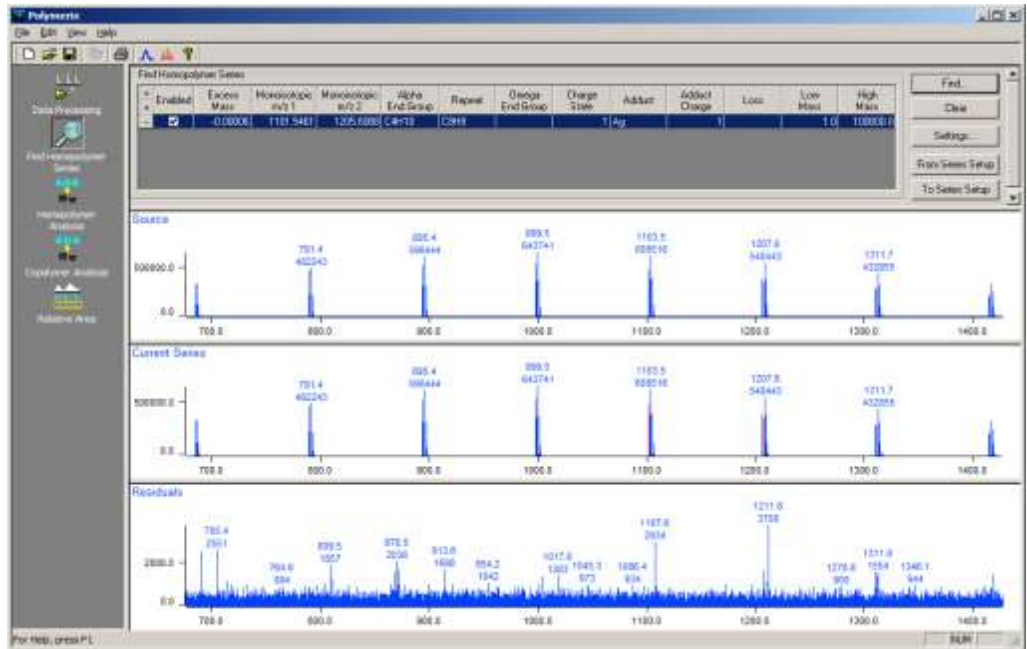

In the topmost pane, series parameters and setup results are displayed. The top spectrum plot shows the spectrum being analyzed. The middle spectrum plot shows a selected series. The bottom plot displays the residual spectrum – in other words, all peaks still not assigned to a series.

Homopolymer Analysis Page

Clicking the **Homopolymer Analysis** button on the workspace control panel switches the right-hand side to the **Homopolymer Analysis** page.

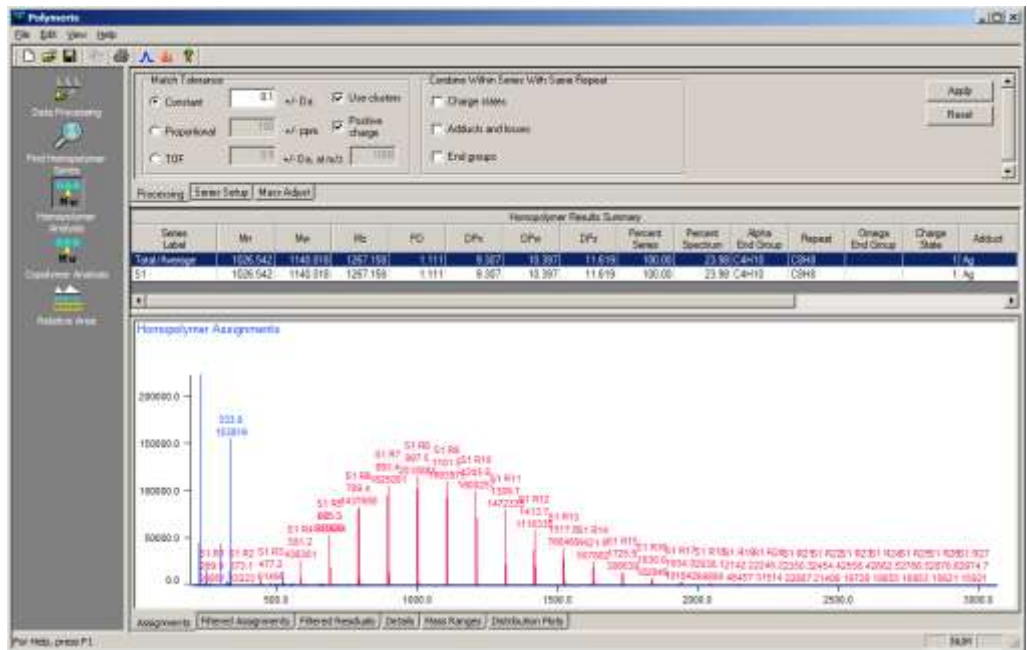


POLYMERIX WORKFLOW AND USER INTERFACE

The top pane of this page contains tabbed views for defining homopolymer series and editing peak matching parameters. The middle pane contains a table of the polymer series molecular weight distribution statistics and relative amounts, and the bottom pane contains a series of tabbed views that display the assignment results in graphical or tabular form.


Copolymer Analysis Page

Clicking the **Copolymer Analysis** button on the workspace control panel switches the right-hand side to the **Copolymer Analysis Page**. It is very similar to the **Homopolymer Analysis** page, but with additional features specific to the two-dimensional nature of copolymer distributions.

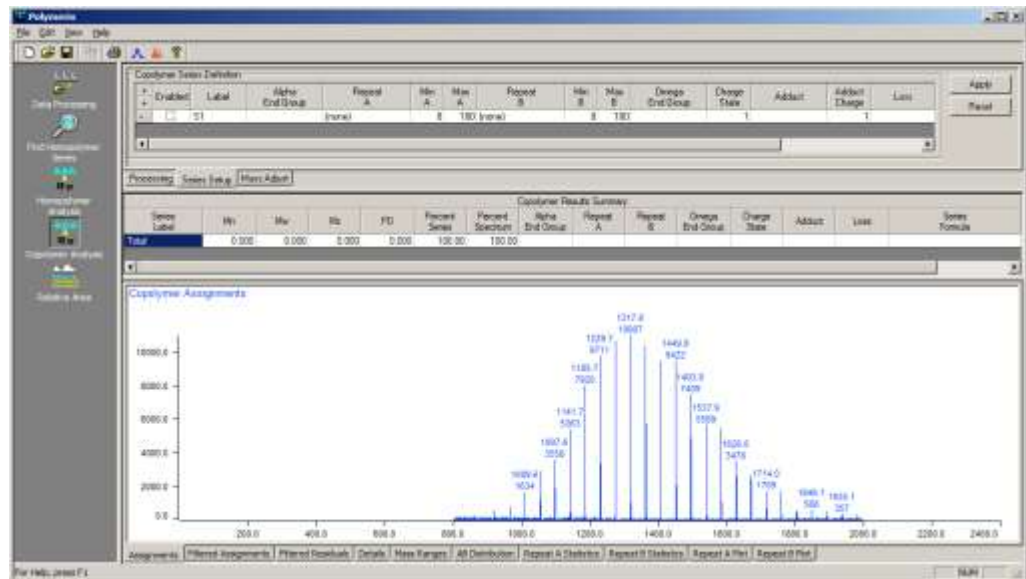

Clicking the **Relative Area** button on the workspace control panel switches the right-hand side to the **Relative Area Page**. This page is used to subdivide the spectrum into two or more regions to measure relative percentages of spectral intensity in each region. This is useful if the polymer has a multimodal distribution or if there are multiple charge states, and you are interested in the relative amount of polymer in each distribution.

Relative Area Page

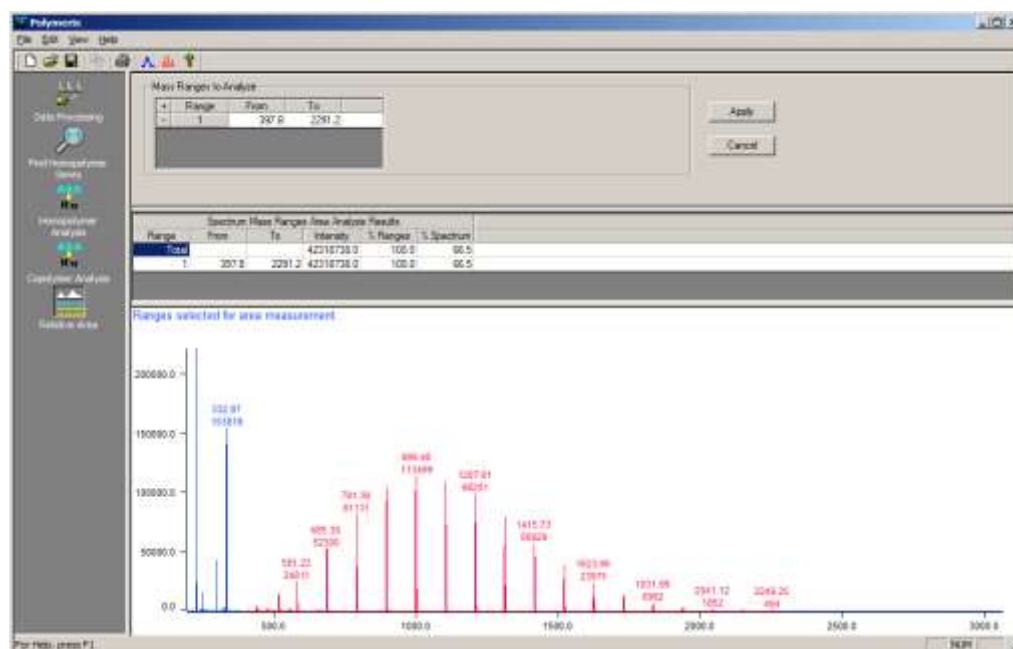

The top panel allows creation of areas by m/z range. The middle panel displays a tabulation of the total relative abundance in each range, and the bottom panel shows a graphical view of each range.

What was covered in this chapter

In this chapter, we

- Described the analytical workflow modeled by Polymerix
- Explained how that workflow is mapped to the user interface, and
- Introduced the major components and views in the user interface.

In the next chapter

We will discuss the first part of the workflow: data import, peak detection, and internal recalibration.

Data Import, Spectrum Selection, and Processing

How to import and reduce raw input spectra to a spectrum appropriate for polymer series assignment

The most critical initial steps in assigning polymer compositions to mass spectral components are selecting spectra to be co-added from a multi-scan data set, accurately processing the input spectrum to create an optional centroid peak list, correcting the mass axis, deisotoping, and applying corrections to accommodate loss of resolution at high m/z .

Goals for this chapter

In this chapter, we will explain how to:

- Use the data browser to select and import raw data from vendor format data files.
- Select a range of spectra to be co-added from a multi-scan data set.
- Select an m/z sub-range to exclude non-polymer peaks
- Set peak detection parameters.
- Apply a mass axis adjustment to reduce mass calibration errors
- Adjust spectral assignment to accommodate resolution loss at high m/z .
- Use deisotoping to merge higher isotopomer peaks into a single monoisotopic peak in each isotope cluster.

Importing data

When you installed the Polymerix software, you selected one or more data format options. This then installed specialized data import software that recognizes files in those vendors' formats and a browser that can display information contained in those files when navigating through your hard disk folder hierarchy. In addition, Polymerix supports import of peak list data contained in ASCII mass, intensity peak list files.




The data import browser

1. Start the software (as usual, by double-clicking the Polymerix icon). Click the **Data Processing** button in the Workspace Control panel.
2. To import vendor format MS data, from the **File** menu click on the **Import Chromatogram** command.




Alternatively, click the **Chromatogram** button in the **Import** section of the **Spectrum Select** tab pane from the Data Processing page.

DATA IMPORT AND PROCESSING

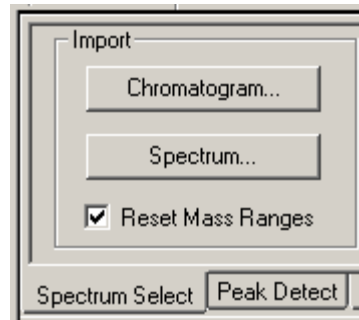

3. The data import browser appears. In the figure below, there are two MS data sets in Waters MassLynx data format shown. Both have been recognized by the browser as raw MS data formats, so they are displayed with a special spectrum icon. If you selected additional formats when you installed Polymerix and those datasets were also in this folder, they would also be shown with spectrum icons.

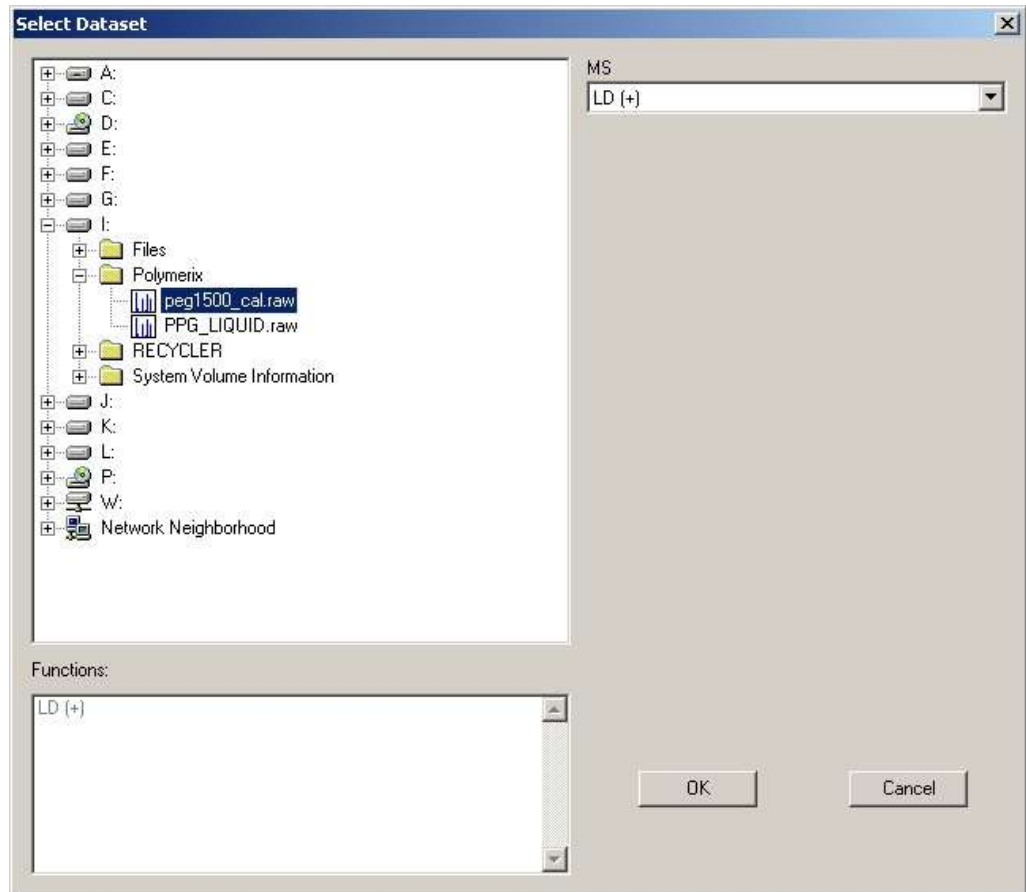


4. Selecting one of the datasets will display a list of the scan functions it holds (using the vendor's nomenclature). In this case, the MassLynx file contains a single scan function in positive ion mode, produced with MALDI (LD) ionization. Click the **OK** button to import the selected

DATA IMPORT AND PROCESSING

data. Click the **Cancel** button to close the browser without importing data.


5. Some datasets may contain multiple spectra, for example where a sample was infused over time or multiple MALDI laser shots were acquired. These are automatically recognized by the data import browser. After selecting the dataset and clicking OK, the total ion chromatogram (TIC) or its equivalent for infusion or other multi-scan acquisitions will be displayed in the TIC pane of the Data Processing page.

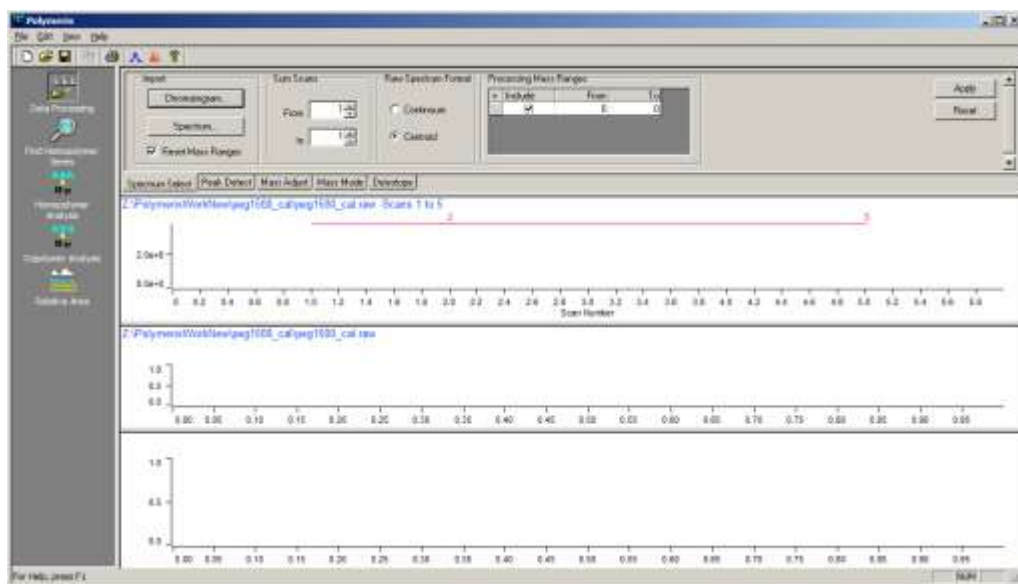

6. To select a single spectrum, simply right-click once in the Total Ion Chromatogram pane at the appropriate scan time or number. To select and co-add a range of spectra, right-click and, while still holding the mouse button down, drag across the desired time range in either direction, then release the mouse button. In each case, the selected scan or scan range will be extracted from the dataset, co-added, and displayed in the middle and bottom panes.
7. To zoom in on a region of the TIC or spectrum, left-click, drag, and release with the mouse. To zoom out, simply left-click once.
8. Each time you right-click and drag to select a spectrum range from the TIC, the input (middle pane) and target (lower pane) spectra will be replaced with the new selection. When you are satisfied with the chosen spectrum, it is time to move onto peak detection and processing.
9. To import a single spectrum from an ASCII peak list file, from the **File** menu click on the **Import Spectrum** command.



10. A standard file browser will be displayed. Select the desired file, then click **OK**. If the file contains valid spectral data, it will be imported and displayed in the middle and bottom spectrum panes. The TIC pane will be empty.
11. By default, Polymerix assumes that data read from peak list files is in centroid mode. If this is not the case, change to profile mode by clicking the **Continuum** button in the **Raw Spectrum Format** section of the **Spectrum Select** page.

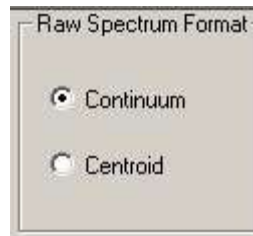


Alternatively, you may edit the ASCII file to add the word *continuum* as the first line in the file to permit the file reader to automatically identify the mode.


12. ASCII peak list files must have the following format: Each line contains a single m/z , intensity pair, separated by a comma, space, or tab delimiter. Most mass spectral data systems will automatically create this format when exporting a peak list.

Restricting the processing mass range


Occasionally a spectrum will contain contaminants or other non-sample peaks such as MALDI matrix ions or chemical noise. If these are in a region of the spectrum that can be isolated from the region containing sample peaks, then you may use the **Processing Mass Ranges** parameters to define one or more spectral ranges to be included for peak assignment.

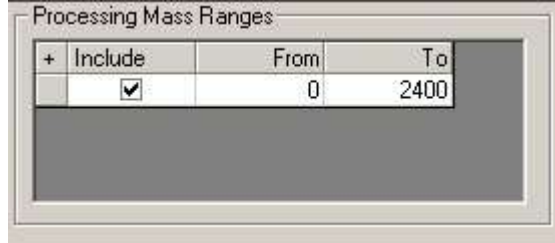


- On the **Spectrum Select** tab, there is a table containing a list of mass ranges that will be included for processing. By default, then entire spectral range is included.

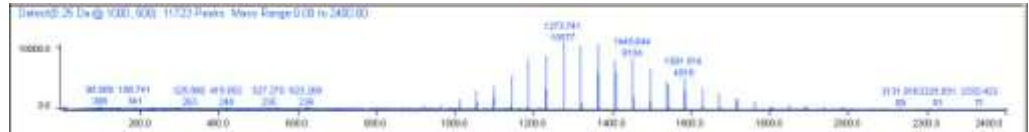

- You chose a restricted mass range in one of two ways: First, by directly editing the mass values in the table, or second, by a right-click and drag operation in the spectrum window. If you edit the mass values manually, click the **Apply** button after setting them. If you right-click and drag, the table and spectrum will be updated automatically. In the spectrum above, there are no useful peaks outside of the 800 – 2000 Da range, so a right-click and drag was used to restrict the processing range, as shown below.

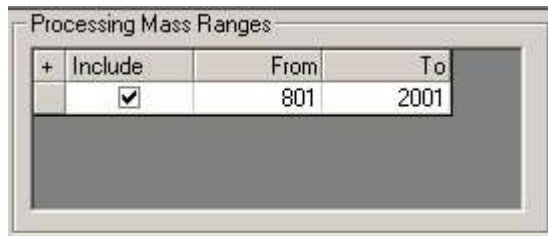


- The processed target spectrum in the bottom pane is also updated to eliminate peaks outside of the chosen range:

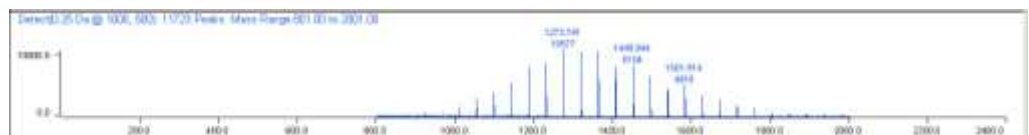

DATA IMPORT AND PROCESSING

11. You may select additional ranges if needed, for example, if a major impurity peak occurs in the midst of sample peaks, and you wish to exclude this peak from processing. In this case, click the **+** button at the upper left of the **Processing Mass Ranges** table to add another row. Select the new sub-range by editing or a right-click and drag in the spectrum. Both the table and processed spectrum will be updated with the new range. In this case, an arbitrary new range was selected for illustration, even though it obvious contains only noise.


+ Include	From	To
<input checked="" type="checkbox"/>	801	2001
<input checked="" type="checkbox"/>	592	759

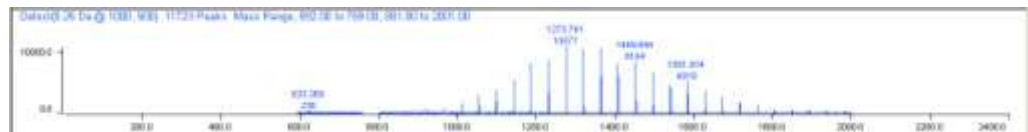

12. To remove a sub-range, click the **-** button at the left edge of the row. You may also temporarily remove a mass range by unchecking the checkbox on that row.
13. When all mass ranges are specified, click the **Apply** button to set them in the parameters list.

Detecting peaks

After importing data and selecting a spectrum range, the Data Processing page should look similar to the figure below.

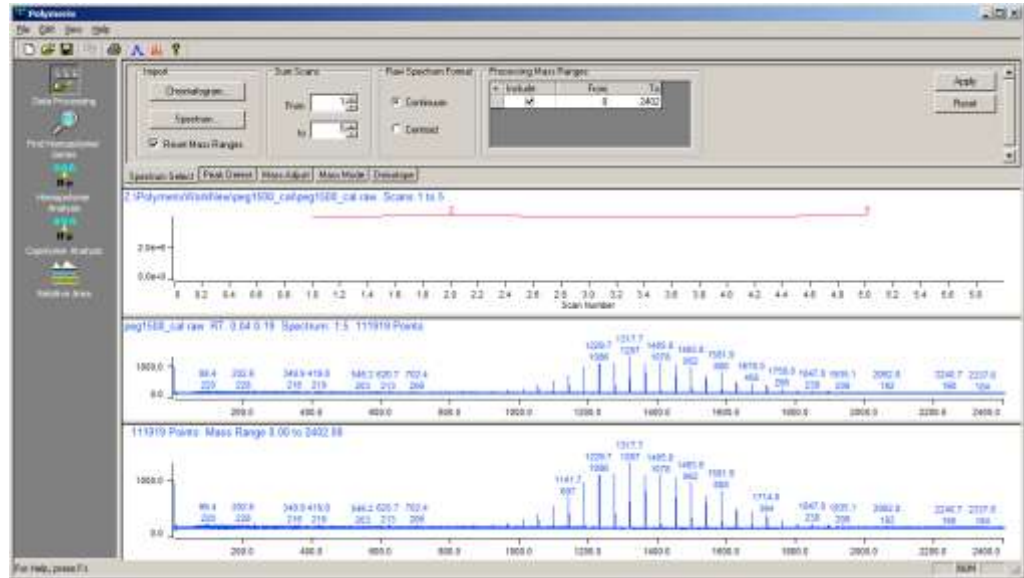

Evaluating peak detection parameters

For most modern instruments, the resolution is high enough to resolve individual isotopic peaks in isotope clusters over most if not all of the m/z range. In this case, it is generally desirable to process data in peak detected (or centroid) mode. For data where peaks are not fully resolved, it is possible to process without centroiding (i.e. in profile mode) with often very satisfactory results.


Polymerix's peak detection parameters allow for three different instrument scan function types:

- Constant (or near constant) width peaks, such as produced by quadrupole or high-resolution FT-ICR or orbitrap instruments;
- Proportional width peaks, such as produced by magnetic sector instruments; and
- TOF-mode peak widths, as produced by time of flight instruments.

Select the mode appropriate for your data type, then click the **Apply** button. Evaluate the results of peak detection by zooming in on a small section of the bottom (target) spectrum, and compare the middle (profile) spectrum with the bottom (peak detected) target spectrum.

DATA IMPORT AND PROCESSING

To zoom out one step, simply click the left mouse button without dragging.

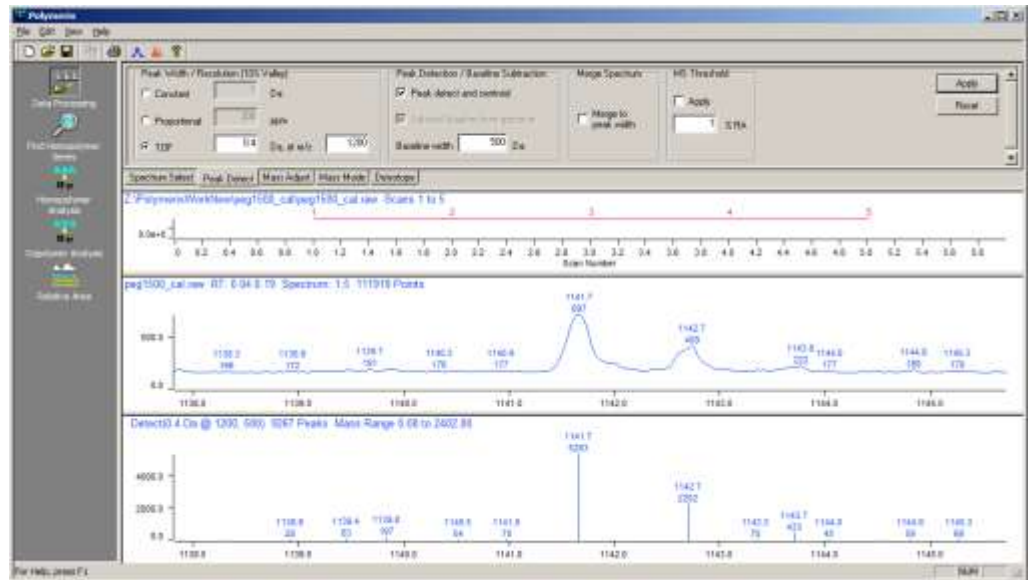


Are there detected peaks in the bottom spectrum where you would expect them to be, based on the profile peaks in the upper spectrum? In this case, the peak detection parameters are probably OK and do not need to be changed.

Are closely spaced profile peaks merged into a single centroid peak? Are profile peaks split into two centroid peaks? Either of these situations indicates that the peak detection parameters are not set correctly.

Setting peak detection parameters

To set peak detection parameters, click the **Peak Detect** tab on the top pane of the **Data Processing** page.

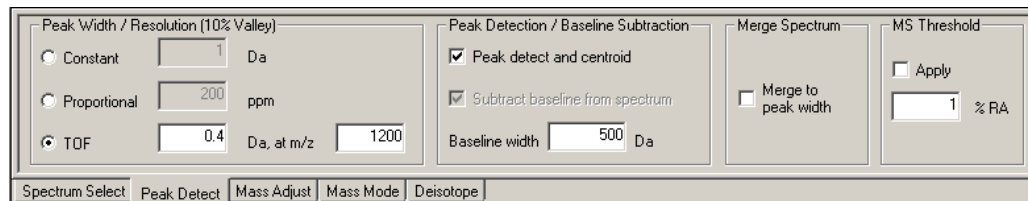

This displays the **Peak Detect** view, shown above.

The scan law determines how the peak width changes as a function of m/z . The peak width value is measured using full width at 5% peak height, or 10% valley. Depending on your instrument type, click the button to the left in the **Peak Width / Resolution** section to select the appropriate type. TOF spectra require two parameters: the peak width (in Da) and the m/z value at which the width was measured. The peak width does not need to be precise – within a factor of two is generally acceptable.

DATA IMPORT AND PROCESSING

1. For very high resolution spectra, such as produced by FT-ICR or Orbitrap instruments, Polymerix's peak detection algorithm works well with a very small peak width, such as a constant width of 0.001 Da. Polymerix, however, will not always detect peaks with a width of less than 4 data points.
2. After choosing the peak width type and changing the width parameter, also select the **Peak detect and centroid** check box in the **Peak Detection / Baseline Correction** section, and then click the **Apply** button. This will process and redisplay the new processed target spectrum in the bottom pane.
3. If you observe peak splitting in the peak detected spectrum (profile peaks are split into two or more centroid peaks), this indicates that the peak width parameter is too small. Missed peaks (or in TOF mode, too many noise peaks or baseline anomalies) indicate a peak width that is too large. Change the value, click **Apply**, and evaluate the result. Repeat if necessary until the centroid spectrum is satisfactory.
4. Polymerix's peak detection algorithm always assigns intensities to the centroided peaks based upon mathematical areas calculated from the spectrum being peak detected.
5. There are three additional sets of parameters on this tab: Baseline correction, peak merging, and peak thresholding.
6. The **Subtract baseline from spectrum** option is used to flatten a sloping baseline to eliminate systematic errors in peak intensity due to a non-zero baseline. It is automatically selected when **Peak detect and centroid** is chosen. If you choose to process in profile mode (i.e. no peak detection), the **Subtract baseline** option is available separately. Using this option is generally recommended. The **Baseline width** parameter determines the width of the spectrum used when averaging to compute the baseline offset at each point. The default value of 500 Da is usually adequate to compensate for broad baseline drifts. If the baseline has more narrow artifacts, using a smaller parameter might provide a better fit.
7. The **Merge to peak width** option helps when the spectra are very noisy, and peak detection results in a cluster of smaller peaks around the true centroid. Selecting this option will merge together each of these peak clusters to yield a single peak using an intensity-weighting algorithm. The peak width mode and peak width parameter selected for peak detection will be used to set the width of the merge window.
8. The **MS Threshold** option is used to filter the smallest peaks out of a peak detected spectrum based upon a relative abundance threshold value.


- Once all parameters have been chosen, click the **Apply** button to set them.

Mass adjustment


Overview of mass adjustment

Occasionally, the m/z assignment read from the data file is incorrect due to inadequate calibration in the instrument software. Polymerix offers a simple way to adjust the mass axis by applying a zero order mass offset, linear slope, or both.


If you suspect a problem in mass assignment, click on the **Mass Adjust** tab of the **Data Processing** parameters pane.

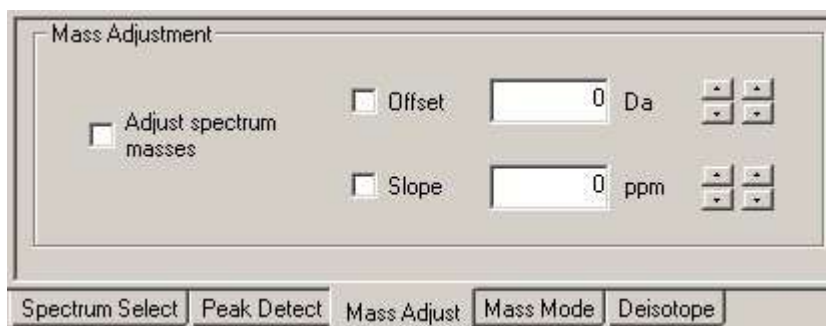

- Click the **Adjust spectrum masses** checkbox to enable mass adjustment mode.
- To apply a mass offset, click the **Offset** checkbox to enable offset adjustment. You may either type an mass adjustment value (in Da) into the edit box, or use the two sets of arrow buttons to interactively change it. The left-hand set of arrow buttons adjusts by 1 Da, while the right-hand set adjusts by 0.1 Da. As the buttons are clicked, the spectrum will shift left or right accordingly, as shown in the pictures below.

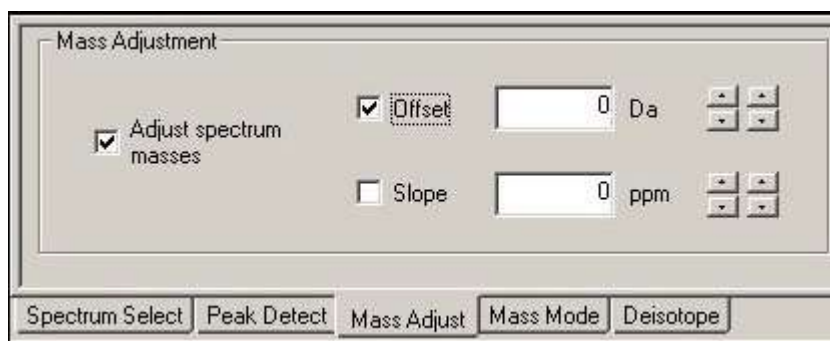

DATA IMPORT AND PROCESSING

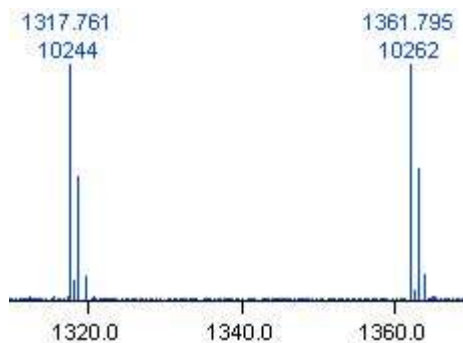


Mass Adjustment

Adjust spectrum masses



Offset Da

Slope ppm

Spectrum Select Peak Detect Mass Adjust Mass Mode Deisotope

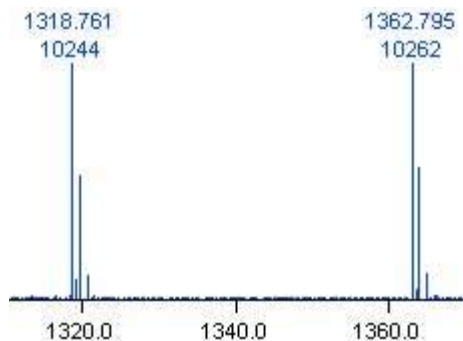

16. To apply a linear slope adjustment, click the **Slope** checkbox. As with offset adjustment, the left-hand arrow buttons adjust the slope in coarse increments of 10 ppm, while the right-hand set adjusts in increments of 1 ppm. Note that in the plot above, the two monoisotopic peaks differ by 44.034 Da, while in the plot below the difference is 44.065 Da. The 10 ppm positive slope results in an increasing mass difference with m/z, whereas an offset alone simply shifts the spectrum while maintaining the same mass differences.

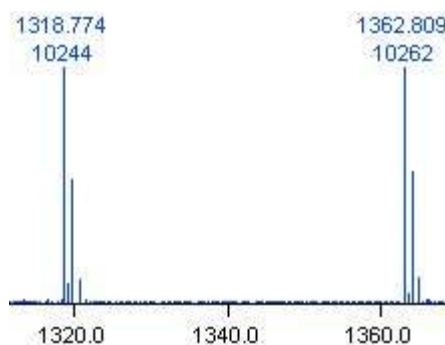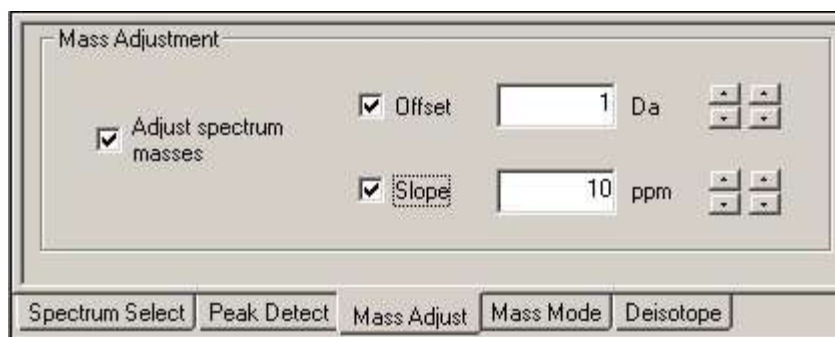

17. Only the processed target spectrum (bottom pane) is changed. This is the convention used throughout the **Data Processing** page – the middle pane shows the raw input spectrum, the bottom panes shows the processed target spectrum that will be used for polymer series assignment.
18. When you are satisfied with the mass adjustment, click the **Apply** button to set the parameters.
19. Often it is necessary to use the mass adjustment in concert with series assignment in an iterative process to ensure that all peaks are assigned. For this purpose, the **Mass Adjust** tab is duplicated on the **Homopolymer Analysis** and **Copolymer Analysis** pages.

Mass mode


In Polymerix, it is most common to do analysis on peak-detected data, with assignments to monoisotopic peaks. For users of instruments with insufficient resolution to resolve isotopic peaks throughout the spectrum, analysis is a bit more difficult when peak detection cannot be uniformly applied. For this purpose, Polymerix permits processing in monoisotopic, chemical (average) or nominal mass modes, and will apply a correction to the matched peaks to convert them to either mode as necessary.

For example, if resolution is degraded above 2000 Da, the detected peak m/z values are the chemical m/z values of the whole cluster, not of the individual isotopic peaks. In this case, you may choose to have molecular weight statistics reported as

DATA IMPORT AND PROCESSING

monoisotopic or chemical mass, and Polymerix will smoothly adjust as the m/z region changes.

The default mass mode is monoisotopic throughout the m/z range. To change the mass mode, click on the **Mass Mode** tab in the **Data Processing** parameters pane.

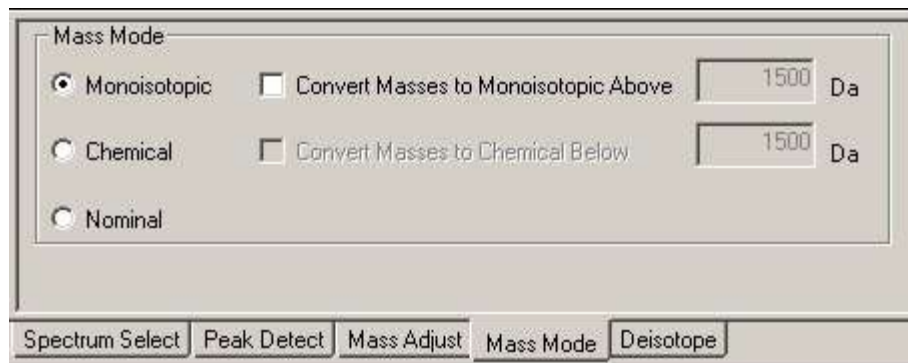

1. As can be seen above, you may select one of three processing modes. If **Monoisotopic** is selected, all results will be reported based on assignments to monoisotopic peaks. Likewise, if **Chemical** or **Nominal** is selected, results will be reported in that mode.
2. For monoisotopic mode, an additional option is to convert unresolved chemical mass peaks to the equivalent monoisotopic peak m/z above the given m/z value. If this option is selected, polymer formulas will be matched to the spectrum based on chemical mass, and the result converted to the equivalent monoisotopic mass above the threshold.
3. For chemical mass mode, the additional option allows matches to monoisotopic peaks to be converted to the equivalent chemical m/z below the given m/z value. If this option is selected, formulas will be matched to monoisotopic peaks based on monoisotopic mass, then converted to the equivalent chemical mass below the threshold.
4. In either of the monoisotopic or chemical mode cases, the optional processing ensures that mass errors due to lack of resolution are minimized, and that mass values are consistently reported across the entire spectral range.

Deisotoping

The final topic in this data processing chapter is deisotoping. This is the process of applying a model chemical formula to each isotope cluster in a centroided spectrum, and merging all higher isotopomeric peaks into a single monoisotopic peak.

The new centroid cluster matching feature which was added to Polymerix in version 3.0 should be used instead of deisotoping. Tools on the Find Homopolymer Series page will only work with non-deisotoped spectra.


In monoisotopic mass mode, Polymerix will match predicted polymer m/z values only to monoisotopic peaks and use the resulting matched intensity in the molecular weight statistics calculations. For polymers with a wide polydispersity (and thus wide mass range), this will result in bias since higher m/z isotope cluster have more abundance in higher isotopes. To eliminate this bias, the deisotoping option is provided.

Note: Deisotoping is only available when the processed spectrum is peak detected. If peak detection is not selected, the deisotoping parameters will be disabled.


To set deisotoping parameters, click the **Deisotope** tab on the **Data Processing** parameters pane.

The screenshot shows a software interface for 'Isotope Correction'. It features a 'Deisotope spectrum' checkbox which is checked. Below it, 'Isotopic Mass Tolerance' is set to '100 ppm'. To the right, there are three input fields: 'Repeat Formula', 'End Group', and 'Adduct', each with a button next to it. At the bottom, a series of tabs includes 'Spectrum Select', 'Peak Detect', 'Mass Adjust', 'Mass Mode', and 'Deisotope', with 'Deisotope' being the active tab.

1. The deisotoping algorithm operates by predicting an isotope cluster from a model chemical formula, matching that to an experimental isotope cluster, then moving the intensity found in the isotopic peaks back into the monoisotopic peak. The model chemical formula is created by adding together an end group composition, adduct, and an appropriate multiple of the repeat formula to yield a model formula with mass close to the experimental mass. The predicted peaks and their relative abundances are then matched to experimental data. This process is repeated for each isotope cluster in the spectrum. To enable deisotoping, click the **Deisotope spectrum** checkbox.
2. Next, you must specify at least the repeat formula. The end group and adduct entries are optional. To specify a repeat formula, click the button next to the **Repeat Formula** edit box.

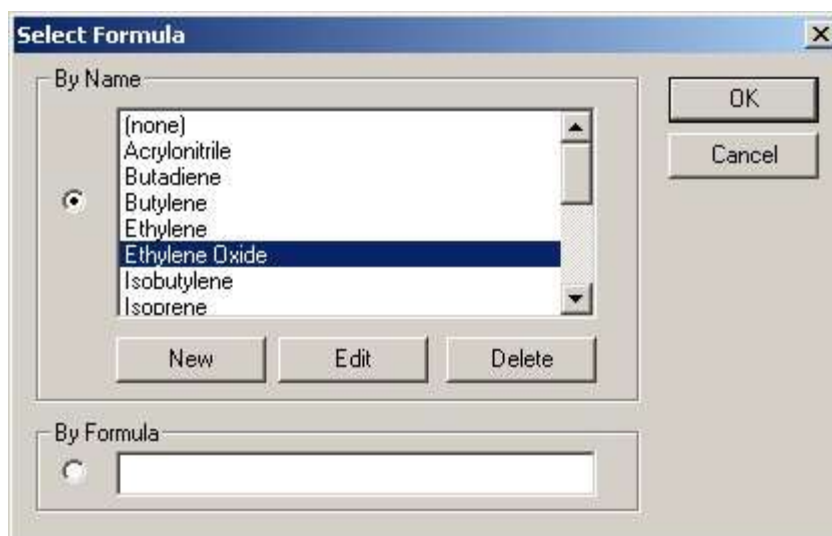


3. This displays the **Select Formula** dialog box. In the top part of the dialog is a list of predefined monomers (repeat units). If your sample includes one of these, you may select it and click the **OK** button.
4. If your monomer is not on the list, you may add a new definition which will be stored for use again in the future. In this case, click the **New** button, and the **Edit Chemical Entity** dialog will be displayed:

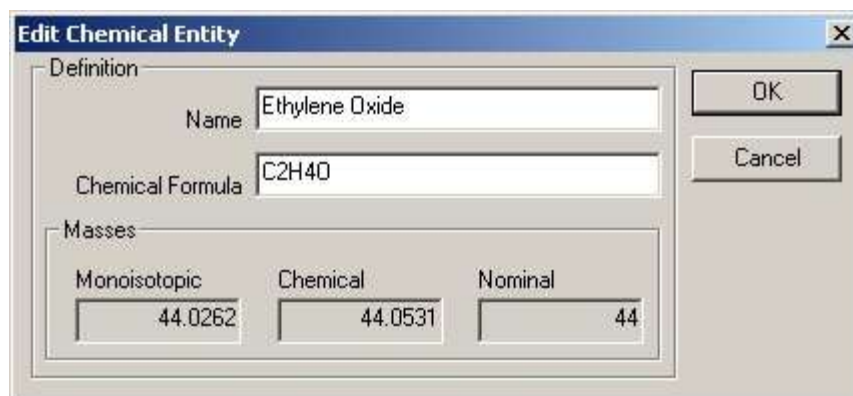

5. Enter a name for the monomer and supply the chemical formula (in proper chemical nomenclature). The formula will be validated, and the masses automatically calculated. Click the **OK** button close the dialog box. The new entry will be added to the list in the **Select Formula** dialog.
6. To specify a repeat formula for a one-time use, type the formula in the **By Formula** edit box. These formula are temporary and are not stored for re-use. Typing into this box will automatically select this option. Click the **OK** button to validate the formula and close the **Select Formula** dialog box.

- After closing the dialog box, the selected entry or formula will be added to the parameters.

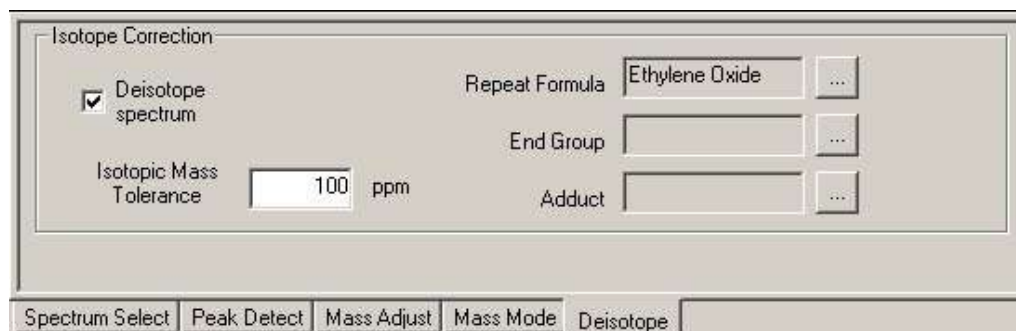

- You may optionally select an end group. This may be important if the end group has a significant contribution to the observed isotope cluster, for example, if it contains chlorine, bromine, silicon, or sulfur. Click the button to the right of the **End Group** edit box to display the **Select End Group** dialog box. As with the repeat formulas, there is a predefined list. You may choose one of these if appropriate, or define a new one using the same process as for repeat formulas. Click the **OK** button to close the dialog and add the formula to the end group parameters.




- Finally, if there is an adduct with an unusual isotopic pattern, such as lithium or silver, you may wish to specify that also. Click the button to the right of the **Adduct** edit box to display the **Select Adduct** dialog box.. As with the previous two, there is a predefined list or you may define a new one. After making your choice, click the **OK** button to close the dialog box and put the adduct in the parameter set.

DATA IMPORT AND PROCESSING



10. Finally, to perform the deisotoping, click the **Apply** button. The effect of deisotoping is shown in the example below. You can see in the zoomed view that the higher isotopes have been completely removed and merged into the monoisotopic peak.

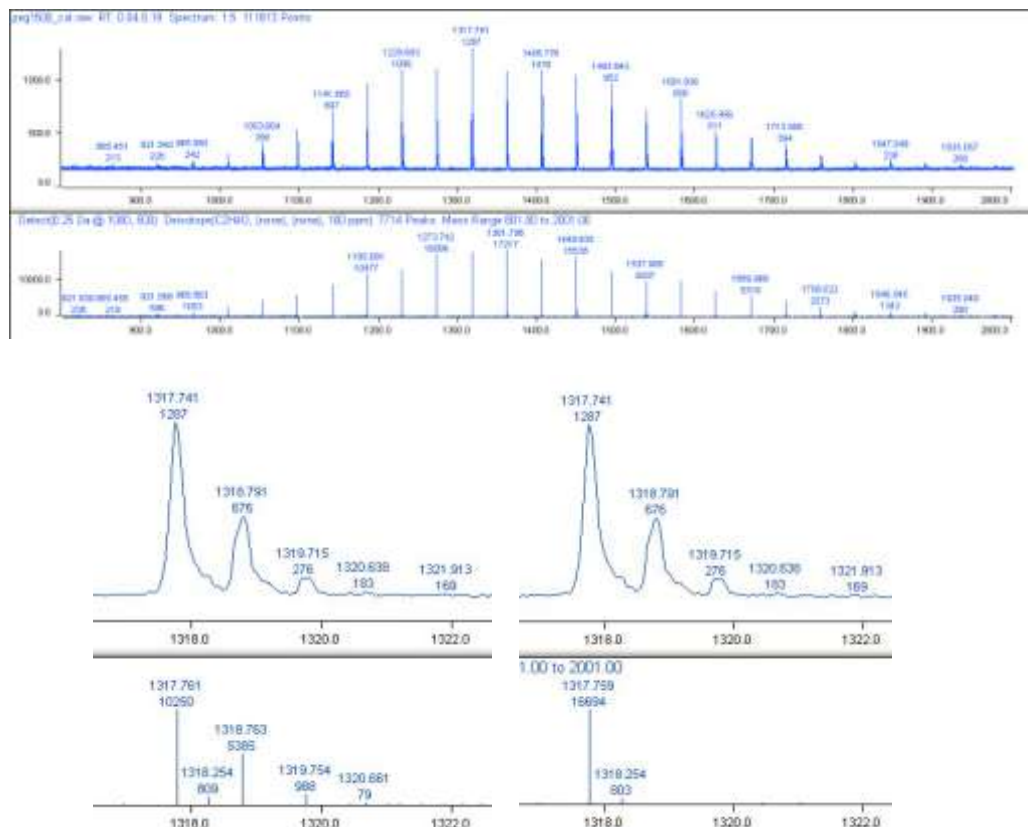


11. It is important to note that the deisotoping algorithm never removes more of an isotopic peak than the model predicts. If there are overlapping isotope clusters where a higher isotopomer of one cluster is isobaric with the monoisotopic peak of another, only as much of the higher isotopomeric peak is removed as is predicted from the formula, leaving a residual peak. This could happen for a polymer


mixture with two different end groups, one containing a C=C double bond and the other a C-C saturated bond. This is a 2 Da difference, and could result in an isobaric overlap between the monoisotopic peak of the saturated species with the M+2 isotope of the unsaturated one.

Saving and loading parameter template files

After all that work of creating appropriate processing parameters, it would be a shame if you had to re-enter all that information every time you used Polymerix on a similar sample. To avoid this, Polymerix allows storing parameters in files as named templates. Polymerix parameter template files use the **.pmt** file extension.

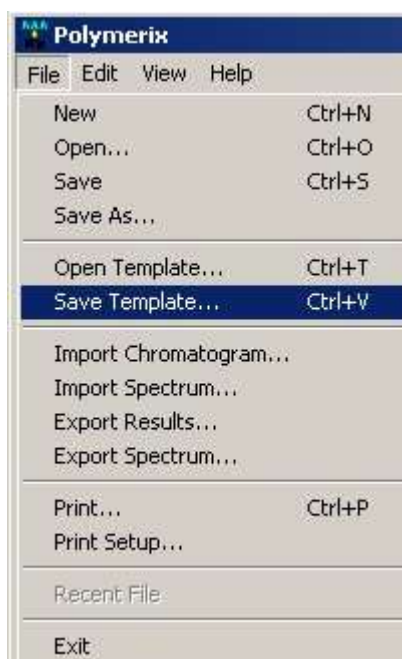

1. To save parameters to a template file, use the **File** → **Save Template...** menu command. This allows you to choose a file name and save the parameters.

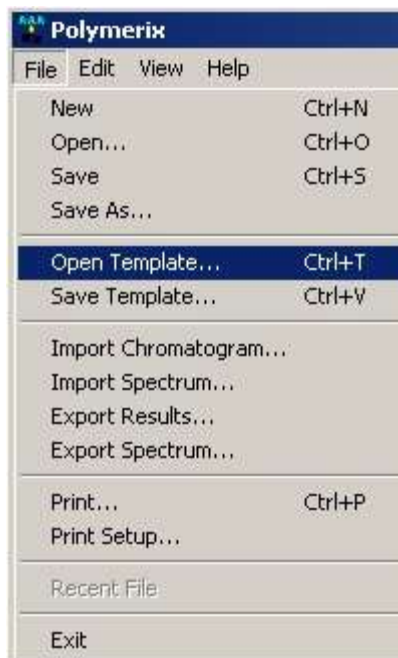

2. To load a previously saved template, use the **File → Open Template...** menu command. This will display a file selection dialog, from which you can choose the desired template file. When you close the file dialog by clicking its **OK** button, the template file will be read in and all of the current parameters replaced with the new ones read from the file.

Parameter template files do not contain any experiment information, only parameters. Therefore parameter templates can be applied to experimental data of any sample of a similar type. They can be transferred from PC to PC, since they also do not contain any absolute folder or file path information.

What was covered in this chapter

In this chapter,

- We presented the Data Import browser and described how to select and import a data file. We also described how, in the case of a multiple-scan dataset, you can select a single spectrum or a range of spectra to be co-added prior to import.
- We described how to select spectral sub-ranges for processing.
- We discussed the peak detection process and how to set peak detection parameters.

DATA IMPORT AND PROCESSING

- We presented ways to overcome inadequate resolution and mass assignment by changing the mass mode and by adjusting the mass axis.
- We discussed deisotoping and how this important algorithm can overcome bias in computing molecular weight distribution statistics.
- And finally, we described how to save and load parameters using Polymerix template files.

In the next chapter

In the next chapter we will describe the Find Homopolymer Series page and how it can be used to assist in determining and specifying the composition of an unknown polymer.

Find Homopolymer Series

How to use the find homopolymer series page to analyze unknown or partially known polymers

Polymerix is normally used to deconvolve polymer series and compute molecular weight distributions for materials of known chemical composition. The **Find Homopolymer Series** page can help a user determine and specify the chemical composition of polymers in a given spectrum.

Goals for this chapter

In this chapter, we will:

- Provide a brief high level overview of the polymer series finding process and its requirements.
- Show how the tools are used with a very simple example.
- Introduce the concept of reference masses and describe how they are used.
- Show details about how to use the tools on the **Find Homopolymer Series** page to construct repeat unit formula, end group formula, and adduct formula specifications.
- Provide tips for solving common problems such as not being able to find any results.

Overview

The spectrum being analyzed must have been peak detected and should not have been deisotoped. The masses of selected monoisotopic peaks from the spectrum are analyzed and candidate lists of the potential chemical components which correspond to those masses are generated. The user chooses from these lists of candidate compositions, generating polymer series definitions which can be used for quantitative analysis. When multiple components are present the process can be repeated.

FIND HOMOPOLYMER SERIES

Spectral data with high mass accuracy will work best with the **Find Homopolymer Series** page. Correct selection of neighboring monoisotopic peaks from the same polymer series is crucial for the correct determination of chemical compositions. Finally, the tools are constrained to determining composition of homopolymer series only, as opposed to copolymer series.

Simplest use case example

The simplest example of using the **Find Homopolymer Series** tool to determine the chemical composition of a homopolymer is as follows:

- Load a dataset and select a peak detected polymer spectrum for analysis.
- Navigate to the **Find Homopolymer Series** page.
- Click the **Find** button.
- Choose a chemical formula for repeat unit from the list provided by Polymerix and click the **OK** button.

#	Repeat Formula	DBE	Charge	m/z	Error mDa	Error mDa	Error ppm
1	C8H8	4.0	1	104.06260	0.035	0.035	0.338
2	C4H8O3	0.0	1	104.04734	15.291	15.291	146.944
3	C7H4O	5.0	1	104.02621	36.421	36.421	349.988
4	C3H4O4	1.0	1	104.01096	51.677	51.677	496.594

Monoisotopic m/z values are 1101.54613 and 1205.60877, repeat m/z is 104.06264, tolerance is 0.1 Da, electron mode is normal, restricted element ratios, DBE range is from -0.5 to 5.0, charge count is 1, maximum result count is 10, C count is from 0 to 10, H count is from 0 to 20, O count is from 0 to 10, N count is from 0 to 3.

Settings...

OK Cancel

- Choose chemical formulas for combined end group and adduct from the list provided by Polymerix and click the **OK** button.

FIND HOMOPOLYMER SERIES

Select end group and adduct for series with monoisotopic peak at $m/z = 1101.54613$

#	Combined End Group	DBE	Adduct	Loss	Computed Series m/z	Error mDa	Error mDa	Error ppm
1	C4H10	0.0	Ag		1101.54620	0.064	-0.064	-0.058
2	C5H20S	5.0	Na		1101.54285	3.287	3.287	2.984
3	C7H18N2O2	0.0	Ag		1101.54217	3.959	3.959	3.594
4	C6H18N4O	0.0	Ag		1101.55341	7.274	-7.274	-6.604
5	H2N2O7	1.0	Na		1101.53882	7.309	7.309	6.636
6	C4H2N2O4	5.0	Na		1101.55408	7.947	-7.947	-7.214
7	C8H18O3	0.0	Ag		1101.53094	15.192	15.192	13.792
8	N6O5	4.0	H		1101.56370	17.563	-17.563	-15.944
9	C4H14N6O	1.0	Ag		1101.52826	17.878	17.878	16.230
10	C5H18N6	0.0	Ag		1101.56464	18.508	-18.508	-16.802

Additional Repeats:

Result End Group:

Monoisotopic m/z is 1101.54613, repeat formula is C8H8, repeat mass is 104.06260, charge count is 1, tolerance is 0.1 Da, electron mode is normal, restricted element ratios, DBE range is from -0.5 to 5.0, minimum end group m/z is 0, maximum end group m/z is 200, maximum result count is 10, C count is from 0 to 20, H count is from 0 to 40, O count is from 0 to 20, N count is from 0 to 6.

Settings...

OK Cancel

In this case Polymerix creates a new polymer series definition using C8H8 (styrene) for its repeat formula, a combined end group of C4H10, and Ag (silver) as the adduct. Peaks from the spectrum being analyzed which match the series definition are extracted and displayed.

Understanding reference peaks

The **Find Homopolymer Series** tool uses pairs of adjacent monoisotopic peaks in a polymer mass spectrum to determine potential repeat compositions and combinations of end group and adduct. Pairs of reference masses are assumed to differ only by the monoisotopic mass of the repeat unit, corresponding to molecular ions whose formulas differ in composition only by a single repeat unit.

Automatic versus manual selection

Monoisotopic reference masses can be selected either automatically or manually. Automatic reference mass selection is accomplished using the **Find** button. Manual reference mass selection is accomplished by right clicking and dragging on spectrum plots.

Excess Mass and errors

Whenever Polymerix has both a reference mass and a hypothetical polymer formula the amount that the reference mass is in excess of the m/z of its closest corresponding molecular ion in the polymer series is calculated and displayed either as “Excess Mass” in mDa (milliDaltons) or as “Error” in both mDa and ppm. Result lists of chemical

FIND HOMOPOLYMER SERIES

compositions corresponding to a reference peak are sorted in order of increasing magnitude of calculated errors.

How Polymerix handles reference peak selection

When the user selects a pair of reference peaks either automatically with the **Find** button or manually by right clicking and dragging on a spectrum plot Polymerix responds as follows:

- The reference peaks are highlighted in red where they are shown in spectrum plots.
- The mass difference between the peaks is computed for use in analyzing the repeat unit's chemical composition. If no repeat unit is currently specified or if the mass difference is incompatible with the currently specified repeat unit then the user is presented with a list of repeat unit chemical formulas to choose from.
- The user is presented with a list of end group and adduct formulas to choose from. Polymerix creates a new polymer series definition from the user's choices.
- Reference peak highlighting is removed where it is shown in spectrum plots.


Automatic reference peak selection

Polymerix uses a heuristic algorithm when attempting to automatically select reference peaks from a spectrum. The algorithm attempts to identify large monoisotopic peaks corresponding to a homopolymer series. It takes account of the range of possibilities for adduct chemical composition when attempting to determine which peak within an isotope cluster is the monoisotopic peak. This accounting is important for higher mass polymers when silver (Ag) is used as an adduct.

Automatic reference peak finding ignores peaks in the spectrum being analyzed which match any existing polymer series specifications.

Because the automatic reference peak finding algorithm is based on a heuristic, it may not always correctly identify monoisotopic reference peaks from the same series, and it may select reference peaks which seem less than ideal. If the wrong peaks are identified the repeat unit selection dialog and the end group selection dialog will often show no potential chemical compositions or only chemical compositions which don't make sense to the user. In such cases, manual selection by right clicking and dragging on spectrum plots may provide a workable alternative.

Find Homopolymer Series page

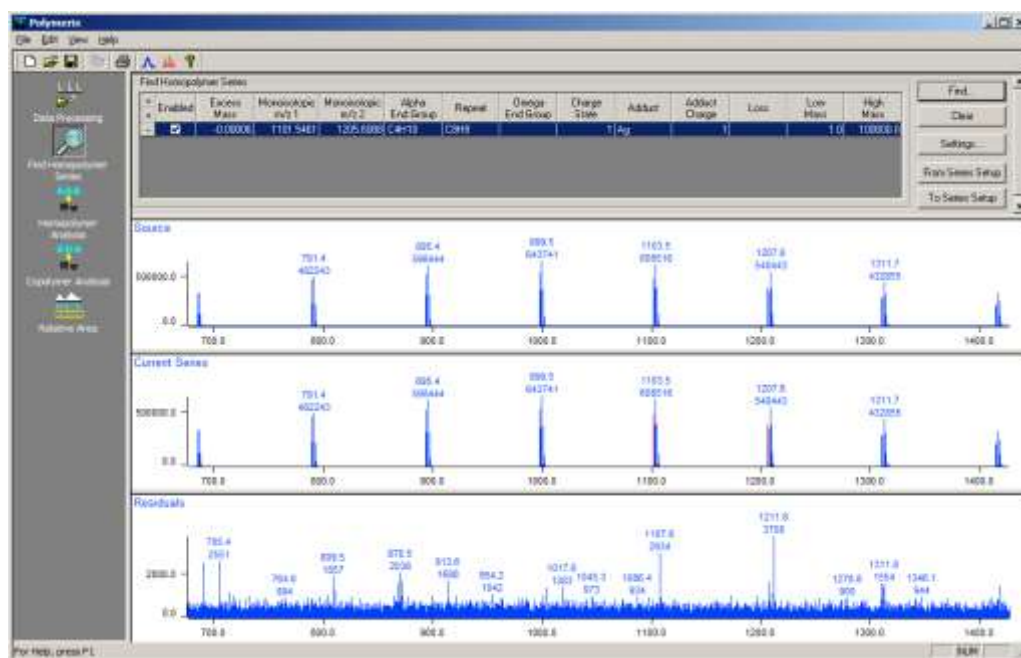

In the topmost panel series specifications and reference masses are displayed. The top spectrum plot shows the complete spectrum being analyzed. The middle spectrum plot shows the currently selected series with its reference peaks highlighted in red. The bottom plot displays the residual spectrum – in other words, all peaks still unassigned. In the above example the residual spectrum mass peaks are all less than 2% RA.

Series list

The Series list shows a list of polymer series specifications which includes a mass range for each series and chemical formulae for repeat unit, end groups, and adduct. When a series is selected as the current series, its row is highlighted and the current series spectrum plot will show peaks which match the series.

The Series list also displays the two reference masses used to determine the chemical composition of the polymer and the “excess mass” at the first reference mass, which indicates the difference between the reference mass and its closest calculated mass from the polymer series which is being specified.

Clicking on an excess mass field when either end group is empty will bring up the **End group and adduct selection** dialog. Clicking on most of the other columns in the series list will allow individual components of the polymer series definition to be entered or edited. The reference masses cannot be entered or edited.

Find button

This invokes the automatic reference peak finder which attempts to find a pair of reference peaks from the residual spectrum which correspond to a polymer series.

Clear button

This clears all polymer definitions in the series list.

FIND HOMOPOLYMER SERIES

Settings button	This brings up the Find Series Settings dialog.
From Series Setup button	This transfers series definitions from the Homopolymer Analysis page. Additionally, some settings such as spectrum match tolerance and charge polarity are transferred from Homopolymer Analysis .
To Series Setup button	This transfers series definitions to the Homopolymer Analysis page for subsequent quantitative analysis. In addition to the series definitions, the spectrum match tolerance and polarity from the Find Series Settings dialog are transferred to Homopolymer Analysis . Finally, the Homopolymer Analysis mass mode is set to “monoistopic” and the use clusters option is enabled.
Source spectrum plot	This shows the spectrum being analyzed. Right clicking and dragging on the plot allows the user to select reference peaks for elemental composition determination.
Current series spectrum plot	This shows peaks from the spectrum being analyzed which match the currently selected series from the series list. Peaks used as reference masses in the series are colored in red. Right clicking and dragging on the plot allows the user to select a mass range for the series.
Residual spectrum plot	This shows the peaks which remain after peaks which match enabled series specifications in the series list have been removed. Right clicking and dragging on the plot allows the user to select reference peaks for elemental composition determination.

Selecting the repeat unit

Repeat unit selection dialog	This dialog is displayed when the user selects a new pair of reference masses. It allows the user to select the polymer’s repeat unit as well as charge state.
-------------------------------------	--

FIND HOMOPOLYMER SERIES

Select repeat and charge for mass difference 104.06264, from m/z 1101.54613 to m/z 1205.60877

#	Repeat Formula	DBE	Charge	m/z	Error mDa	Error mDa	Error ppm
1	C8H8	4.0	1	104.06260	0.035	0.035	0.338
2	C4H8O3	0.0	1	104.04734	15.291	15.291	146.944
3	C7H4O	5.0	1	104.02621	36.421	36.421	349.988
4	C3H4O4	1.0	1	104.01096	51.677	51.677	496.594

Monoisotopic m/z values are 1101.54613 and 1205.60877, repeat m/z is 104.06264, tolerance is 0.1 Da, electron mode is normal, restricted element ratios, DBE range is from -0.5 to 5.0, charge count is 1, maximum result count is 10, C count is from 0 to 10, H count is from 0 to 20, O count is from 0 to 10, N count is from 0 to 3.

Settings...

OK Cancel

The currently selected row represents the user's selection of repeat unit composition. In the bottom panel, all parameters that were in effect are shown. These parameters can be adjusted using the **Settings** button, which brings up the **Find Series Settings** dialog.

Selecting end groups and adducts

End group and adduct selection dialog

Polymerix displays this dialog when reference peaks are selected, or when the "Excess Mass" column for an entry in the Series list is clicked and either the "Alpha End Group" or "Omega End Group" column for the entry is empty.

FIND HOMOPOLYMER SERIES

Select end group and adduct for series with monoisotopic peak at m/z = 1101.54613

#	Combined End Group	DBE	Adduct	Loss	Computed Series m/z	Error mDa	Error mDa	Error ppm
1	C4H10	0.0	Ag		1101.54620	0.064	-0.064	-0.058
2	C5H205	5.0	Na		1101.54285	3.287	3.287	2.984
3	C7H18N2O2	0.0	Ag		1101.54217	3.959	3.959	3.594
4	C6H18N4O	0.0	Ag		1101.55341	7.274	-7.274	-6.604
5	H2N2O7	1.0	Na		1101.53882	7.309	7.309	6.636
6	C4H2N2O4	5.0	Na		1101.55408	7.947	-7.947	-7.214
7	C8H18O3	0.0	Ag		1101.53094	15.192	15.192	13.792
8	N6O5	4.0	H		1101.56370	17.563	-17.563	-15.944
9	C4H14N6O	1.0	Ag		1101.52826	17.878	17.878	16.230
10	C5H18N6	0.0	Ag		1101.56464	18.508	-18.508	-16.802

Additional Repeats: None 1
 Result End Group: C4H10 C12H18

Monoisotopic m/z is 1101.54613, repeat formula is C8H8, repeat mass is 104.06260, charge count is 1, tolerance is 0.1 Da, electron mode is normal, restricted element ratios, DBE range is from -0.5 to 5.0, minimum end group m/z is 0, maximum end group m/z is 200, maximum result count is 10, C count is from 0 to 20, H count is from 0 to 40, O count is from 0 to 20, N count is from 0 to 6.

Settings... OK Cancel

This is actually two lists each of which has a highlighted selection.

In the top list, chemical compositions which differ only by an integral number of repeat units are eliminated. Additionally, the columns which are presented in the top list will vary slightly depending upon whether the end group being determined is an Alpha or Omega endgroup versus determining a combined end group.

The bottom list corresponds to the composition which is currently selected in the top list. When the End group maximum mass from Find Series Settings is large enough then the bottom list will present multiple choices of end group which differ by integral numbers of repeat units.

In the bottom panel, all parameters that were in effect are shown. These parameters can be adjusted using the **Settings** button, which brings up the **Find Series Settings** dialog.

Setting constraints on chemical composition

Find Series Settings dialog

The **Find Series Settings** dialog is used to set constraints on the chemical compositions considered when attempting to match reference peaks. The available chemical elements used for constructing repeat units, end groups, and adducts are specified here.

FIND HOMOPOLYMER SERIES

The **Find Series Settings** are made persistent on the user's local computer so that when Polymerix is next started the same settings will be in effect. None of the **Find Series Settings** are saved with Polymerix document files. The screen shot below shows the default **Find Series Settings** for Polymerix.

Find Series Settings

Repeat Tolerance: mDa ppm

End Group Tolerance: mDa ppm

Maximum Result Count:

End Group Mass: Minimum Maximum

Charge: Minimum Maximum

DBE: Minimum Maximum

Electron Mode:

Element Ratio Restrictions

Positive Charge Mixed adducts

Spectrum Match Tolerance:

Constant +/- Da

Proportional +/- ppm

TOF +/- Da, at m/z

Repeat Unit Elements Table

	Enable	Atomic Symbol	Min. Count	Max. Count
-	<input checked="" type="checkbox"/>	C	0	10
-	<input checked="" type="checkbox"/>	H	0	20
-	<input checked="" type="checkbox"/>	O	0	10
-	<input type="checkbox"/>	N	0	3

End Group Elements Table

	Enable	Atomic Symbol	Min. Count	Max. Count
-	<input checked="" type="checkbox"/>	C	0	20
-	<input checked="" type="checkbox"/>	H	0	40
-	<input checked="" type="checkbox"/>	O	0	20
-	<input checked="" type="checkbox"/>	N	0	6

Adducts

	Enable	Adduct	Loss
-	<input checked="" type="checkbox"/>	H	
-	<input checked="" type="checkbox"/>	Na	
-	<input type="checkbox"/>	K	
-	<input type="checkbox"/>		H

Element tables

When considering possible chemical formulas Polymerix restricts the formulas under consideration using element tables. There are three independent tables which correspond to repeat units, end groups, and adducts. These tables are shown on the right side of the **Find Series Settings** dialog.

For repeat units and end groups, the number of atoms of a particular element can be numerically constrained for the repeat unit or end group.

Change an entry by selecting and modifying its text. Add a new entry to an element table by clicking on the “+” in column one. Remove an entry by clicking on the “-” in column one of the entry's element table line.

FIND HOMOPOLYMER SERIES

Repeat tolerance	Repeat Tolerance is the maximum mass tolerance allowed when computing potential repeat formulas. Can be specified in mDa (milliDaltons) or ppm.
End group tolerance	End Group Tolerance is the maximum mass tolerance allowed in computing end group formulas. Can be specified in mDa (milliDaltons) or ppm.
Maximum result count	Maximum Result Count is the maximum number of repeat formulas or end group formulas that are displayed for user selection. If more than this number of potential formulas satisfy the end group or repeat tolerance then only those formulas which best match the reference mass are shown.
End group mass range	The End Group Mass Minimum and Maximum are mass limits for the mass of the end group whose composition is being determined. All end group formulas are computed using the minimum and maximum masses provided.
Charge range	Charge Minimum and Maximum define the charge count range to be used for the computations. To enable multiply charged ions Maximum Charge must be set higher than its default value of 1.
DBE range	DBE Minimum and Maximum is the allowed double bond equivalent range to be considered for both repeat formulas and end group formulas. In the case of a polystyrene polymer, a maximum of at least 4 must be used, since the styrene repeat formula (C ₈ H ₈) has a DBE value of 4.
Electron mode	This sets another restriction on DBE values, limiting whether integral and/or non-integral values are allowed. For repeat units, selecting an Electron Mode of “Normal” will limit result chemical compositions to those with integral DBE values and selecting “Reversed” will limit result compositions to those with non-integral DBE values. For end groups, composition results are limited in such a way that the sum of end group DBE values is integral when “Normal” is selected and non-integral when “Reversed” is selected. Selecting “None” disables electron mode restrictions.
Element ratio restrictions	Enabling Element Ratio Restrictions requires that the formulas for repeat and end group are “reasonable”, for example H ₁₄ won’t be allowed.
Positive charge	Specifies ion polarity. Polymerix uses this to account for electron masses when determining elemental composition. The To Series Setup button transfers this setting to homopolymer analysis and the From Series Setup button sets this setting using the corresponding value from Homopolymer Analysis.
Mixed adducts	Mixed Adducts is only used for multiply charged data. It specifies that different adducts may be combined with each other when forming multiply charged ions. For instance, a doubly charge species with H ⁺ and NH ₄ ⁺ would be allowed.
Spectrum match tolerance	The Spectrum Match Tolerance values are the same as in Homopolymer Analysis .

FIND HOMOPOLYMER SERIES

The **To Series Setup** button transfers this setting to homopolymer analysis and the **From Series Setup** button sets this setting using the corresponding value from **Homopolymer Analysis**.

Reset button

This restores all the parameters in the **Find Series Settings** dialog to default values.

Troubleshooting Tips

With high mass accuracy data and chemical composition constraints which have been properly set up, Polymerix should be able to correctly determine chemical compositions. If Polymerix is showing no results or only results which don't make chemical sense, the problem is usually that Polymerix has done a poor job of automatically selecting reference peaks for use in determining chemical compositions.

- Restricting mass ranges using the Data Processing page may help automatic reference peak selection work. Once a series is found the mass range restriction may be removed.
- Double check the results of automatic reference peak selection. Manual reference peak selection, even of a different pair of reference peaks from the same series, may solve the problem.
- Double check the chemical constraints being used. If multiply charged ion series aren't being recognized make sure the Charge Maximum setting is set higher than its default value of one.

What was covered in this chapter

In this chapter we

- Showed how to use Find Homopolymer Series page to easily set up homopolymer series definitions by selecting from candidate lists of chemical compositions generated from polymer spectral data.
- Described reference masses and how they are used by the Find Homopolymer Series page.
- Showed how to manipulate the various settings which are used when constructing candidate chemical compositions.

In the next chapter

In the next chapter, we will present the procedure for homopolymer analysis and describe the tables and graphics produced as a result of homopolymer series assignment.

Homopolymer Analysis

Defining and reviewing homopolymer series assignments

After Polymerix has applied all of the data processing parameters to produce a processed target spectrum, the time has come in the case of a homopolymer sample, to set up the definitions of the homopolymer series and assign compositions to spectral peaks..

Goals for this chapter

In this chapter, we will

- Describe the molecular weight distribution statistics calculated by Polymerix
- Illustrate in detail the user interface for creating homopolymer series definitions.
- Explain how to browse the series assignment results in tables and graphics.

Molecular weight distribution statistics

One of the primary goals of the Polymerix software is to produce molecular weight distribution statistics on the total sample as well as each of the polymer series within it. These statistics are commonly used in polymer chemistry to describe a polymer. The statistics calculated by Polymerix are:

- M_n : Number average molecular weight

$$M_n \equiv \frac{\sum_i (M_i \cdot n_i)}{\sum_i n_i}$$

- M_w : Weight average molecular weight

$$M_w \equiv \frac{\sum_i (M_i^2 \cdot n_i)}{\sum_i (M_i \cdot n_i)}$$

Molecular weight averages

HOMOPOLYMER ANALYSIS

- M_z : Third order molecular weight

$$M_z \equiv \sum_i (M_i^3 \cdot n_i) / \sum_i (M_i^2 \cdot n_i)$$

Polydispersity

- **PD**: Polydispersity

$$PD \equiv M_w/M_n$$

Degrees of polymerization

- DP_n : Number average degree of polymerization

$$DP_n \equiv M_n/R$$

- DP_w : Weight average degree of polymerization

$$DP_w \equiv M_w/R$$

- DP_z : Third order average degree of polymerization

$$DP_z \equiv M_z/R$$

Relative abundances

- $\%S_i$: Percentage of series of i-th entry

$$\%S_i \equiv 100 \cdot n_i / \sum_i n_i$$

- $\%S_i$: Percentage of spectrum of i-th series entry

$$\%S_i \equiv 100 \cdot n_i / \sum_j \left(\sum_i n_i \right)_j$$

- $\%W_i$: Weight percentage of series of i-th entry

$$\%W_i \equiv 100 \cdot (n_i \cdot M_i) / \sum_i (n_i \cdot M_i)$$

- $\%W_i$: Weight percentage of spectrum of i-th series entry

$$\%W_i \equiv 100 \cdot (n_i \cdot M_i) / \sum_j \left(\sum_i (n_i \cdot M_i) \right)_j$$

Miscellaneous

- Formula and mass of i-th series entry

HOMOPOLYMER ANALYSIS

- Monoisotopic peak error for i-th series entry, the difference between the m/z value where the most abundant matched peak occurs and its calculated value
- Cluster error for i-th series entry, the difference between the intensity weighted mass average for a matched entry and its calculated value
- Totals by series and by spectrum

Where:


M_i = mass of i-th series member

n_i = moles (\approx relative abundance) of i-th series peak

R = mass of repeat unit

The Homopolymer Analysis page user interface

Switch to the **Homopolymer Analysis** page by clicking the **Homopolymer Analysis** button on the **Workspace Control Panel**. The **Homopolymer Analysis** page is shown below after the data processing steps have been completed. For the PEG example spectrum shown, the data processing options included limiting the mass range to between 800 and 2000 Da and applying peak detection and baseline correction.

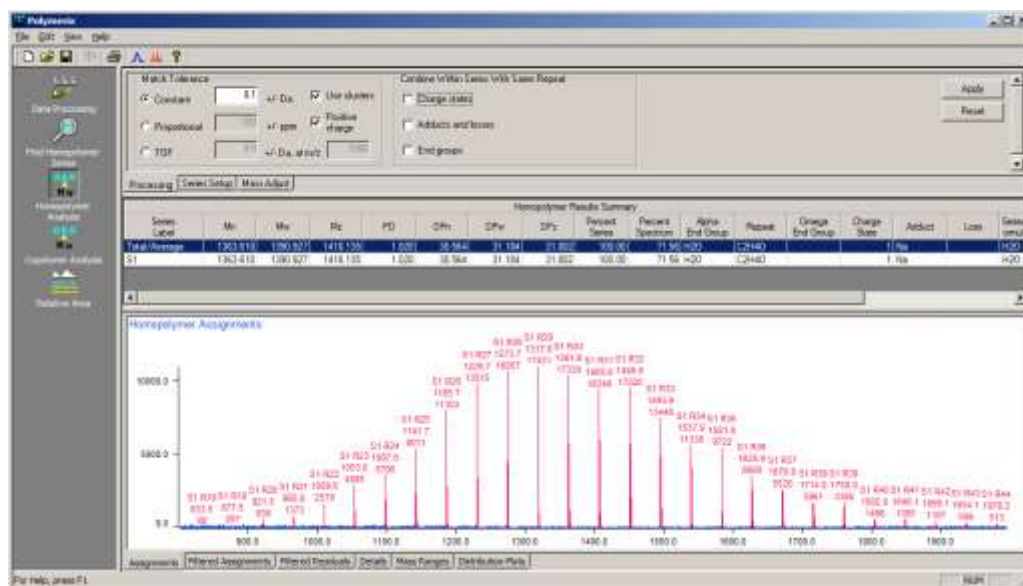

At the top of the page are the parameter setting tabs. There are three of these, a **Processing** tab for setting peak matching tolerances and options for combining related homopolymer series, a **Series Setup** tab used to define the homopolymer homologous series that will be matched as part on the analysis, and a copy of the **Mass Adjust** tab described previously in the **Data Processing** chapter.

In the middle of the page is a table of the homopolymer molecular weight distribution statistics. These are calculated for the sample as a whole and for each series individually.

At the bottom of the page there is a series of tabbed panes showing the assignment results in various graphical and tabular formats. Each of these will be described in more detail in the sections below.

Setting homopolymer processing parameters

Homopolymer analysis parameters are spread among three tabs at the top of the **Homopolymer Analysis** page.


The screenshot shows a software interface with two main panels. The left panel is titled 'Match Tolerance' and contains three radio button options: 'Constant' (selected) with a value of 0.1 and unit '+/- Da', 'Proportional' with a value of 100 and unit '+/- ppm', and 'TOF' with a value of 0.5 and unit '+/- Da, at m/z' and a sub-value of 1000. There are also two checked checkboxes: 'Use clusters' and 'Positive charge'. The right panel is titled 'Combine Within Series With Same Repeat' and contains three unchecked checkboxes: 'Charge states', 'Adducts and losses', and 'End groups'. At the bottom, there are three tabs: 'Processing' (selected), 'Series Setup', and 'Mass Adjust'.

Match tolerance

1. The first tab, **Processing**, contains parameters to control the tolerance for matching m/z values predicted from the homopolymer series definition with experimental peaks in the target spectrum.
 2. First, the **Match Tolerance** parameter should be set to correspond to the instrument scan function type. As in peak detection, these choices are:
 - Constant, used for quadrupole and ion trap instruments where the peak width is constant or nearly so over the mass range of interest. This choice also works well for very high resolution data such as is produced by FT-ICR or Orbitrap instruments,
 - Proportional, used for magnetic sector instruments, and
 - TOF, used for time of flight instruments.
20. Select the appropriate scan type, and then set the matching tolerance appropriately. In general, for well-resolved, peak detected spectra where there are now very closely-spaced peaks, the match tolerance can be set to any reasonable value that would not result in a tolerance window so wide as to cause mismatches. In the example spectrum, peaks are separated by 44 Da, so any window narrower than that is acceptable. Typically, a value of 0.5 Da or less (or the equivalent in ppm) is acceptable.

HOMOPOLYMER ANALYSIS

21. The **Use clusters** option should be selected when the target spectrum has not been deisotoped or has not been peak detected and it is desirable to match all peaks belonging to the target isotope cluster. With this option selected and continuum data, the monoisotopic peak will be matched, then the window will be broadened to match the additional peaks in the isotope cluster. With this option selected and centroided data, individual peaks within computed isotope clusters will be matched using the tolerance. In the example below an isotope cluster is shown. The match tolerance is set to 0.1 Da, which matches the isotope cluster peaks but not the peaks which lie between isotope cluster peaks.

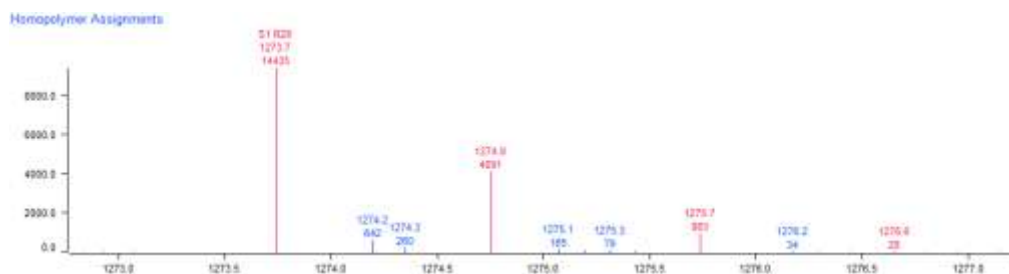

22. The **Positive charge** option should be selected to match the polarity of the spectrum being analyzed. It controls whether Polymerix adds or subtracts electron masses when computing the m/z values for a polymer series.

Combining series


23. **Combine Within Series With Same Repeat:** Next are optional choices for combining homopolymer series. In many cases, multiple series are found in the spectrum simply due to sampling or ionization artifacts. For example, the appearance of multiple charge states in ESI, or of both proton and sodium adducts is simply a result of the ionization process and not a true reflection of the sample. In this case, Polymerix provides options to combine the results for these different series into a composite result. In all cases though, the series to be combined must have the same repeat unit. The choices are:

- **Charge states:** By selecting this option, results for series with the same repeat unit, end group, and adduct that differ only by charge state will be combined into a single composite series result.
- **Adducts and losses:** By selecting this option, results for series with the same end group, repeat unit, and charge state that differ only by an adduct or a neutral loss will be combined into a single series results.
- **End groups:** Finally, by selecting this option, results for series that differ only by end group are combined into a single composite result.

24. Any combination of these options can be chosen, depending on the nature of the sample and ionization artifacts. With nothing selected, all series will be assigned and reported separately.
25. Once the match tolerance and series combination options have been chosen, click the **Apply** button to set them.

Defining homopolymer series

A homopolymer series consists of three parts:

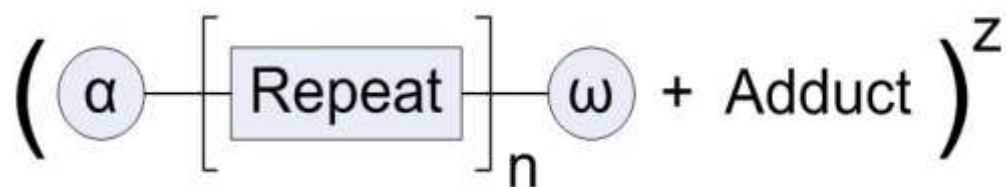

- An *alpha end group*, which starts the polymer chain;
- Some number of *repeat units* or *monomers*; and
- An *omega* end group that terminates the chain.


In addition, the process of sample introduction and ionization may result in modification of this core structure, including addition of one or more adducts and multiple charge states, or a neutral loss.

For each combination of end group, repeat unit, adduct, loss and charge state, Polymerix builds a list of target species to be matched against the processed target spectrum. In the nomenclature used by Polymerix this list of targets is called a *series*.

Polymerix does not know in advance what target lists to build; there is no built-in database of polymers and their chemical and mass spectral properties. You must specify everything Polymerix is to look for in the spectrum, by defining homopolymer series.

The **Find Homopolymer Series** page introduced in version 3.0 of Polymerix provides an alternative means of defining homopolymer series in which the user can specify the details of a series definition by selecting from lists of chemical compositions generated by Polymerix based upon polymer spectral data. Resultant series definitions can be transferred from the **Find Homopolymer Series** page. Using the **Find Homopolymer Series** page it takes five mouse clicks and no keystrokes to set up the example which follows.

HOMOPOLYMER ANALYSIS

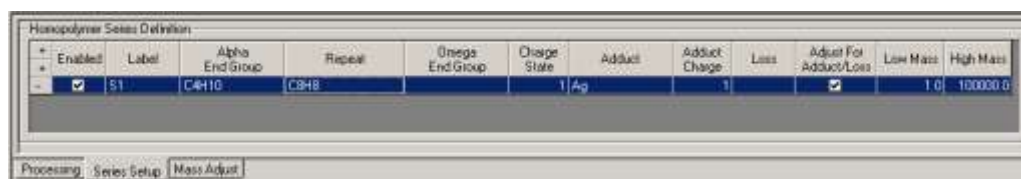

1. Click on the **Series Setup** tab of the **Homopolymer Analysis** parameters pane. With no series defined, it should look as shown above.
26. To start, *enable* the first series by selecting the **Enabled** check box to the left of the first row. Next, if desired, change the name of the new series by clicking in the **Label** box and typing a new name. By default, Polymerix entitles each series as S1, S2, S3, and so forth. You may change these to more descriptive names.




27. Next, click in the **Alpha End Group** box to display the **Specify End Group** dialog box, shown above. If the desired alpha end group appears in the list, you may select it, and then click the **OK** button.
28. Otherwise, you may define a new end group by clicking the **New** button. This will display the **Edit End Group** dialog box, which was described in an earlier chapter. End groups defined in this way will be added to the list for future use.
29. Alternatively, you may temporarily specify the end group either by chemical formula or by mass. Using a chemical formula is **strongly** preferred, because then Polymerix can calculate monoisotopic, chemical,

HOMOPOLYMER ANALYSIS

or nominal mass to high precision. If you do not know a chemical formula and must enter a mass, do so to the highest precision possible (as many decimal places as can be experimentally determined).

30. Click the **OK** button to close the dialog box and place the selected end group information in the series definition.

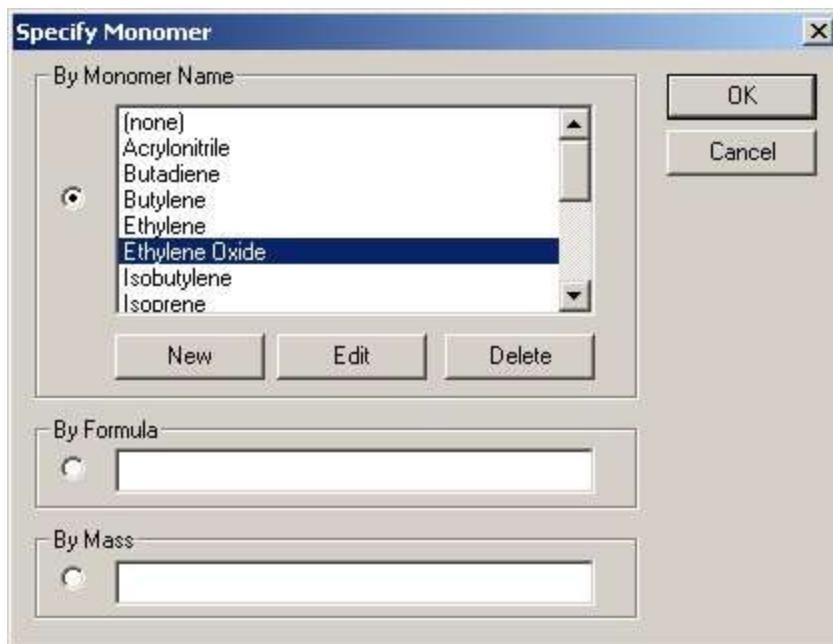


31. Next, click in the **Repeat** box to display the **Specify Monomer** dialog box. As with end groups, there is a predefined list, an option to create a new monomer definition, and options to enter a temporary definition by either chemical formula or mass. If you will be using Polymerix frequently on the same class of polymer, it is very convenient to add new monomer definitions so they will be available for future use. After defining the monomer, click the **OK** button to close the dialog box and copy the monomer into the series definition.
32. For the **Omega End Group**, follow the same procedure as for the alpha end group selection. If you know only the total end group, you may specify this as either an alpha or omega end group and leave the other entry empty.
33. Next, specify the charge state (a signed quantity) by typing into the **Charge State** box.. If more than one charge state is observed, you must define additional homopolymer series to match each of those.
34. Next, specify an adduct, if any. Click in the **Adduct** box to show a button; click the button to display the **Specify Adduct** dialog box. As with other chemical entities, there is a predefined list of common adducts, an option



HOMOPOLYMER ANALYSIS

to define new ones, and options to temporarily specify the adduct by formula or mass. After choosing the adduct, click the **OK** button to copy it into the definition.

35. Specify **Adduct Charge**, the charge state corresponding to the adduct. Polymerix uses **Adduct Charge** and **Charge State** to compute the number of adducts used to create a molecular ion.
36. A few more options remain in the series definition. If a loss occurs in sampling or ionization, specify that by formula or mass by typing in the **Loss** box. If you wish to have the molecular weight distribution statistics adjusted for the mass difference of the adduct or loss, click the **Adjust For Adduct / Loss** check box. Finally, to restrict the mass range searched for the newly-defined series edit the range by typing into the **Low Mass** or **High Mass** boxes.
37. This completes the definition of the first series. Click the **Apply** button.
38. To define additional series, click the **+** button next to the **Enabled** box. This will make a copy of the last series onto a new line. Make whatever changes are necessary, then click the **Apply** button again. Continue until all needed series are defined.
39. To remove a series definition, click the **-** button in the column to the left of the **Enabled** box.
40. At this point, it might be a good idea to save the definitions into a template file for later use.

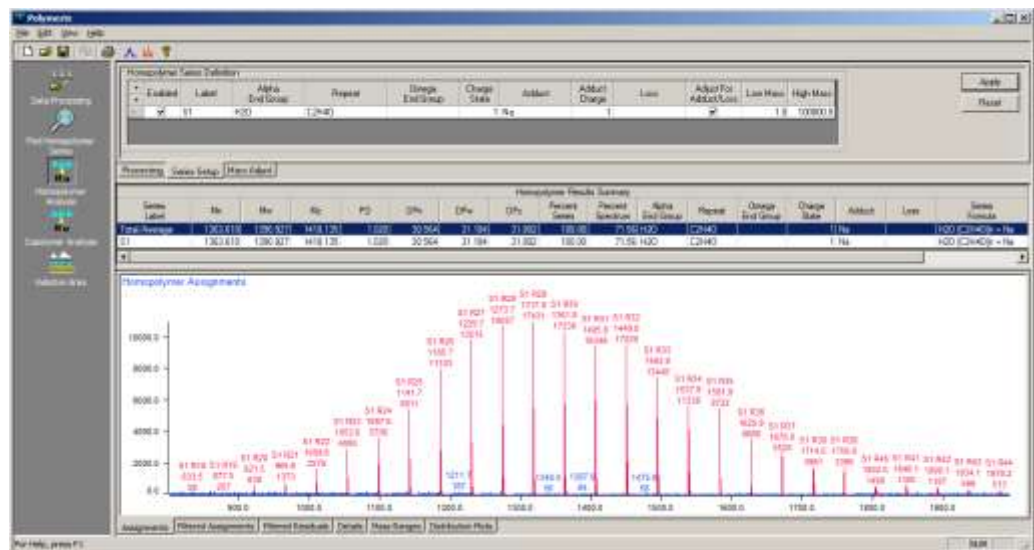

HOMOPOLYMER ANALYSIS

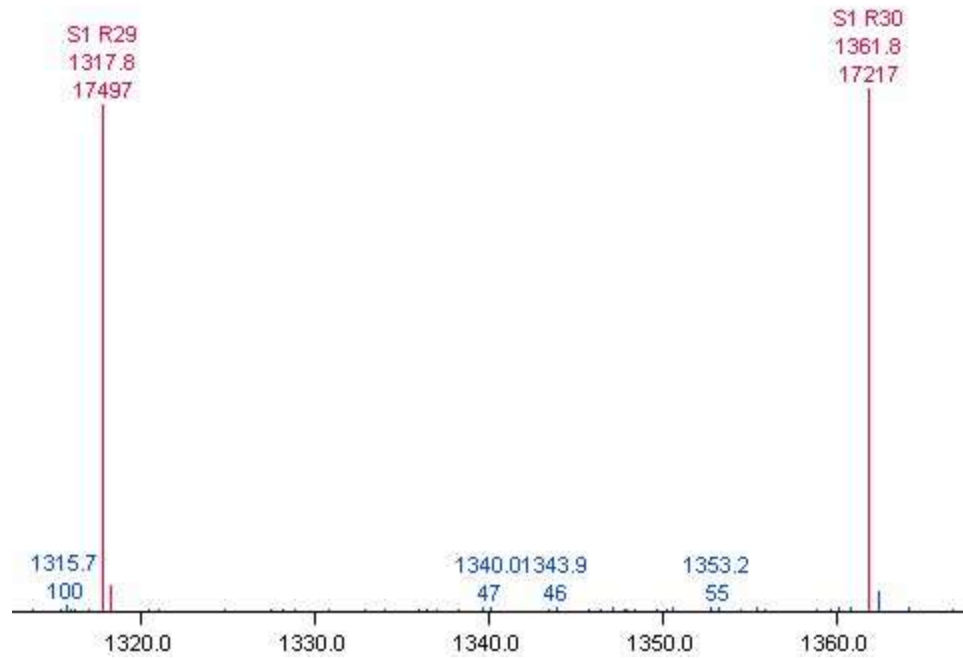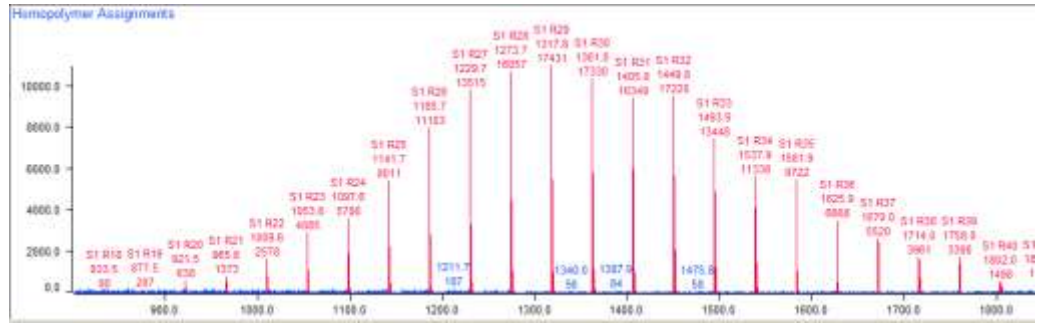

With one series defined, our example looks like the screen shots above. The series definition has been expanded into a list of targets appropriate for the 800 – 2000 Da mass range chosen, and these have been matched against the spectral features. Matched peaks have been labeled and colored. The labels show the series name, (S1), repeat count (Rn), m/z, and intensity.

Homopolymer molecular weight statistics

When homopolymer series are defined and the **Apply** button clicked, Polymerix immediately calculates molecular weight statistics for the whole spectrum as well as for each series. These are displayed in the middle pane of the **Homopolymer Analysis** page.

Series Label	Rn	Rm	Mt	RD	CPn	DPn	DPm	DPw	Percent Series	Percent Structure	Alpha End Group	Repeat	Charge End Group	Charge Size	Added	Loss	Series Formula
Total Average	183.33	138.48	1413.28	1.03	30.48	25.02	21.03	100.00	68.84	24	C2H43	01					(C2H43)n + 8a
S1	183.33	138.48	1413.28	1.03	30.48	25.02	21.03	100.00	68.84	24	C2H43	01					(C2H43)n + 8a

HOMOPOLYMER ANALYSIS

Series Label	Mn	Mw	Mz	PD	DPn	DPw	DPz
Total/Average	1359.331	1386.488	1413.386	1.020	30.466	31.083	31.694
S1	1359.331	1386.488	1413.386	1.020	30.466	31.083	31.694

Percent Series	Percent Spectrum	Alpha End Group	Repeat	Omega End Group	Charge State	Adduct	Loss	Series Formula
100.00	68.84	OH	C2H4O	H	1	Na		OH [C2H4O] _n H + Na
100.00	68.84	OH	C2H4O	H	1	Na		OH [C2H4O] _n H + Na


This table starts with a summary row for the entire sample, followed by a row for each series. If one or more of the options to combine series was selected during setup, the combined series will be shown instead of individual series (e.g. S1 + S2). The left side of the table presents the statistics, and the right side summarizes the series definition.

Visualizing homopolymer series assignments

There are four tabs in the bottom pane of the **Homopolymer Analysis** page, each presenting the assignment details in a different way.

Assignments


The **Assignments** tab presents the target spectrum with all matched series peaks labeled. This can be a cluttered plot if there are many series with many assigned peaks. However, it shows at a glance if there are major unassigned peaks in the spectrum. Here we have defined an additional series, with a proton adduct.

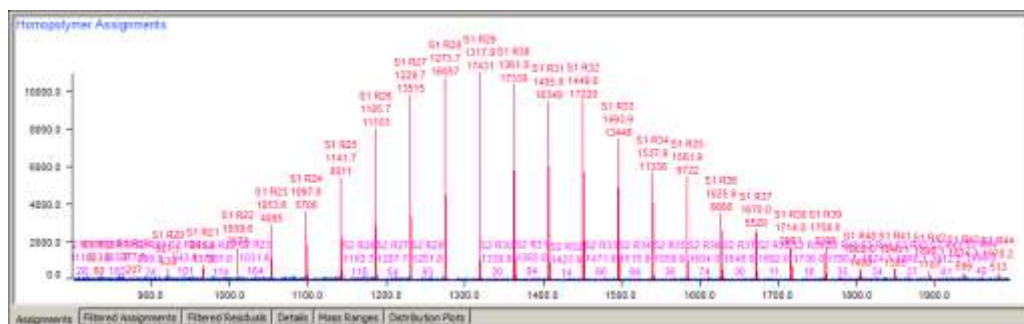

Filtered Assignments


The **Filtered Assignments** tab shows only the assigned peaks in the spectrum.

HOMOPOLYMER ANALYSIS

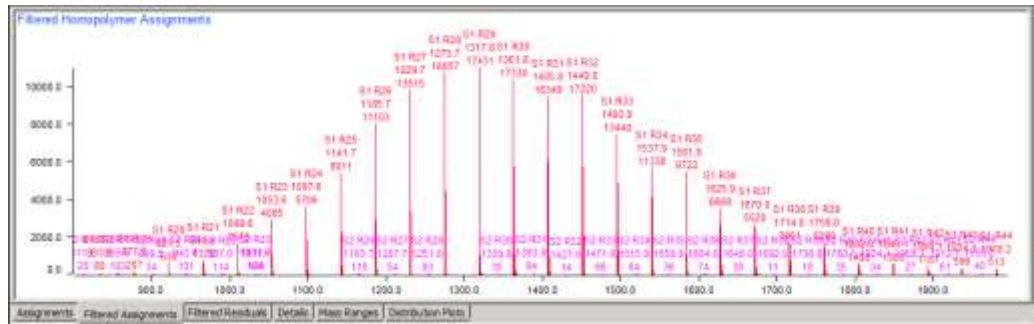

Filtered Residuals

The **Filtered Residuals** tab shows only the peaks in the spectrum which have not been assigned to a series. In the current example, since the peaks corresponding to the polymer have been matched, all that remains in the spectrum is low intensity noise.

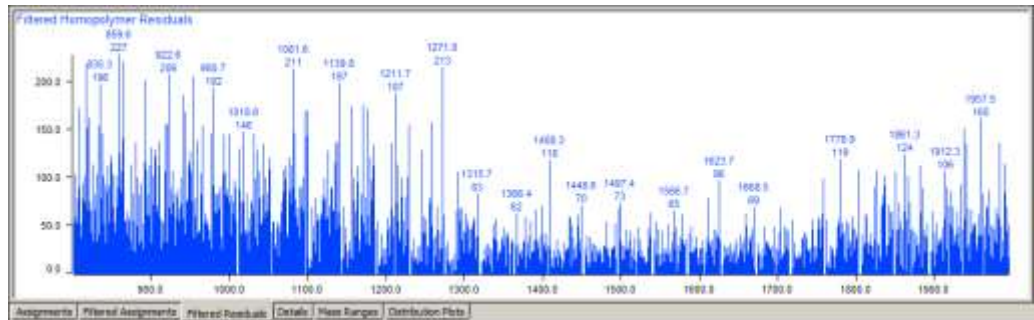

Details

This presents a table containing quantitative details of every peak assignment for all series. The left side of the table summarizes assignments for all series, followed by additional groups of columns giving assignment details for each series, by repeat unit count. Depending upon the data being analyzed and the parameters being used for quantitative analysis, columns for “Homopolymer m/z” and “Cluster m/z” may or may not be shown.

Homopolymer Analysis Details										Homopolymer Analysis Details									
Totals										S1									
Repeat Count	Found Intensity	% Found Intensity	% Found Weight	Percent Spectrum	Percent Series	S Series Weight	Percent Spectrum	Monomeric m/z			Cluster m/z			Found Intensity	Percent Series	% Series Weight	Percent Spectrum		
								Computed	Found	Err (mDa)	Computed	Found	Err (mDa)						
18	338.78	0.2825	0.0337	0.0408	0.3421	0.3251	0.0332	823.472			833.538	834.507	-0.969	38.46	1.5732	1.1303	0.0108		
19	389.56	0.32625	0.1271	0.1485	2.3352	0.2678	0.074	877.488	-11.554	877.995	877.896	-0.099	206.51	12.1903	8.1692	0.0896			
20	472.01	0.3982	0.2504	0.2918	0.3343	0.2282	0.2290	923.524	821.572	-101.952	923.038	923.976	-0.938	437.83	2.7398	1.6122	0.0128		
21	1473.75	0.7689	0.5380	0.5524	0.7190	0.4973	0.5145	365.956	365.961	0.005	366.089	366.091	0.002	1172.81	6.6854	4.9607	0.0379		
22	2692.32	1.3992	1.0134	1.0592	1.3505	0.8776	0.9864	1009.577	1009.609	0.032	1010.139	1010.044	-0.095	2576.22	7.6461	5.8771	0.0428		
23	4248.57	2.2079	1.6793	1.9326	2.1397	1.6177	1.6312	1953.600	1953.689	0.089	1954.189	1954.348	0.159	4384.97	10.6201	8.7998	0.0613		
24	5796.30	2.9655	2.2381	2.7392	2.9890	2.3946	2.1390	1987.629	1987.525	-0.104	1988.239	1988.078	-0.161	5706.30					
25	9816.88	4.9520	3.6183	4.0527	4.7963	3.4435	3.0927	1541.658	1541.554	-0.104	1542.268	1542.093	-0.175	8816.89					
26	11220.26	5.6111	4.2751	4.2258	8.8184	4.9652	4.1617	1185.881	1185.884	0.003	1186.337	1186.229	-0.108	11102.52	7.7994	7.1378	0.0441		
27	15669.09	7.8347	6.2454	6.6863	7.8791	6.2642	5.0859	1229.708	1229.713	0.005	1230.386	1230.325	-0.061	12514.77	5.9322	3.4218	0.0204		
28	16150.18	8.0931	7.7052	6.8538	8.4113	7.7174	6.0190	1273.754	1273.741	-0.013	1274.426	1274.154	-0.272	16027.30	6.1380	6.0536	0.0348		
29	17430.93	8.9587	8.0090	6.5338	9.1304	8.6725	6.5339	1317.762	1317.761	-0.001	1318.518	1318.207	-0.311	17430.93					
30	17360.06	8.8219	8.0553	6.5272	9.0776	8.9158	6.4961	1361.796	1361.799	0.003	1362.565	1362.292	-0.273	17330.96	1.8846	2.0894	0.0112		
31	16433.97	6.5490	6.6893	6.1803	8.9543	8.8870	6.1325	1405.812	1405.817	0.005	1406.620	1406.318	-0.302	16349.50	5.9885	6.1904	0.0317		

Mass Ranges

The **Mass Ranges** table displays a list of the calculated m/z ranges used for determining series assignments ordered by m/z value. The table lists the m/z window


HOMOPOLYMER ANALYSIS

(as defined by the matching constraints), series name, repeat unit count, and the number of isobaric overlaps (collisions) if any.

Low Mass	High Mass	Series	Repeat	Collisions	
811.3957	874.5304	S2	18		
833.3717	836.5304	S1	18		
855.4155	858.5247	S2	19		
877.3979	880.8986	S1	19		
899.4421	882.6509	S2	20		
921.4341	924.8329	S1	20		
943.4604	946.6772	S2	21		
965.4903	968.8591	S1	21		
987.4946	990.7034	S2	22		
1009.4765	1012.8854	S1	22		
1031.5208	1034.7297	S2	23		
1053.5027	1056.7116	S1	23		
1075.5470	1078.7958	S2	24		
1097.5289	1100.7379	S1	24		
1119.5732	1122.7822	S2	25		
1141.5552	1144.7641	S1	25		
1163.5994	1166.8084	S2	26		

Distribution Plots

This graphical view is linked to the **Homopolymer Results Summary** table. Selecting a row in the table displays the peaks assigned to the series defined in that row. In the first illustration below, the first series (S1) is selected; in the second series (S2) is selected.

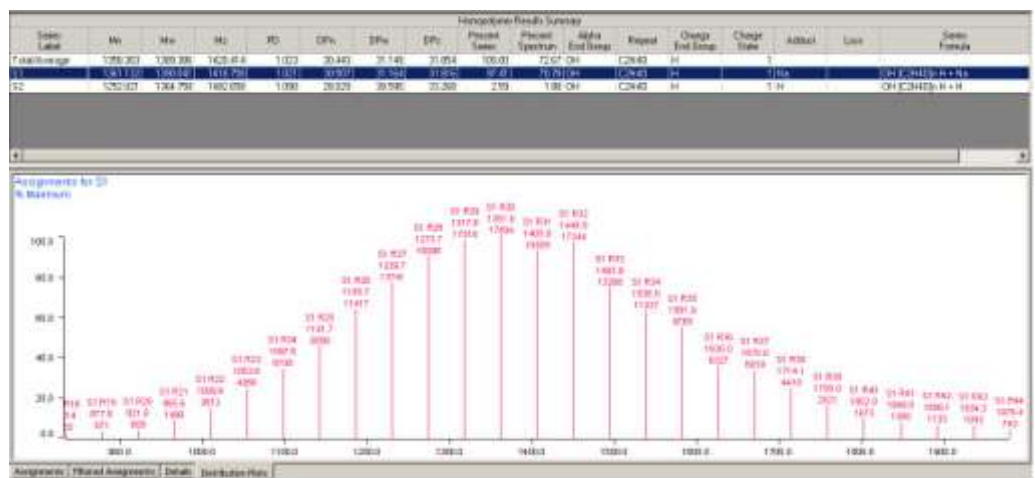

Assignment and matching details

The algorithms which Polymerix uses for quantitative analysis vary depending upon the data being analyzed and the settings being used for analysis. When multiple series have been defined and there are series overlaps, the different algorithms assign intensity values to series differently.

At its highest level, mass matching can be divided into matching based upon single masses versus matching based upon computed isotope clusters. When computed isotope clusters are being used for matching, centroided data is matched differently than continuum data.

Matching based upon single masses and matching of continuum data based upon isotope clusters share the same algorithm for making intensity assignments and resolving intensity assignments in overlapping series. Isotope cluster based matching of centroid data uses a different algorithm for resolving intensity assignments in overlapping series.

The simple matching algorithm

This algorithm is used when the **Use clusters** option is not selected, or when **Use clusters** is selected and the data being analyzed is continuum data. The steps in the algorithm are as follows:

1. Use polymer series definitions and the match tolerance to generate mass ranges corresponding to each series. Each mass range corresponds to a distinct repeat count or, for copolymers, a combination of repeat A and repeat B counts.
2. Partition the mass ranges into those without m/z collisions and those with m/z collisions.
3. Assign intensity values from the spectrum being analyzed to the collision-less mass ranges.
4. Construct absolute intensity estimates for the mass ranges which have not yet been assigned using the assignments determined in the previous and linear interpolation and extrapolation.
5. Assign intensity values from the spectrum being analyzed to the collided mass ranges. The intensity estimates from the previous step are used to determine a proportional (to the estimates) intensity distribution to each mass range for a spectrum data point which matches multiple mass ranges.

Cluster matching

When the **Use clusters** option is enabled, Polymerix will calculate theoretical isotope clusters corresponding to a series, and use these theoretical clusters together with the match tolerance to construct the mass ranges described above. The low m/z value for


HOMOPOLYMER ANALYSIS

a mass range corresponds to the monoisotopic peak of its isotope cluster. The high m/z value for a mass range corresponds to the highest mass peak in the isotope cluster for which the peak's intensity is greater than or equal to 2% relative abundance.

Centroid cluster matching

When the **Use clusters** option is enabled and centroided data is being analyzed, the assignment process is similar to that used for simple matching, but with an important difference – the individual peaks in theoretical isotope clusters are used for matching, in addition to the mass ranges described above. With small match tolerance values, this excludes some intensity values from being assigned to a series.

The example below shows an isotope cluster. Because the match tolerance is set to 0.1 Da, only the isotope cluster peaks are matched. Match contributions from data between the isotope cluster peaks are filtered out.

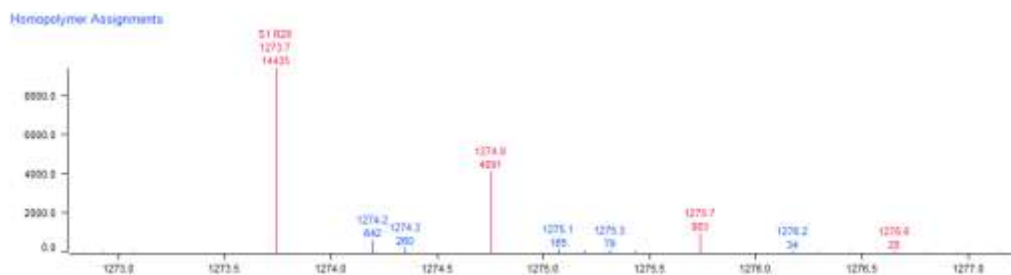

Centroid cluster matching algorithm

The steps used for centroid cluster matching are as follows.

1. Determine mass ranges as described above.
2. Assign intensity values from the spectrum being analyzed to different series, based upon mass ranges, theoretical isotope clusters, and the match tolerance. When a data point in the spectrum being analyzed matches multiple series, its intensity values are equally distributed to each of the matching series.

Centroid cluster matching also does special accounting for theoretical isotope cluster monoisotopic peaks. The m/z value for the most abundant peak which matches each monoisotopic peak is determined and used for calculating error values.

When series assignments are shown in data plots, the series annotations are positioned at m/z values corresponding to computed monoisotopic masses.

What was covered in this chapter

In this chapter, we

HOMOPOLYMER ANALYSIS

- Described the statistics calculated by Polymerix to describe homopolymer series.
- Showed how to define a homopolymer series and set matching parameters.
- Explained the result tables and plots produced by Polymerix in the series assignment process.
- Explained the algorithms used for quantitative analysis.

In the next chapter

In the next chapter, we will present the procedures for defining copolymer series and visualizing copolymer assignment results.

Copolymer Analysis

Defining and reviewing copolymer series assignments


The definition and analysis of copolymer series is very similar to homopolymer analysis, with one crucial difference. Instead of one repeating unit, there are two, and as a consequence the assignments for a series form a two-dimensional grid. Keeping the count of one repeat unit fixed, the second repeat unit typically has a repeat distribution similar to a homopolymer. Increment the first repeat unit count, there is another distribution for the second repeat unit, and so forth.

Goals for this chapter

In this chapter, we will:

- Describe Polymerix's definition of copolymers
- Present the **Copolymer Analysis** page and how to use it.
- Show how to define copolymer series.
- Explain the tables and plots produced by Polymerix during copolymer series assignment.

Copolymer definition

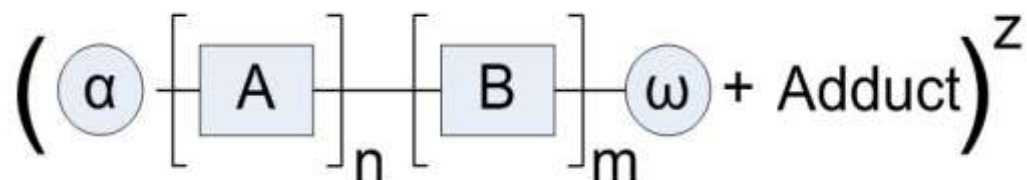


COMPOLYMER ANALYSIS

In the Polymerix world, a copolymer series is a polymer containing two different repeat units, which we label as **A** and **B** in the schematic above. Like a homopolymer, a copolymer series also has alpha and omega end groups, an adduct and/or neutral loss, and a charge state. In a mass spectrometric determination of a copolymer molecular weight distribution, there is information only about the total m/z value of a given peak. Nothing can be inferred about the copolymer sequence responsible for the observed m/z value. It could be a block, alternate, or random copolymer or something else. All that is known is that, after subtracting out the end group and adduct masses, there are n counts of repeat unit A and m counts of repeat unit B.

However, as will be shown in the description of copolymer series setup below, you can add your knowledge of the chemistry to help guide the assignment process, by putting constraints on the type of copolymer and ranges of A and B counts. This may serve to eliminate any possibility of isobaric overlaps or at least minimize them to compositions that are chemically meaningful.

The Copolymer Analysis page user interface

Switch to the **Copolymer Analysis** page by clicking the **Copolymer Analysis** button on the **Workspace Control Panel**. The **Copolymer Analysis** page is shown below, after the data processing steps have been completed.

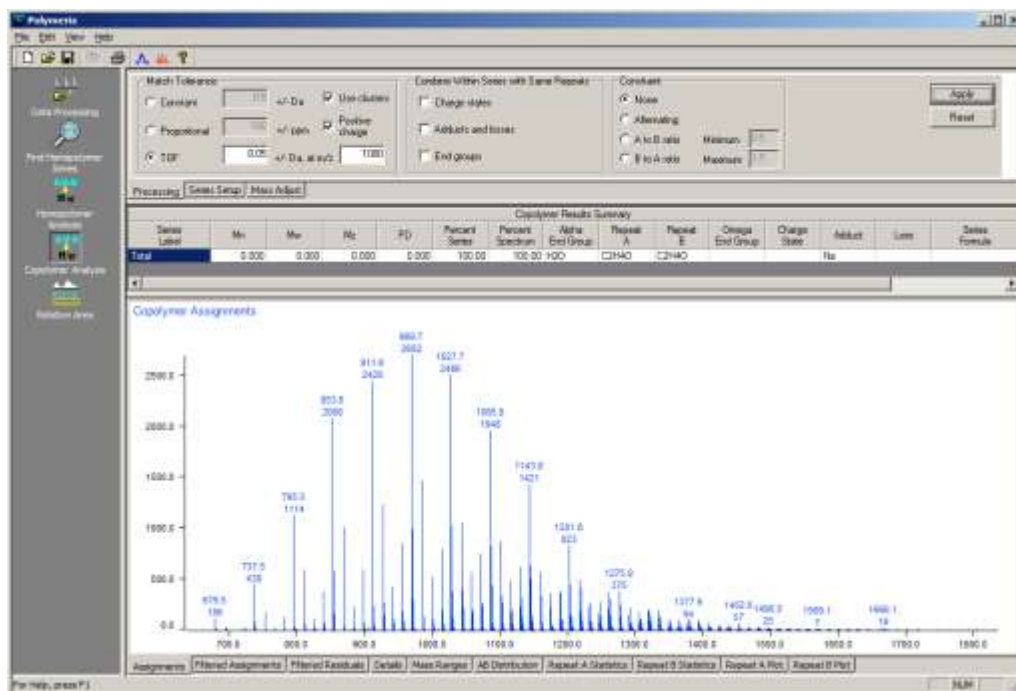

As with the **Homopolymer Analysis** page, there are three panes; the upper pane contains several tabs of copolymer analysis parameters, the middle pane displays molecular weight distribution statistics, and the bottom pane is a tabbed display of plots and tables with different views onto the assignment results.

The analysis parameters pane contains a copy of the **Mass Adjust** tab. This is a copy of the same tab in the **Data Processing** page and will not be described further here.


Setting copolymer processing parameters

In addition to the same match tolerance and series combination options as are available for homopolymer analysis, there is an additional parameter set used to define the type of copolymer and constrain the number of A and B combinations.

- To set a constraint on the relative counts of A and B units, click the **Processing** tab. On the right side is a new section, **Constraint**. There are four constraints:
 - None:** The A and B monomer units may occur in any combination.
 - Alternating:** The copolymer is strictly alternate, and the A count = B count ± 1 for all assignments. This allows for one extra A or B unit at the end of the chain. If the A and B counts must be strictly equal, use the ratio constraint below instead.
 - A to B ratio:** The ratio of A count to B count must fall within a specified range. When either this or the next option is selected, the **Minimum** and **Maximum** edit boxes are enabled. Enter the minimum and maximum values of the A/B ratio.
 - B to A ratio:** This is simply the inverse of the A to B ratio and is provided for convenience.
- After setting the desired matching, combination, and constraint parameters, click the **Apply** button.

Defining copolymer series

As was mentioned above, the major difference between a homopolymer and a copolymer series definition is the presence of a second repeat unit.

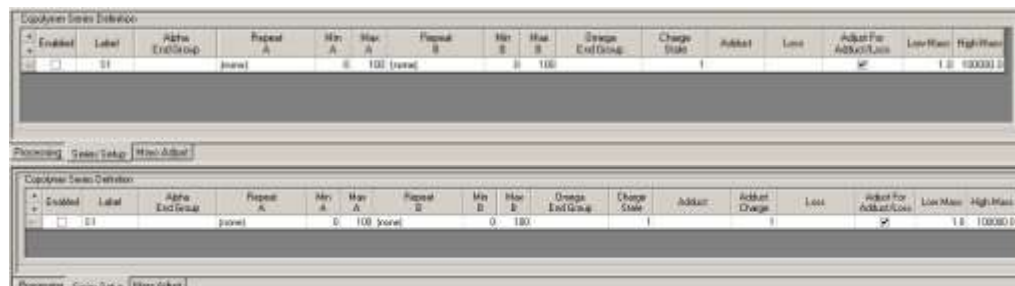

1. Click the **Series Setup** tab to view the copolymer series definition table.

Copolymer Series Definition						
+	Enabled	Label	Alpha End Group	Repeat A	Min A	Max A
--	<input type="checkbox"/>	S1		(none)	0	100

Repeat B	Min B	Max B	Omega End Group	Charge State	Adduct	Adduct Charge	Loss
(none)	0	100		1		1	

2. Most of the editable components in this table are the same as for homopolymer series setup, and behave in the same way: click in a box, click the button, and then select or define a chemical entity from the dialog.
3. As can be seen from the screen shots above, there are a few new items as well. First, instead of a single repeat unit, there are now separate selections for A and B repeats. Next, there are lower and upper limits on the A repeat count and B repeat count (set by typing into the **Min A**, **Max A**, **Min B**, and **Max B** edit boxes). These repeat count constraints are used in conjunction with the ratio constraints on the **Processing** tab to limit the number of combinations of A and B count generated for matching to the spectrum.
4. For example, suppose the **A to B ratio** constraint was chosen with minimum and maximum ratio both set to 2. This means that for every B, there must be two A units. So, if the A and B count limits are each set from 1 – 100, then the only (A,B) counts that will actually be considered are (2,1), (4,2), (6, 3),..., (100,50) since no others satisfy both the ratio and count range limits.

COMPOLYMER ANALYSIS

5. After defining the necessary series, click the **Apply** button to assign the spectrum.

Copolymer molecular weight statistics

The table of copolymer molecular weight distribution statistics is identical to that created for homopolymers, with the substitution of **Repeat A** and **Repeat B** columns for the single **Repeat** column.


Series Label	M _n	M _w	M _v	PD	Percent Series	Percent Spectrum	Alpha	Repeat A	Repeat B	Charge	Charge State	AdSet	Lot	Series Formula
Total Average	1038.026	1059.088	1073.097	1.020	100.00	95.77	0.4	C2H4O	C3H6O	H		1 Na		CH ₂ (C2H4O) _x (C3H6O) _y H + Na
BT	1038.026	1059.088	1073.097	1.020	100.00	95.77	0.4	C2H4O	C3H6O	H		1 Na		CH ₂ (C2H4O) _x (C3H6O) _y H + Na

The statistics are computed in the same way as for homopolymers.

Visualizing copolymer series assignments


Because copolymer assignments are a two-dimensional result, the results are presented in a two-dimensional manner. Since assignments are deconvolved into separate distributions by A and B count, some additional tables and plots are provided to show the distribution of each repeat unit as a composite and separately.

Assignments

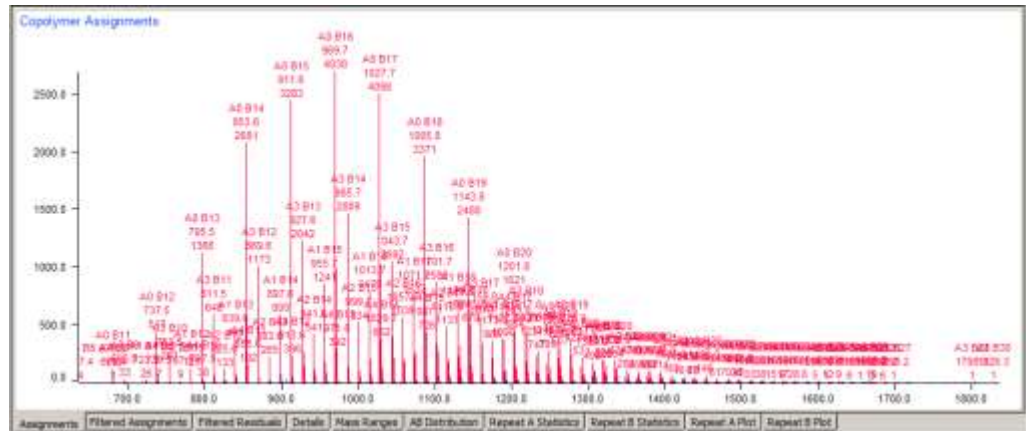


This is basically the same plot as for homopolymer analysis, showing the entire target spectrum with all assigned peaks labeled. Instead of an R_n repeat count label, there is an A_nB_m label. A detail view of the screen shot above illustrates this:

COMPOLYMER ANALYSIS

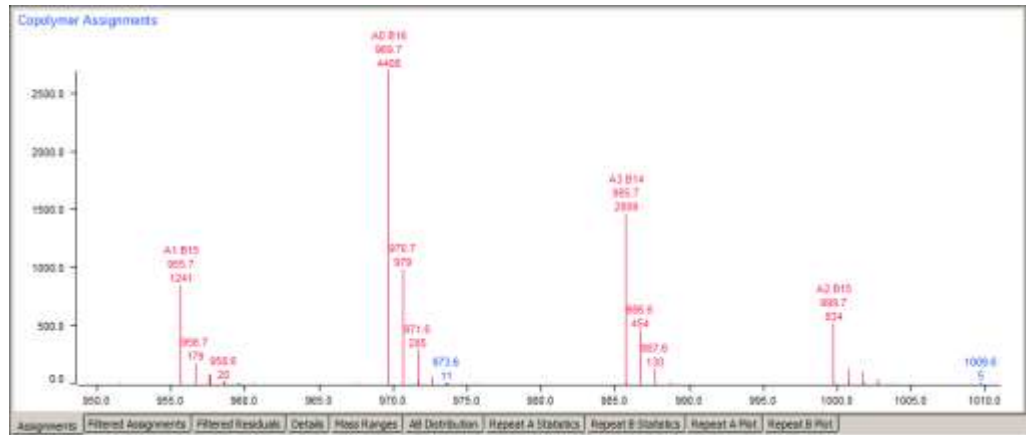

Filtered Assignments


This plot is also the same as for homopolymer analysis; only the assigned peaks are shown in the plot.

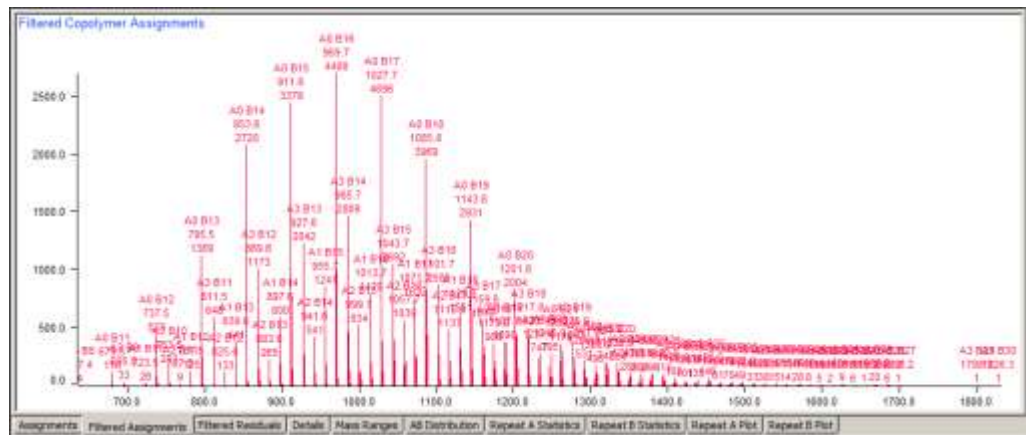

Filtered Residuals

This plot is also the same as for homopolymer analysis; only peaks which have not been assigned to any series are shown in the plot.

COMPOLYMER ANALYSIS

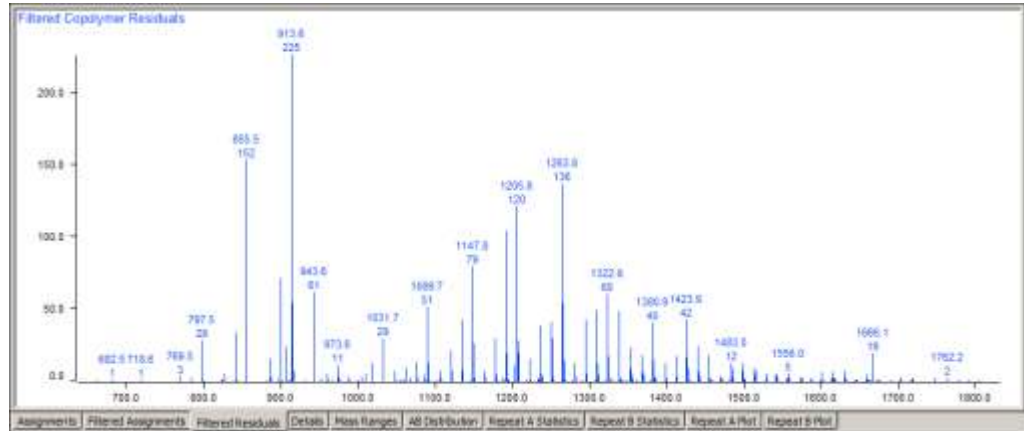

Details

The **Details** table groups the assignments by B count within A count. That is, the first group contains the range of B counts matched for the minimum A count, the next group contain the B counts for the next A count, and so forth.

Copolymer Analysis Details										Copolymer Analysis Details									
Totals										S1									
Repeat A Count	Repeat B Count	Found Intensity	% Found Intensity	% Found Weight	Percent Spectrum	Percent Series	% Series Weight	Percent Spectrum	Monoisotopic m/z			Cluster m/z			Found Intensity				
									Computed	Found	Error (mDa)	Computed	Found	Error (mDa)					
0	11	116.00	0.1831	0.1161	0.1737	0.1831	0.1161	0.1737	679.480	679.481	0.681	679.839	679.548	291.461	116.00				
0	12	523.00	0.8257	0.5698	0.7832	0.8257	0.5698	0.7832	737.502	737.503	0.826	737.946	737.690	205.993	523.00				
0	13	1369.00	2.1612	1.6128	2.0500	2.1612	1.6128	2.0500	795.544	795.544	0.123	795.623	795.736	287.440	1369.00				
0	14	2725.00	4.2940	3.4453	4.0731	4.2940	3.4453	4.0731	853.586	853.586	0.119	854.100	853.862	237.589	2725.00				
0	15	3379.00	5.3344	4.5794	5.0600	5.3344	4.5794	5.0600	911.620	911.620	0.654	912.176	911.967	208.603	3379.00				
0	16	4408.00	6.9589	6.3656	6.6009	6.9588	6.3656	6.6009	969.670	969.670	0.470	970.252	970.250	-1.361	4408.00				
0	17	4696.00	7.4135	7.1984	7.0322	7.4135	7.1984	7.0322	1027.711	1027.712	0.761	1028.327	1028.480	153.024	4696.00				
0	18	3969.00	6.2658	6.4258	5.9435	6.2658	6.4398	5.9435	1085.753	1085.754	0.741	1086.402	1086.623	220.848	3969.00				
0	19	2931.00	4.6271	5.0121	4.3891	4.6271	5.0121	4.3891	1143.795	1143.796	0.336	1144.476	1144.684	207.875	2931.00				
0	20	2064.00	3.1637	3.6047	3.0009	3.1637	3.6047	3.0009	1201.837	1201.837	-0.061	1202.951	1202.886	304.767	2064.00				
0	21	1176.00	1.8471	2.2088	1.7520	1.8471	2.2088	1.7520	1259.879	1259.880	0.977	1260.658	1261.332	674.382	1176.00				
0	22	747.00	1.1793	1.4768	1.1186	1.1793	1.4768	1.1186	1317.921	1317.921	0.244	1318.735	1319.746	1011.219	747.00				
0	23	296.00	0.4673	0.6116	0.4433	0.4673	0.6116	0.4433	1375.963	1375.965	2.617	1376.812	1378.242	1430.568	296.00				
0	24	135.00	0.2131	0.2909	0.2022	0.2131	0.2909	0.2022	1434.005	1434.004	-0.163	1434.899	1436.581	1691.991	135.00				
0	25	49.00	0.0774	0.1100	0.0734	0.0774	0.1100	0.0734	1492.046	1492.052	5.927	1492.965	1494.926	1969.365	49.00				
0	26	14.00	0.0221	0.0325	0.0210	0.0221	0.0325	0.0210	1550.088	-	-	1551.942	1552.366	1324.990	14.00				
0	27	2.00	0.0032	0.0048	0.0030	0.0032	0.0048	0.0030	1608.130	-	-	1609.118	1612.117	2968.899	2.00				
0	28	29.00	0.0336	0.0501	0.0299	0.0336	0.0501	0.0299	1666.172	1666.202	29.508	1667.193	1668.468	726.621	29.00				
1	11	26.00	0.0410	0.0276	0.0389	0.0410	0.0276	0.0389	723.487	723.488	1.967	723.886	723.726	-166.367	26.00				

Mass Ranges

The **Mass Ranges** table displays a list of the calculated m/z ranges used for determining series assignments ordered by m/z value. The table lists the m/z window (as defined by the matching constraints), series name, A and B unit counts, and the number of isobaric overlaps (collisions) if any.

COMPOLYMER ANALYSIS

Copolymer Mass Ranges					
Low Mass	High Mass	Series	Repeat A	Repeat B	Collisions
637.3734	639.4593	S1	3	8	
651.3886	653.4754	S1	2	9	
665.4039	667.4916	S1	1	10	
679.4191	681.5077	S1	0	11	
695.4135	697.5030	S1	3	9	
709.4287	711.5191	S1	2	10	
723.4440	725.5352	S1	1	11	
737.4592	740.5540	S1	0	12	
753.4537	756.5494	S1	3	10	
767.4689	770.5655	S1	2	11	
781.4842	784.5815	S1	1	12	
795.4994	798.5976	S1	0	13	
811.4939	814.5929	S1	3	11	
825.5092	828.6090	S1	2	12	
839.5244	842.6251	S1	1	13	
853.5397	856.6411	S1	0	14	
869.5342	872.6364	S1	3	12	
883.5495	886.6525	S1	2	13	
897.5647	900.6686	S1	1	14	
911.5800	914.6846	S1	0	15	
927.5745	930.6799	S1	3	13	
941.5898	944.6959	S1	2	14	
955.6051	958.7120	S1	1	15	

Assignments | Filtered Assignments | Filtered Residuals | Details | Mass Ranges | AB Distr

AB Distribution

This interesting table is a “top down” view of the A vs. B distribution. A repeat unit counts increase downward on the Y axis, while B counts increase along the X axis. For every AB combination where a matching peak was found, the table cell contains the relative abundance on the peak. In addition, the cells are color-coded; the more intense the color, the higher the relative abundance.

		Copolymer Distribution																																			
		Repeat B																																			
		8	9	10	11	12	13	14	15	16	17	18	19	20	21	22	23	24	25	26	27	28															
Repeat A	0				2.47	11.14	21.15	57.82	71.95	56.87	109.89	94.52	62.41	42.67	24.91	15.91	6.30	2.97	1.04																		
	1					2.84	9.39	17.04	26.43	31.49	34.75	33.03	23.21	16.93	10.96	6.09	2.15	1.49																			
	2						2.83	5.64	11.52	17.76	22.13	24.21	21.00	15.91	11.31	6.05	3.47	1.30																			
3				4.28	13.83	24.58	43.48	55.82	57.33	52.07	39.55	34.77	22.64	12.01	5.82	3.11																					

Assignments | Filtered Assignments | Filtered Residuals | Details | Mass Ranges | AB Distribution | Repeat A Statistics | Repeat B Statistics | Repeat A Plot | Repeat B Plot

COMPOLYMER ANALYSIS

Repeat A Statistics

On this tab is a table of molecular weight distribution statistics for all A unit series by fixed B count. That is, keeping the B count fixed, statistics are computed for every peak in the series with that B count and variable A count. This is repeated for the next B count, and so forth. This is the equivalent of taking each column in the AB distribution table and computing statistics for it.

Repeat B	Repeat A Statistics For Constant Repeat B Counts						
	Mn	Mw	Mz	PD	DPn	DPw	DPz
8	614.404	614.404	614.404	1.000	3.000	3.000	3.000
9	672.627	672.627	672.628	1.000	3.004	3.004	3.004
10	730.645	730.646	730.646	1.000	3.003	3.003	3.003
11	766.114	769.114	771.768	1.004	2.490	2.559	2.619
12	802.770	806.958	810.891	1.005	2.005	2.100	2.189
13	848.690	852.938	857.061	1.005	1.729	1.826	1.919
14	896.973	900.996	904.986	1.004	1.508	1.599	1.690
15	948.458	951.977	955.517	1.004	1.359	1.439	1.519
16	999.691	1002.842	1006.049	1.003	1.204	1.276	1.348
17	1050.523	1053.164	1055.880	1.003	1.040	1.100	1.162
18	1109.290	1111.758	1114.290	1.002	1.057	1.113	1.170
19	1165.494	1167.771	1170.110	1.002	1.015	1.067	1.120
20	1220.322	1222.280	1224.294	1.002	0.942	0.986	1.032
21	1277.573	1279.383	1281.247	1.001	0.924	0.965	1.008
22	1331.481	1333.054	1334.678	1.001	0.830	0.866	0.903
23	1384.301	1385.534	1386.810	1.001	0.712	0.740	0.769
24	1451.112	1452.491	1453.905	1.001	0.911	0.942	0.974
25	1503.499	1504.873	1506.293	1.001	0.782	0.814	0.846
26	1579.473	1581.214	1582.976	1.001	1.190	1.229	1.269
27	1621.900	1622.706	1623.519	1.000	0.835	0.853	0.872
28	1649.941	1650.448	1650.992	1.000	0.154	0.165	0.177
29	0.000	0.000	0.000	0.000	0.000	0.000	0.000
30	1807.244	1807.244	1807.244	1.000	1.090	1.090	1.090

Assignments | Filtered Assignments | Filtered Residuals | Details | Mass Ranges | AB Distribution | Repeat A Statistics | Repeat B S

Repeat B Statistics


Likewise, by slicing the AB distribution table by rows, the molecular weight distribution over B counts for each fixed A count can be computed.

Repeat B Statistics For Constant Repeat A Counts							
Repeat A	Mn	Mw	Mz	PD	DPn	DPw	DPz
0	1004.688	1025.574	1046.607	1.021	16.999	17.359	17.722
1	1060.729	1080.050	1099.410	1.018	17.206	17.539	17.873
2	1099.642	1117.677	1135.749	1.016	17.118	17.429	17.740
3	1046.184	1067.114	1088.205	1.020	15.439	15.799	16.163

Assignments | Filtered Assignments | Filtered Residuals | Details | Mass Ranges | AB Distribution | Repeat A Statistics | Repeat B Statistics

Repeat A Plot

More useful than tables of numbers, even with color coding, is to display spectra filtered by series and by A and B count. In the **Repeat A Plot** tab, the vertical slices taken through the **AB Distribution** table are displayed as spectra.

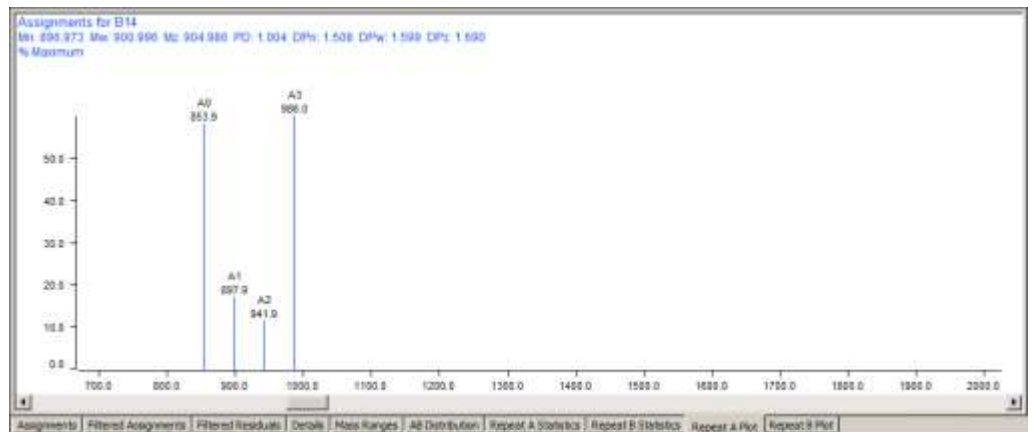

The screen shot above shows the A unit spectrum for a fixed B count of 14.



The scroll bar at the bottom of the plot is used to step through each B unit count. Clicking to the right increments the B count, clicking to the left decrements it. Each

COMPOLYMER ANALYSIS

plot is initially scaled to actual relative abundance on a 0 – 100 scale; left-click with the mouse in the plot window to autoscale it to the full height of the plot.

Repeat B Plot

Likewise, horizontal slices through the **AB Distribution** table can create plots of the B unit spectrum for fixed A counts. Similarly, the scroll bar is used to step through the A count increments. Two steps in A count are shown in the screen shots below. The first is the B spectrum for A0, the next is the B spectrum for A1.

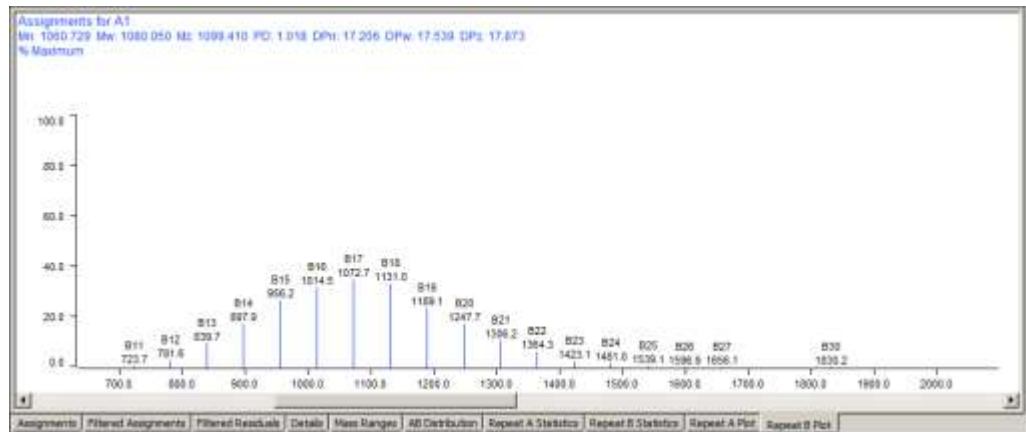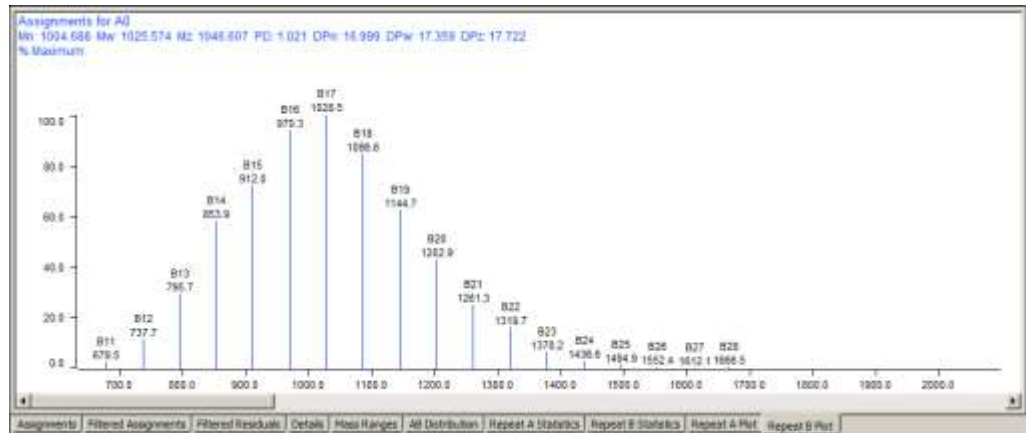

Here it can be seen that the B distribution is smooth for each of the A counts.

What was covered in this chapter

In this chapter we

- Described the definition of a copolymer used by Polymerix.
- Explained how to set processing parameters and define copolymer series.

COMPOLYMER ANALYSIS

- Presented the various result plots and tables and illustrated the utility of each one.

In the next chapter

In the next chapter we will describe the relative area tool.

The Relative Area Tool

Using the relative area tool

The Relative Area tool is just that, a simple tool that gives an estimate of the total relative abundance due to spectral features in selected parts of the spectrum.


Goals for this chapter

In this chapter, we will:

- Describe the **Relative Area** page
- Explain how to select spectral regions for relative area measurement

The Relative Area page

Like other Polymerix pages, the **Relative Area** page consists of three panels.

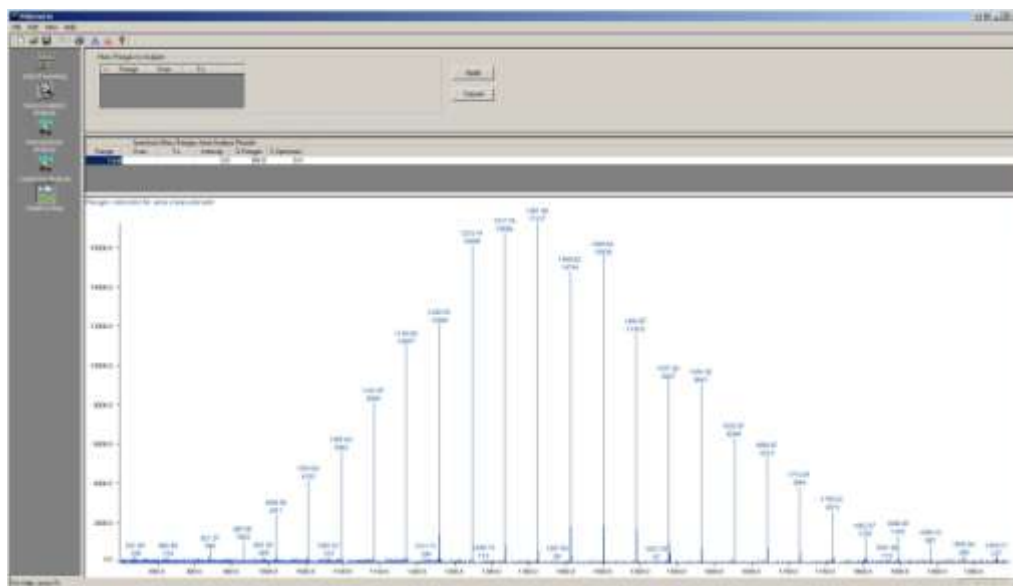


RELATIVE AREA

At the top of the page are parameters for specifying the relative area regions. In the middle is a pane containing a table of the relative area ranges and relative abundance percentages. Finally, the bottom pane displays the processed target spectrum with each area highlighted in color.

Specifying ranges

There are two ways to specify relative area ranges.

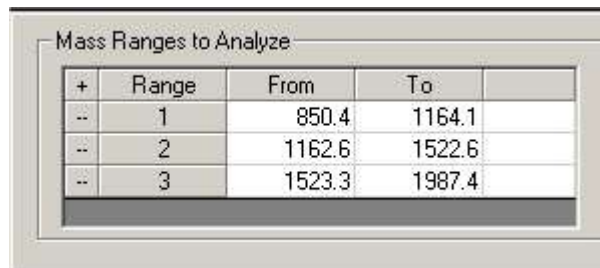

+	Range	From	To
--	1	850.4	1164.1
--	2	1162.6	1522.6
--	3	1523.3	1987.4

1. Click the **+** button in the upper left corner of the **Mass Ranges to Analyze** table. This adds a new row to the table. Type the m/z range start and end values. Continue until all ranges have been defined, and then click the **Apply** button.
2. Alternatively, you may right-click and drag in the spectrum window. Right-click at the beginning of the range, and then while holding the mouse button down, drag to the end of the range. Repeat this for each new range. The ranges will automatically be added to the table. If you wish, you may fine-tune the range setting by editing the table afterward. Click the **Apply** button to calculate the percentages.

Relative Area results

The relative areas and their percentages are displayed in the table in the middle pane.

Spectrum Mass Ranges Area Analysis Results					
Range	From	To	Intensity	% Ranges	% Spectrum
Total			264411.3	100.0	96.5
1	850.4	1164.1	66333.0	25.1	24.2
2	1162.6	1522.6	137556.4	52.0	50.2
3	1523.3	1987.4	60521.9	22.9	22.1

For each range, Polymerix computes the total integrated intensity, the percentage that intensity reflects of all of the defined ranges and of the spectrum as a whole (which may include area outside of defined ranges).

In the next chapter

In the next chapter, we will describe the autocorrelation analysis process and how it can assist in determining the composition of an unknown polymer. Autocorrelation analysis has been superseded by the functionality provided by the Find Homopolymer Series page, so the chapter can be skipped.

Autocorrelation

How to use autocorrelation to analyze unknown polymers

Polymerix is normally used to deconvolve polymer series and compute molecular weight distributions for materials of known composition. What can you do when you have an unknown polymer, or a polymer where only partial information is known?

Goals for this chapter

In this chapter, we will:

- Describe how to enable and access the **Autocorrelation** page
- Describe the autocorrelation algorithms and what the resulting spectra mean, and
- Show how to use the information in the autocorrelation and end group spectra to assist in determining repeat unit or end group composition.

Accessing the Autocorrelation page

The **Autocorrelation Analysis** page and its button are not displayed by default, because **Find Homopolymer Series** supersedes the functionality of autocorrelation analysis in most cases. To enable and show the button, select the **View** menu's **Autocorrelation...** item and then exit and restart Polymerix. Once Polymerix has been restarted there will be a new Autocorrelation icon not previously shown. Clicking on the icon will access the **Autocorrelation Analysis** page.

Summary of the algorithm

The autocorrelation spectrum

AUTOCORRELATION

Autocorrelation is a method by which all m/z differences between peaks in a mass spectrum can be computed. The peaks in an autocorrelation spectrum indicate a frequently-occurring m/z difference where the mass spectral peaks have high relative abundance.

The autocorrelation spectrum can be computed by a brute-force process: For every point in a spectrum, compute the m/z difference between it and every other point in the spectrum, and assign to each difference an intensity weighted by the intensity of the two points. Create a difference spectrum by summing together the differences computed for every point. For a large spectrum with many points, this can be a very compute-intensive process.

There is a simpler way, based on the Fourier transform. Without going into great detail, the autocorrelation spectrum can be computed simply by taking the Fourier transform of the original mass spectrum, convoluting it with itself, and then transforming the result back.

What does the autocorrelation spectrum look like? Take for example a singly-charged ethylene oxide (EO) homopolymer spectrum. Regardless of the polymer molecular weight, there will be a series of peaks separated by the EO monomer mass, or 44 Da. There may be more than one EO series present, possibly due to different end groups or adducts, but each of these series will have a spacing of 44 Da as well.

The autocorrelation spectrum will therefore have a strong peak at 44 Da. It will also have peaks of lower intensity at 88 Da, 132 Da, and higher multiples of 44 Da. These *overtones* are due to the fact that for every group of peaks with a difference of 44 Da, the first and third peaks, second and fourth peaks, and so on are separated by 88 Da, while the first and fourth, second and fifth, etc. has a 132 Da spacing, etc. However, there are fewer pairs of peaks with each higher multiple of the base separation, so these overtones form a progression of decreasing abundance in the autocorrelation spectrum.

In addition, if there is more than one homologous series, there will be autocorrelation peaks corresponding to the mass shifts between the series. For example, if a sample contains both protonated and sodiated series, there will be an autocorrelation peak at 22 Da, the difference in m/z between H (1 Da) and Na (23 Da). Likewise, if there are two or more end groups, there will be autocorrelation peaks corresponding to the mass *differences* between them.

Thus, the peaks in the autocorrelation spectrum give clues to the nature of the repeating unit chemical composition. If the original spectrum is singly-charged, then the peaks in the autocorrelation spectrum that are due to real chemical differences can be assigned a chemical composition.

The end group spectrum

Computing the end group spectrum

Once an autocorrelation spectrum has been computed, it is possible to get an estimate of the end groups by computing an end group spectrum. The process is as follows: From the autocorrelation spectrum, pick the base peak and assign it as the mass of the repeat unit (monomer). Then, for each peak in the experimental spectrum, subtract from it the highest integer multiple possible of the repeat mass. Insert these remainder mass values as peaks in the end group spectrum. If the repeat mass was correctly chosen, there will be clusters of peaks in the end group spectrum at positions corresponding to each of the total end group plus adduct masses.

To give a concrete example, suppose the experimental spectrum has peaks at 63, 85, 107, 129, 151, 173, 195, and 217 Da. The autocorrelation spectrum shows, among other differences, a strong peak at 44 Da. If we use 44 as a repeat unit, then we compute an end group spectrum according to the results shown in the table below:

Expt. m/z	Max # of 44 repeats	Max # x 44	Remainder
63	1	44	19
85	1	44	41
107	2	88	19
129	2	88	41
151	3	132	19
173	3	132	41
195	4	176	19
217	4	176	41

As should be clear from examining the table, the end group spectrum will have only two peaks, at 19 Da and 41 Da. The difference between these two is 22 Da, leading to the inference that there are two homologous series in the sample, one with proton adduct and the other with sodium adduct. So, the peaks at 19 Da must be due to the total of the end groups plus a proton, and likewise the peaks at 41 Da are due to the total of the end groups plus a sodium ion. Subtracting 1 Da from the 19Da peak and 23 Da from the 41 Da peak leaves 18 Da in both cases. This leads to the conclusion that the total end group is water, H₂O.

With this information and some knowledge of polymer chemistry, it seems likely that the sample is an ethylene oxide polymer with H and OH end groups.

A real example would undoubtedly be more complex and subject to errors in calculation of the repeat mass and remainder masses. In particular, if the total end group plus adduct mass is larger than the repeat mass, then the algorithm described above will subtract too many repeat units, leading to masses in the end group spectrum that are too small by one or more repeat masses. This will lead to nonsensical


AUTOCORRELATION

compositions. In such cases, you may need to add one or more repeat masses into the peak values to see if that makes more sense as a total end group mass.


Nevertheless, autocorrelation analysis can often give valuable insights into composition of an unknown material.

User interface and parameters


The autocorrelation and end group spectra are computed automatically from the process target spectrum. To see these spectra, click the **Autocorrelation Analysis** button on the **Workspace Control Panel** to switch to the **Autocorrelation Analysis** page.

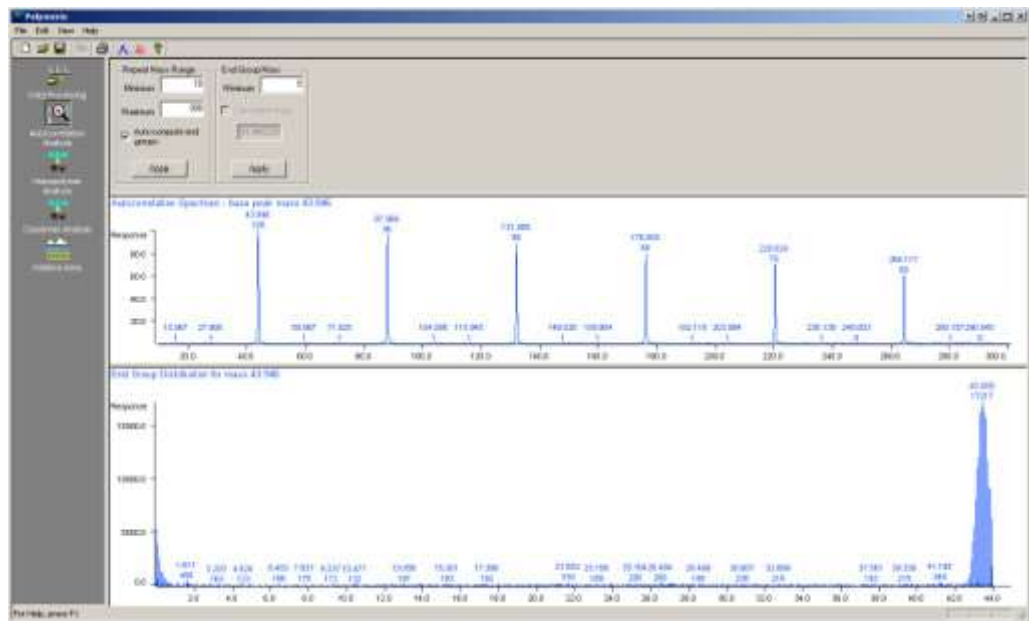

A magnified view of portions of these two spectra is shown below:

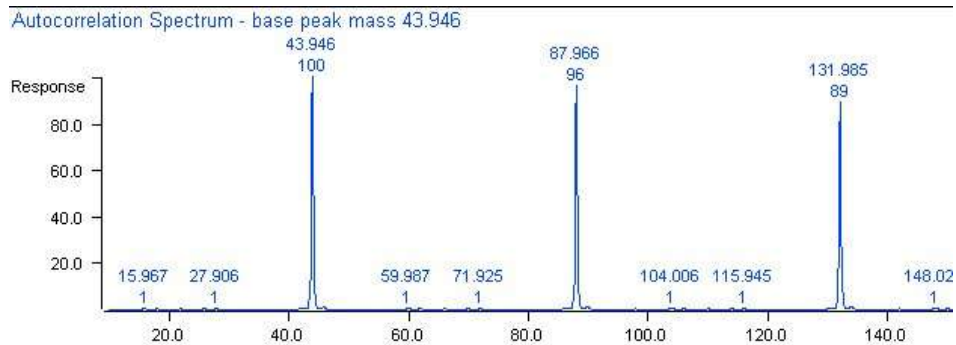

AUTOCORRELATION

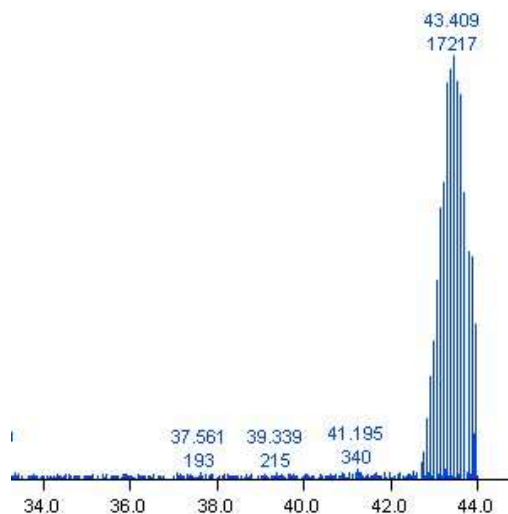

1. There are two sets of parameters used in the autocorrelation analysis. The first controls the computation of the autocorrelation spectrum itself, and whether the repeat mass determined automatically from it is used to compute the end group spectrum.

A dialog box titled "Repeat Mass Range" with the following controls:

- Minimum:
- Maximum:
- Auto-compute end groups
-

2. The default repeat mass range for the autocorrelation spectrum is from 10 to 300 Da. This will cover a large number of common monomers. If observation of the experimental spectrum gives evidence for a larger repeat mass, adjust these parameters, and then click the **Apply** button.
3. The second set of parameters controls how the end group spectrum is computed. By default, the mass of the maximum intensity peak in the autocorrelation spectrum is chosen as the repeat mass. In this case, the only end group parameter available is the minimum end group mass, which is usually set to zero. To set a different end group minimum mass, change the value, then click the **Apply** button.

AUTOCORRELATION




End Group Mass

Minimum

Use repeat mass:

- In the illustration above, the repeat mass value 43.946 Da was chosen, using the mass of the maximum abundance point in the autocorrelation spectrum. Applying this to compute the end group spectrum leads to a cluster of peaks near 43 Da.
- However, 43.946 Da does not correspond to any meaningful chemical formula for a monomeric repeat unit. A better, but still close value is 44.0262 Da, corresponding to the mass of the ethylene oxide monomer with formula C_2H_4O . Particularly in TOF spectra, peak tailing due to detector dead time can result in intensity-dependent variations in peak centroids. This leads to broadening and shifts in the autocorrelation spectrum. So, let's make the assumption that the actual repeat mass is 44.0262 Da and recompute the end group spectrum accordingly.
- In the **Repeat Mass Range** section, uncheck the **Auto-compute end groups** checkbox. In the **End Group Mass** section, check the **Use repeat mass** checkbox, and enter the value **44.0262** in the edit box. Click the **Apply** button. The end group spectrum will be recomputed, this time using 44.0262 Da as the repeat mass.

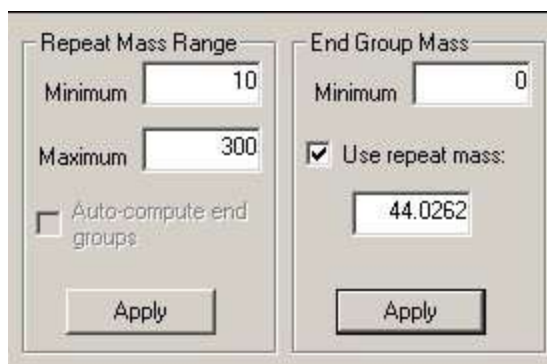

Repeat Mass Range

Minimum

Maximum

Auto-compute end groups

End Group Mass


Minimum

Use repeat mass:

- Now, as can be seen from the portion of the end group spectrum below, it consists of just a single peak at 41.000 Da. This implies a sodium adduct. Subtracting the mass of Na (22.990 Da) yields a

AUTOCORRELATION

remainder of 18.010 Da, very close to the exact mass for H₂O of 18.0105 Da.

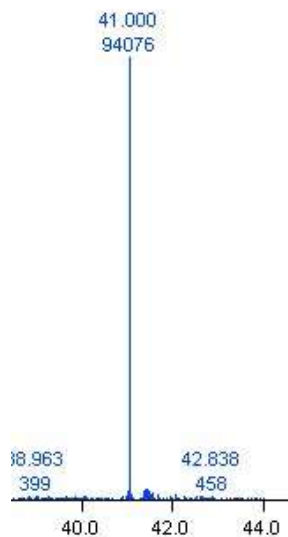

What was covered in this chapter

In this chapter, we

- Showed how to access the autocorrelation analysis page.
- Described the process for creating autocorrelation and end group spectra.
- Demonstrated how these spectra might be used to help identify an unknown polymer.

In the next chapter

the next chapter we will describe how to get all results out of Polymerix and into other documents, files, or printed reports and how to set some plot customization options.

Creating and Customizing Output

How to do get creative with Polymerix results

After spending a lot of time and thought setting up polymer analysis parameters, series, and analyzing data, it would be nice to be able to do something with it besides look at it on screen. In this chapter we will present all of the various options for getting information out of Polymerix and into your other documents, reports, and files.

Goals for this chapter


In this chapter, we will:

- Describe how to customize the appearance of Polymerix's plots
- Explain how to copy plots, spectral data, and tabular results to the Windows Clipboard.
- Show how to save and export Polymerix results

Polymerix options

There are few options that can be set to customize plot colors and labeling.


Plot labeling options

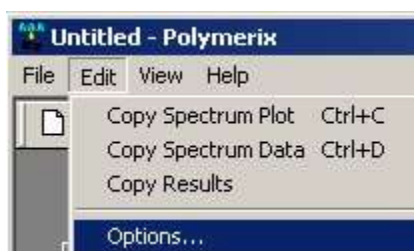

1. To set options, select the **Edit → Options...** menu command. This will display the **Polymerix Options** dialog box.




2. The **Display** tab allows you to specify the X-axis labeling preference for the chromatogram (TIC) plot as either scan number or retention time, and to set options for labeling of peaks in mass spectra. To include a label on the spectra, click the checkbox to enable it. Optionally change the number of decimals of precision in the label.

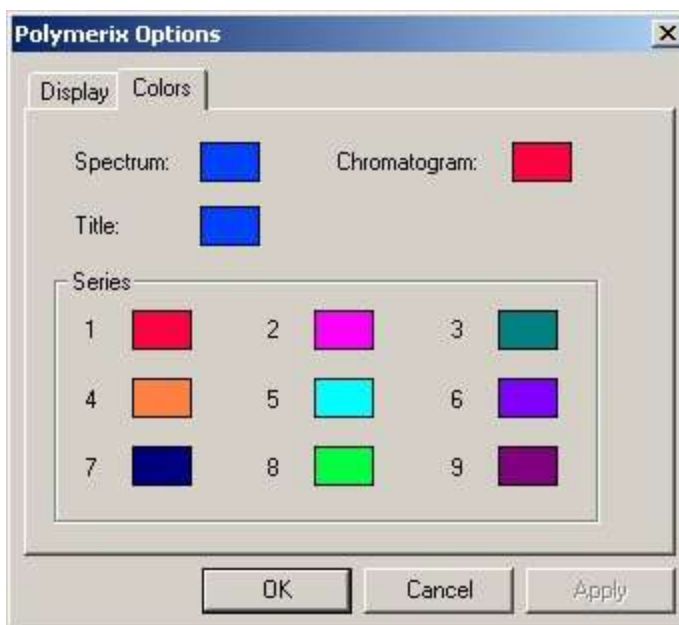

Plot color options

3. The **Colors** tab allows you to specify the colors used for the TIC, spectrum, plot titles, and assigned series. The colored rectangles to the right of each item are buttons. Clicking the button displays a **Color**

selection dialog box, from which you may choose a basic color or define a custom color.

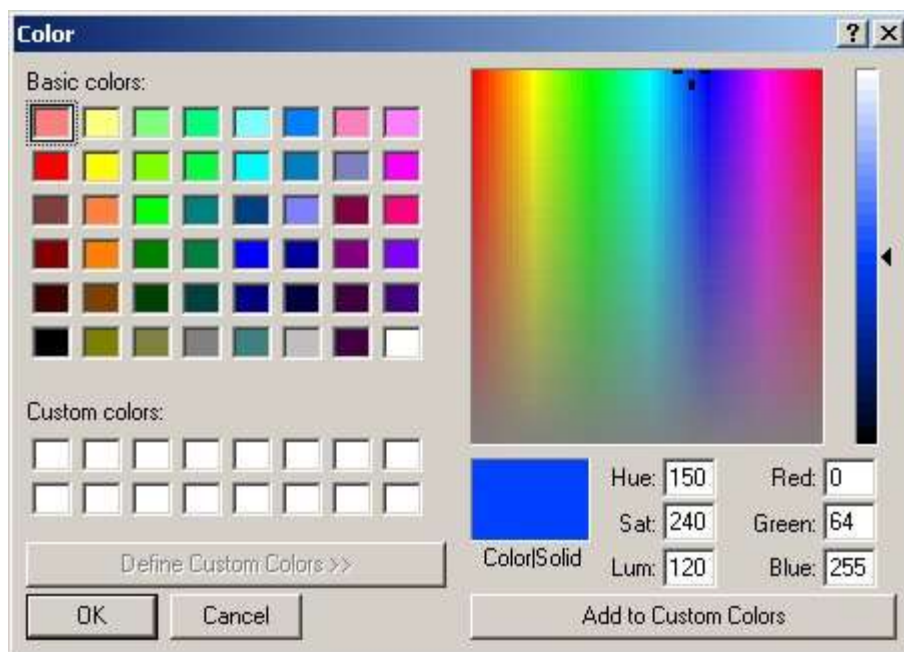

4. Click **OK** to close the **Polymerix Options** dialog box. The user interface will be updated immediately to reflect any changes.


Copying to the clipboard

The graphics in any of the spectrum panes may be copied to the Windows clipboard as a bitmap picture, or the spectrum details shown in that plot can be copied as a spectrum peak list.


Copying a spectrum

The currently-selected spectrum pane is always the source for a copy operation. The current spectrum pane is the one in which the last mouse click operation was performed. To copy the pane, do the following:

1. Left-click on the spectrum to select it. Zoom in if desired – the copied image will exactly as seen on screen.
2. To copy the plot graphics, select the **Edit → Copy Spectrum Plot** menu command (or press the **Ctrl + C** keys) to copy the spectrum image to the Clipboard. The graphics can then be pasted into another application, such as Microsoft Word or PowerPoint.

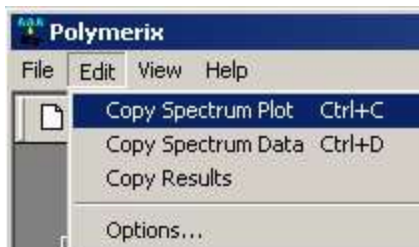


3. To copy the spectrum data as a peak list, select the **Edit → Copy Spectrum Data** menu command (or press the **Ctrl + D** keys). This places a list of the spectrum peaks on the Clipboard. These may be pasted into Microsoft Word or Excel as a table.

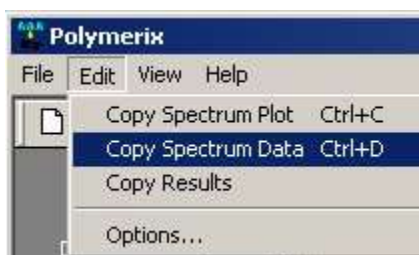

Copying results


The series definitions, statistics, and detailed assignments can be copied to the clipboard as a formatted table.

1. Click on the **Details** tab to bring it to focus..
2. Select the **Edit → Copy Results** menu command to place a formatted table on the Clipboard. This may be pasted into Microsoft Word or Excel as a table.

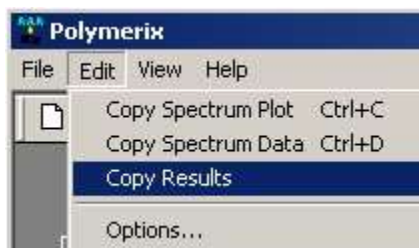

Saving and exporting results

Polymerix results and the input (unprocessed) spectrum can also be saved as ASCII files for import into Microsoft Excel or other software.

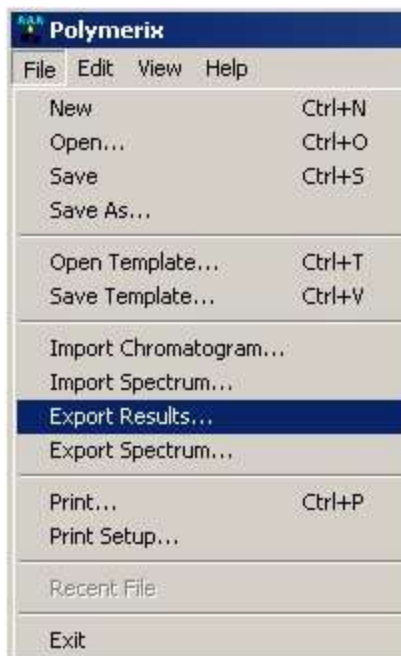

Exporting results

1. To save Polymerix results, select the **File → Export Results...** menu command. A file dialog will appear to allow the choice of file name and location.



Exporting the input spectrum


42. To export the input spectrum as an ASCII peak list, select the **File → Export Spectrum...** menu command. A file dialog will be displayed to allow file name and location selection. The spectrum will be saved as a tab delimited file, with one (m/z, intensity) pair per line.

Printing

Polymerix produces a preformatted printed report as output. There are a number of choices for the items to include in the report.



1. Select the **File** → **Print...** menu command. An enhanced **Print** dialog will be displayed.

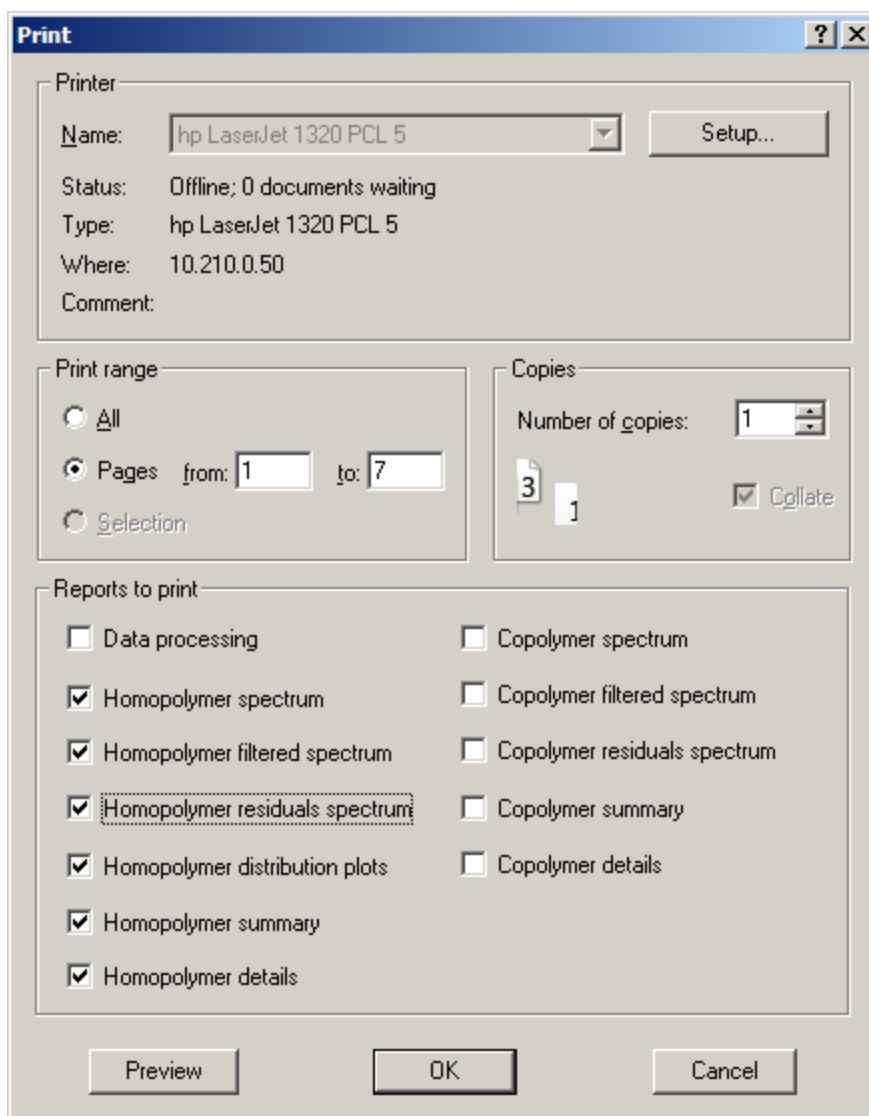

2. In addition to the normal printing options at the top of the dialog, the lower part of the dialog contains Polymerix-specific printing options. Selecting an option will create one or more pages in the output containing the chosen results. These are:
 - **Data Processing:** Produces a page containing data file header information, the raw input spectrum plot, and the processed target spectrum plot. Note that the plots will be displayed in the report at their current on-screen zoom state.
 - **Homopolymer spectrum:** Produces a page containing the plot shown on the **Assignments** tab of the **Homopolymer Analysis** page.

- **Homopolymer filtered spectrum:** Produces a page containing the plot shown on the **Filtered Assignments** tab of the **Homopolymer Analysis** page.
- **Homopolymer residuals spectrum:** Produces a page containing the plot shown on the **Filtered Residuals** tab of the **Homopolymer Analysis** page.
- **Homopolymer distribution plots:** Produces multiple pages, one containing a plot of the total series spectrum, and followed by additional pages, each containing an individual series spectrum plot.
- **Homopolymer summary:** Produces a page containing the **Homopolymer Results Summary**, as seen in the middle pane of the **Homopolymer Analysis** page.
- **Homopolymer details:** Produces one or mages of the tabulated homopolymer assignment details, as seen on the **Details** tab.
- **Copolymer spectrum:** Produces a one page plot of the copolymer **Assignments** spectrum.
- **Copolymer filtered spectrum:** Produces a one-page plot containing the copolymer **Filtered Assignments** spectrum.
- **Copolymer residuals spectrum:** Produces a one-page plot containing the copolymer **Filtered Residuals** spectrum.
- **Copolymer summary:** Produces a page containing the **Copolymer Results Summary** as seen in the middle pane of the **Copolymer Analysis** page.
- **Copolymer details:** Produces one or more pages tabulating the copolymer assignment details, sorted by series, and by increasing B count within A count.
- **Autocorrelation results:** This check box is only shown if autocorrelation has been enabled by use of the **View** menu's **Autocorrelation...** item. Produces a page with the autocorrelation and end group distribution plots, at their current on-screen zoom level.

An example homopolymer analysis report is included in the Appendix.

What was covered in this chapter

In this chapter, we

- Described how to customize plot display options
- Showed how to copy plots, spectral data, and results to the Clipboard
- Explained how to save and export Polymerix results.

In the next chapter

There is no next chapter... you've reached the end of the manual! It's time to go to work and analyze some polymers.



Release Notes

Version 3.0

Major new features

The major new features improve polymer analysis of centroided data with high mass accuracy.

- The new **centroid cluster matching** feature matches individual peaks in each isotope cluster for a polymer series. This allows matching of centroided data without deisotoping. For data with high mass accuracy, the new matching feature allows use of much smaller match tolerance values. Use of smaller match tolerance values can reduce or eliminate mass overlaps between different series.
- The **Find Homopolymer Series** feature can help a user select chemical compositions for a homopolymer series from lists based upon pairs of reference peaks in the spectrum being analyzed. Selecting compositions using **Find Homopolymer Series** is often easier and quicker than manually setting up polymer series definitions in **Homopolymer Series Setup**. **Find Homopolymer Series** uses centroid cluster matching.

Updates, bug fixes, and new features

- User interaction is essentially unchanged from Version 2.0. Some superfluous mouse clicks have been eliminated.
- Added support for multiply charged data with mixed adducts.
- Added a Filtered Residuals tab and plot to polymer analysis results.
- Added a Mass Ranges tab and table to homopolymer analysis results.
- Added error metrics to **Homopolymer Details** and **Copolymer Details**.
- Added percent of series molecular weight metric for series components.
- Changed the position of series annotations in plots so that their x positions always correspond to calculated monoisotopic peaks for the series.
- Depreciated the Autocorrelation tool, but added a menu item to the **View** menu so that it is still accessible.

RELEASE NOTES

- Account for electron mass when computing mass ranges for use in matching. Added a **Positive charge** checkbox to both **Homopolymer Setup** and **Copolymer Setup**.
- Fixed a problem with the calculations of intensity weighted mass averages. Negative intensity values are now ignored in these computations.
- New data file format. Data files created by earlier versions of Polymerix can be loaded by Polymerix version 3.0, but data files created by Polymerix version 3.0 cannot be used with earlier versions of Polymerix.
- Fixed problem with multiple charge states combined with a loss.
- Added % RA threshold to MS peak processing.
- Added the **Reset Mass Ranges** checkbox to the **Data Processing** page.
- Improved peak detection.

Version 2.0

Updates, bug fixes, and new features

- User interaction is essentially unchanged from Version 1.0.
- Moved copolymer processing from Polymerix V2 beta to the 2.0 Release version including full copolymer reporting.
- Implemented improved mass spectral peak detection.
- Fixed known printing and report generation problems, including missing lines, blank pages, incorrect page count when switching between landscape/portrait, incorrect sizing of collate icons, blank copolymer assignments spectrum.
- Fixed bug in copolymer details display which used to manifest when copolymer series other than the last were disabled.
- Result summaries now include entities specified by mass and consolidate shared formulas.
- Polymer statistics now have zero values when intensities don't match.
- Fixed heap corruption bug in Homopolymer spectrum view when attempting to color unassigned masses.
- Fixed homopolymer assignment bug where assignments were generated for masses of zero.
- Added "Filtered Assignments" spectrum view.
- Added homopolymer distribution plots.
- Added intensity weighted mass metrics.
- Changed MS data access to use ordinary DLL instead of COM DLL. (If low-level data access tools supplied by the vendor use COM, this is not affected by this change. Only Sierra's high-level data access is changed).
- Updated installation script to use new DLLs.

RELEASE NOTES

- Changed licensing to use a new scheme that does not depend on MAC address. Two methods are provided:
 - “fixed” license (licensed to a single PC) or
 - “floating” license using a designated portable USB device (licensed up to 4 PCs). Copies using a floating license must have the USB device installed at all times when the software is in use. The software will exit if the USB device is removed.



Example Report

An example of a homopolymer analysis report

On the following pages is an example of a full report for a PEG homopolymer analysis. On the **Print** dialog, the **Data Processing** and all **Homopolymer** report options were selected.


RELEASE NOTES

Polymerix Processing Report
Data Processing Results
Printed: Wed Mar 06 2013 15:42:07


Page 1 of 8

Data file: C:\Polymerix\peg1500_cal.raw
Processing mass range 0.0 to 2402.0
Peak width TOF, 0.400 Da at 1000.000 m/z
Peak detected, area intensity
Baseline subtraction width 500.0 Da


C:\2.0e+6\erix\peg1500_cal.raw Scans 1 to 5

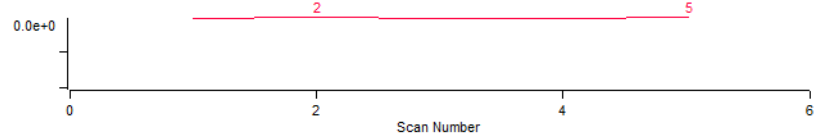

peg1500_cal.raw RT: 0.04:0.19 Spectrum: 1:5 111919 Points

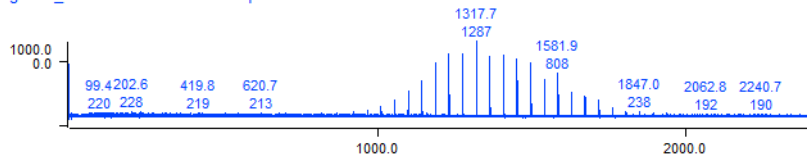


Detect(0.4 Da @ 1000, 500) 8590 Peaks Mass Range 0.00 to 2402.00

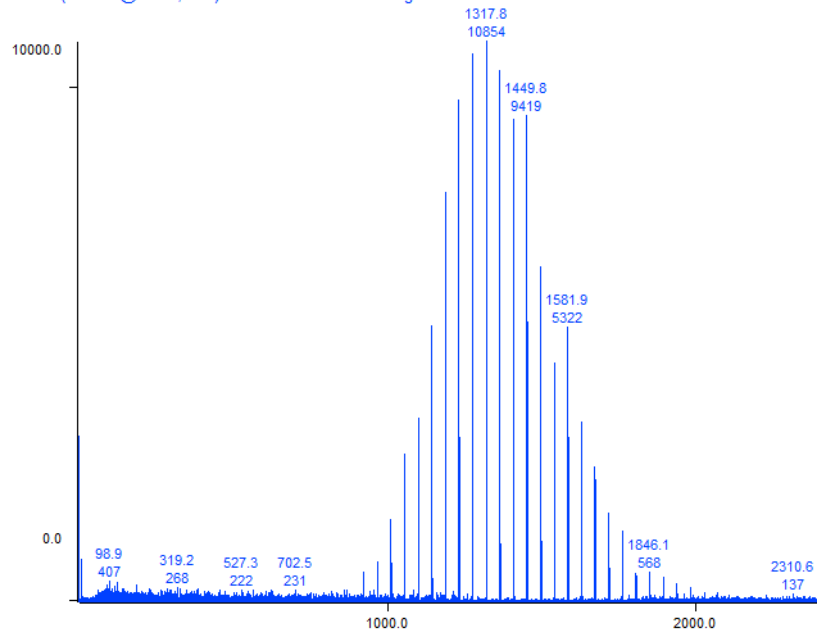


Data file: C:\Polymerix\peg1500_cal.raw
Monoisotopic mass mode
Match tolerance +/- 0.100 Da

Homopolymer Assignments

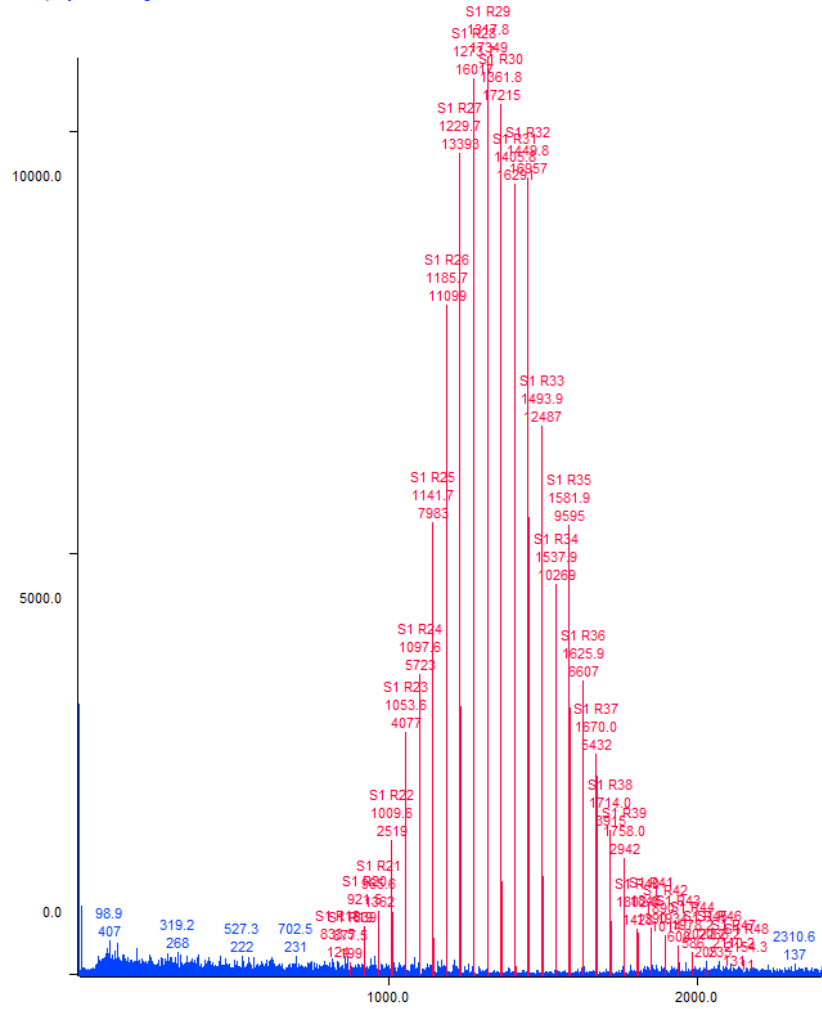


Data file: C:\Polymerix\peg1500_cal.raw
Monoisotopic mass mode
Match tolerance +/- 0.100 Da

Filtered Homopolymer Assignments

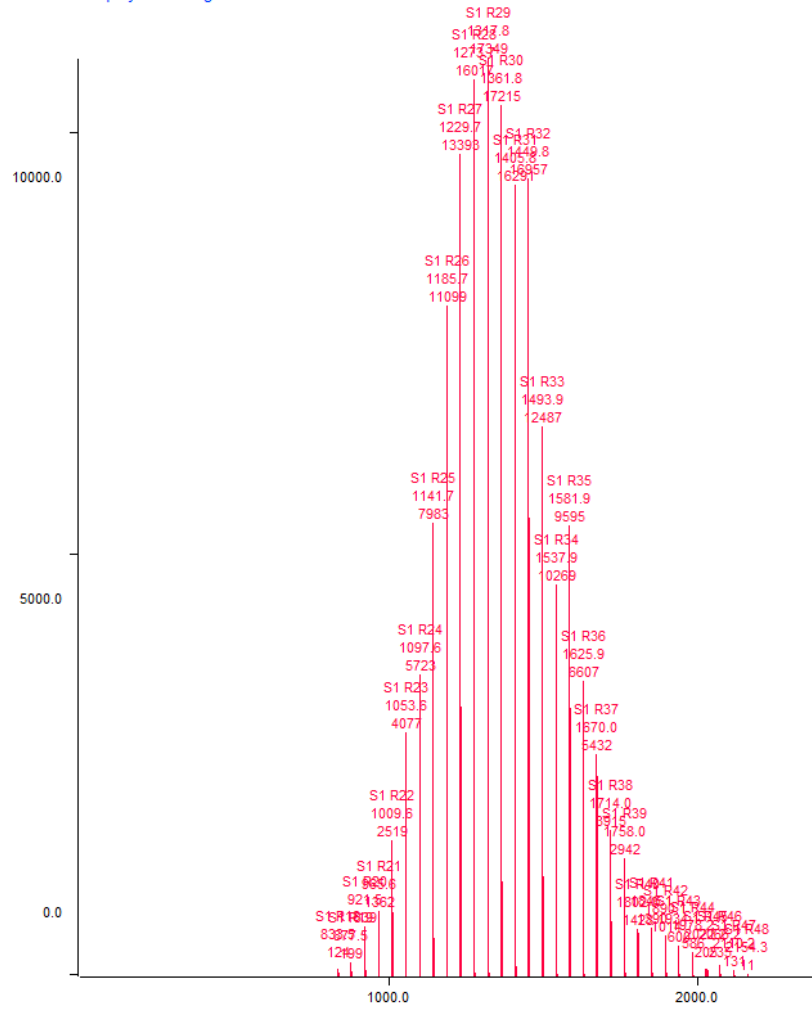


Data file: C:\Polymerix\peg1500_cal.raw
Monoisotopic mass mode
Match tolerance +/- 0.100 Da

Filtered Homopolymer Residuals

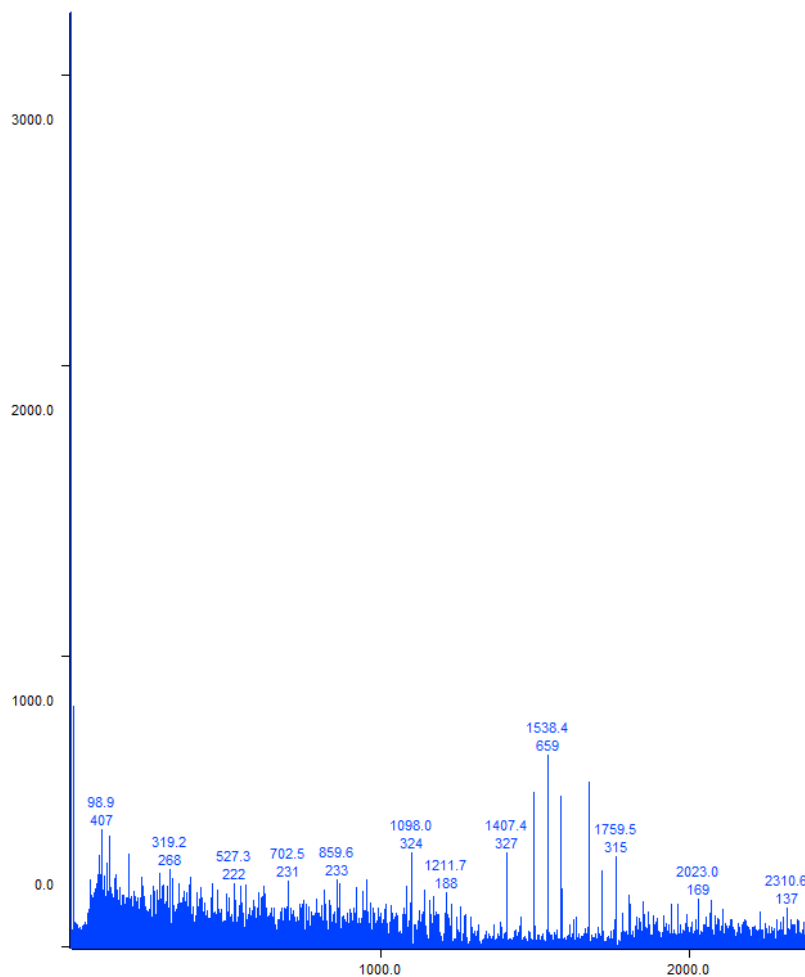


Data file: C:\Polymerix\peg1500_cal.raw
Monoisotopic mass mode
Match tolerance +/- 0.100 Da

Assignments for Total/Average
% Maximum

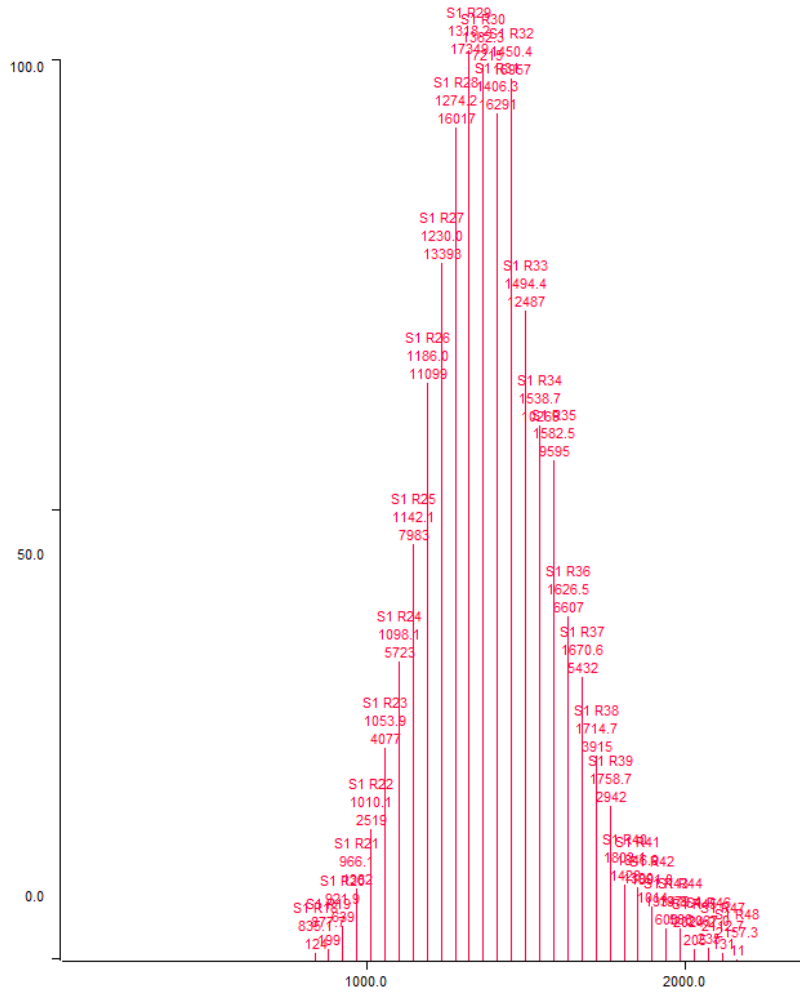

Data file: C:\Polymerix\peg1500_cal.raw
Monoisotopic mass mode
Match tolerance +/- 0.100 Da

Assignments for S1
% Maximum

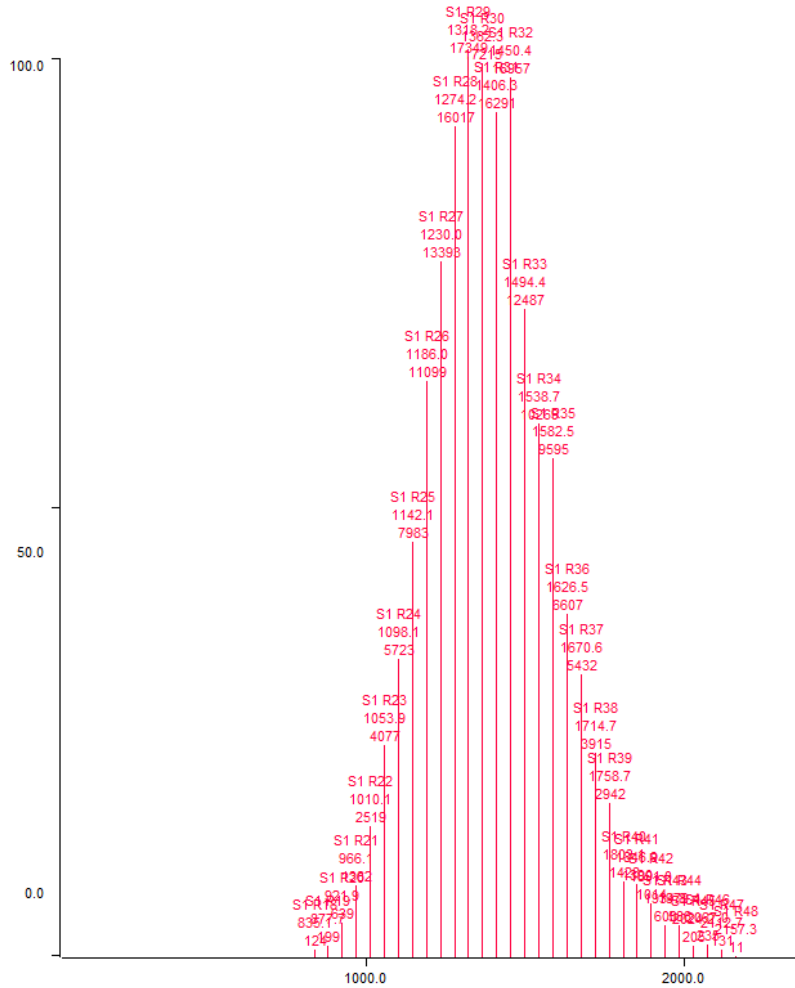

RELEASE NOTES

Polymerix Processing Report
 Homopolymer Analysis Results
 Printed: Wed Mar 06 2013 15:42:07

Page 7 of 8

Data file: C:\Polymerix\peg1500_cal.raw
 Monoisotopic mass mode
 Match tolerance +/- 0.100 Da

Homopolymer Results Summary

Series Label	Mn	Mr	Mz	PD	DPn	DPv	DPz	Percent Series	Percent Spectrum
Total/Average	1362.742	1390.981	1419.490	1.021	30.544	31.185	31.833	100.00	34.83
S1	1362.742	1390.981	1419.490	1.021	30.544	31.185	31.833	100.00	34.83

Series Label	Alpha End Group	Repeat	Omega End Group
Total/Average	H2O	C2H4O	
S1	H2O	C2H4O	

Series Label	Charge State	Adduct	Adduct Charge	Loss	Series Formula
Total/Average	1	Na	1		H2O [C2H4O] _n + Na
S1	1	Na	1		H2O [C2H4O] _n + Na

Homopolymer Analysis Details

Series Label: S1									
Repeat Count	Percent Series	% Series Weight	Percent Spectrum	Computed Mono. m/z	Found Mono. m/z	Error mDa	Found Intensity		
18	0.07	0.04	0.02	833.472	0.000	0.000	124.29		
19	0.11	0.07	0.04	877.498	877.486	-11.547	199.43		
20	0.34	0.22	0.12	921.524	921.572	48.418	639.41		
21	0.73	0.50	0.25	965.550	965.561	10.767	1361.97		
22	1.34	0.97	0.47	1009.577	1009.608	31.269	2518.73		
23	2.17	1.64	0.76	1053.603	1053.608	5.570	4076.74		
24	3.05	2.40	1.06	1097.629	1097.625	-4.319	5723.39		
25	4.25	3.49	1.48	1141.655	1141.654	-1.438	7983.43		
26	5.91	5.04	2.06	1185.681	1185.684	2.684	11098.52		
27	7.13	6.32	2.48	1229.708	1229.710	2.653	13392.53		
28	8.53	7.83	2.97	1273.734	1273.741	6.968	16017.27		
29	9.24	8.78	3.22	1317.760	1317.761	0.552	17349.32		
30	9.17	9.01	3.19	1361.786	1361.795	9.234	17214.57		
31	8.67	8.81	3.02	1405.812	1405.817	4.322	16291.07		
32	9.03	9.46	3.14	1449.839	1449.844	5.413	16957.00		
33	6.65	7.18	2.32	1493.865	1493.867	2.578	12487.15		
34	5.47	6.08	1.90	1537.891	1537.893	1.436	10269.04		
35	5.11	5.85	1.78	1581.917	1581.914	-3.718	9598.10		
36	3.52	4.14	1.23	1625.944	1625.961	37.130	6606.95		
37	2.89	3.50	1.01	1669.970	1669.976	5.817	5431.56		
38	2.08	2.59	0.73	1713.996	1713.995	-0.870	3914.74		
39	1.57	2.00	0.55	1758.022	1758.043	21.251	2942.41		
40	0.76	0.99	0.26	1802.048	1802.108	59.407	1427.70		
41	0.74	0.99	0.26	1846.075	1846.059	-16.036	1389.92		
42	0.54	0.74	0.19	1890.101	1890.118	17.454	1013.69		
43	0.32	0.45	0.11	1934.127	1934.105	-22.173	600.23		
44	0.31	0.45	0.11	1978.153	1978.118	-34.847	585.97		
45	0.11	0.16	0.04	2022.179	2022.100	-79.001	205.46		
46	0.13	0.19	0.04	2066.206	2066.187	-18.865	234.87		
47	0.07	0.11	0.02	2110.232	0.000	0.000	131.16		
48	0.01	0.01	0.00	2154.258	0.000	0.000	10.70		
Repeat Count	Percent Series	% Series Weight	Percent Spectrum	Computed Cluster m/z	Found Cluster m/z	Error mDa	Found Intensity		
18	0.07	0.04	0.02	833.938	835.106	1167.264	124.29		
19	0.11	0.07	0.04	877.989	877.713	-276.026	199.43		
20	0.34	0.22	0.12	922.039	921.867	-172.183	639.41		
21	0.73	0.50	0.25	966.089	966.063	-26.874	1361.97		
22	1.34	0.97	0.47	1010.139	1010.052	-87.239	2518.73		
23	2.17	1.64	0.76	1054.189	1053.946	-243.443	4076.74		
24	3.05	2.40	1.06	1098.239	1098.085	-154.035	5723.39		

RELEASE NOTES

Polymerix Processing Report
Homopolymer Analysis Results
Printed: Wed Mar 06 2013 15:42:07

Page 8 of 8

25	4.25	3.49	1.48	1142.288	1142.052	-236.025	7983.43
26	5.91	5.04	2.06	1186.337	1186.023	-313.878	11098.52
27	7.13	6.32	2.48	1230.386	1230.018	-367.631	13392.53
28	8.53	7.83	2.97	1274.435	1274.151	-284.208	16017.27
29	9.24	8.78	3.22	1318.518	1318.206	-312.340	17349.32
30	9.17	9.01	3.19	1362.569	1362.279	-290.354	17214.57
31	8.67	8.81	3.02	1406.620	1406.317	-303.756	16291.07
32	9.03	9.46	3.14	1450.671	1450.418	-253.130	16957.00
33	6.65	7.18	2.32	1494.722	1494.443	-278.712	12487.15
34	5.47	6.08	1.90	1538.773	1538.653	-119.716	10269.04
35	5.11	5.85	1.78	1582.823	1582.490	-332.725	9595.10
36	3.52	4.14	1.23	1626.873	1626.529	-344.434	6606.95
37	2.89	3.50	1.01	1670.924	1670.589	-334.780	5431.56
38	2.08	2.59	0.73	1714.974	1714.741	-232.861	3914.74
39	1.57	2.00	0.55	1759.024	1758.697	-327.147	2942.41
40	0.76	0.99	0.26	1803.073	1803.079	5.266	1427.70
41	0.74	0.99	0.26	1847.123	1846.917	-206.618	1389.92
42	0.54	0.74	0.19	1891.203	1891.022	-180.390	1013.69
43	0.32	0.45	0.11	1935.254	1934.763	-491.404	600.23
44	0.31	0.45	0.11	1979.305	1979.412	107.008	585.97
45	0.11	0.16	0.04	2023.357	2024.193	836.846	205.46
46	0.13	0.19	0.04	2067.408	2066.971	-436.807	234.87
47	0.07	0.11	0.02	2111.459	2112.748	1289.486	131.16
48	0.01	0.01	0.00	2155.510	2157.319	1808.680	10.70